



Instruction Sheet

ULT Plus Series Battery Replacement Low Battery Light On

See page 4 for applicable models
October 23, 2014

Power off the unit in this sequence:

- First, power off “battery” switch. This is located on the front at the bottom of the unit below the door in behind a small pull out panel (see photo below).

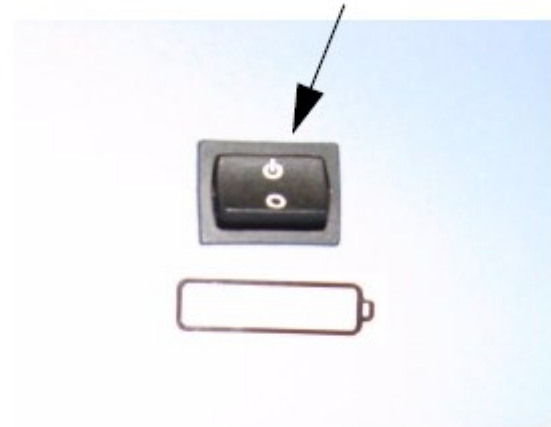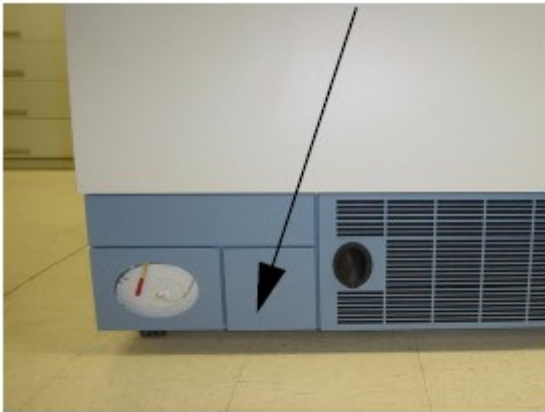

Front Panel, Battery Enable Switch

- Next, power off the “key” switch on the front of the display.



- Last, power off the “power/breaker” switch on the back of the unit below.



Remove the six screws that secure the lower left side panel. See photo below. Set the panel aside.



Locate the battery on the right side of the relay enclosure below.



Battery location (above)

- Disconnect the two red and black leads from the battery.
- Remove the screws securing the battery. Discard the battery, according to proper disposal procedures.
- Install new battery. Connect the wires = **Red** is the positive lead and black is the negative.
- Reinstall panel and restart in this order Battery Switch on, Breaker Switch on, Key switch on.

IMPORTANT: Allow 4-6 days for the battery to fully charge. Monitor unit for proper operation.

This process is applicable to the following models:

Revco Plus -86C Freezers; All Sizes & Voltages

Ultima Plus

Value Plus

Elite Plus

ULT1786-4-6-10

ULT2186-4-6-10

ULT2586-4-6-10