


TSX 505

Slider Basket – Installation Instructions

Description here: How to install Slider Baskets in a TSX505 refrigerator

TSX505
Revision Date: June 20, 2018
Created by: Wayne Anders

1. Remove all cold storage products from refrigerator.
2. Open the door and unplug the power cord from the wall and the power receptacle on the back of the refrigerator.
3. Carefully remove shelves and clips that you will be replacing with the slider baskets.
4. Decide the height you want your slider basket and identify the 4 Philips head screws you will need to remove. Note: The slider basket can be installed with either the screw holes on the bottom or top depending on your preferred storage configuration. When installing the bottom shelf, the bracket should be installed so that the screw hole is below the bracket.
5. Using a Philips head screwdriver unscrew the 4 screws.
6. Place the slider bar in the refrigerator aligning the screw holes on the cabinet and the slider bracket. Note that regardless of the top or bottom installation the stopper on slider bracket must face the rear as shown.
7. Tighten all the screws to secure the bracket to the cabinet.

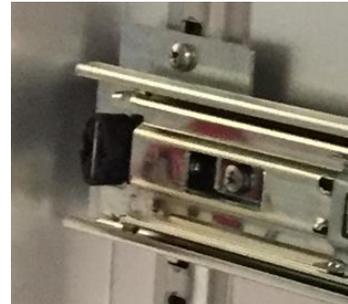

8. Repeat the slider bracket installation steps on the opposite side of the cabinet.
9. Install the basket by aligning and inserting the basket into the slider bracket. Note: The basket will only fit into the bracket with the open end of the basket rail facing the rear of the refrigerator.
10. Once installed, if you need to remove a basket from the slider bracket simply pull the basket out as far as possible. This will expose a plastic lever on the inside of each basket rail. Push these levers one up and one down to release the basket.
11. Restore power to the refrigerator by installing the power cord.
12. Wait until the chamber has reached set point, before loading any products.

