



Setting Up Qtegra for ICP-OES on a New Windows 10 PC

Backing Up Qtegra From Old Computer	2
Capturing Qtegra Configurations	2
New Instrument License Seat	5
Setting Up the New Computer.....	6
Viewing File Extensions.....	6
Network Setup	6
Windows Defender and Firewall.....	9
Create a System Restore Point	9
Installing Qtegra on New Computer	10
Installing Backed Up Files.....	10
Create Configurations	11
Restore Plugin Settings	11

Backing Up Qtegra From Old Computer

Back up the entire **Qtegra** folder located in the following location:
C:\ProgramData\Thermo on the old computer.

Capturing Qtegra Configurations

- 1) Open  Configurator
- 2) Navigate to  Experiment Configurator
- 3) Record the list of configurations

Note: One easy way to do this is to take a screen capture as in [Figure 1](#). It is helpful to expand the configurations before taking the screen capture so that the plugins contained within each configuration can be noted.

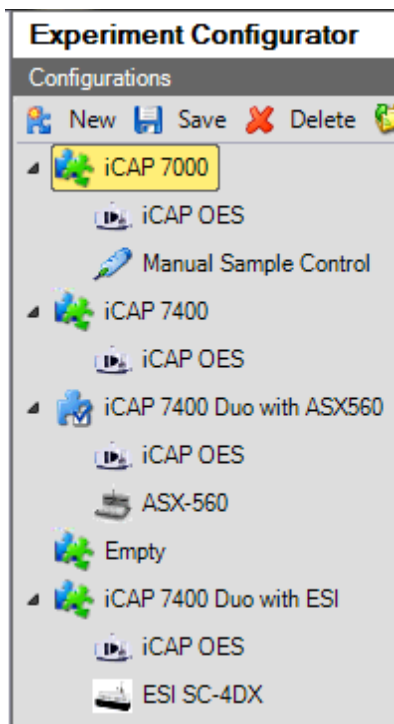


Figure 1: Experiment Configurations

- 4) Find the active configuration (evident by black checkmark in icon) as demonstrated in [Figure 2](#).

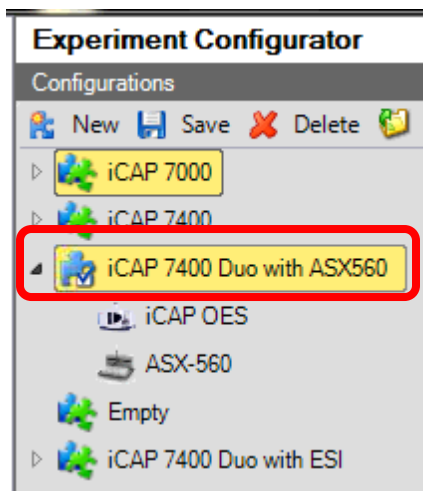


Figure 2: Active Configuration

- 5) Right-click on iCAPQ plugin and select Edit settings as shown in [Figure 3](#).

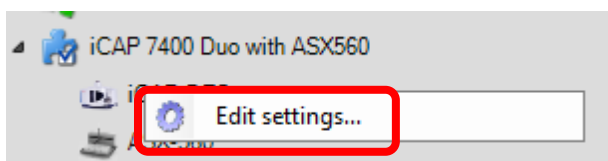


Figure 3: Edit the plugin settings

- 6) Note the Settings for that particular plugin. One easy way to do this is to take a screen capture of the settings for each plugin like that seen in [Figure 4](#).

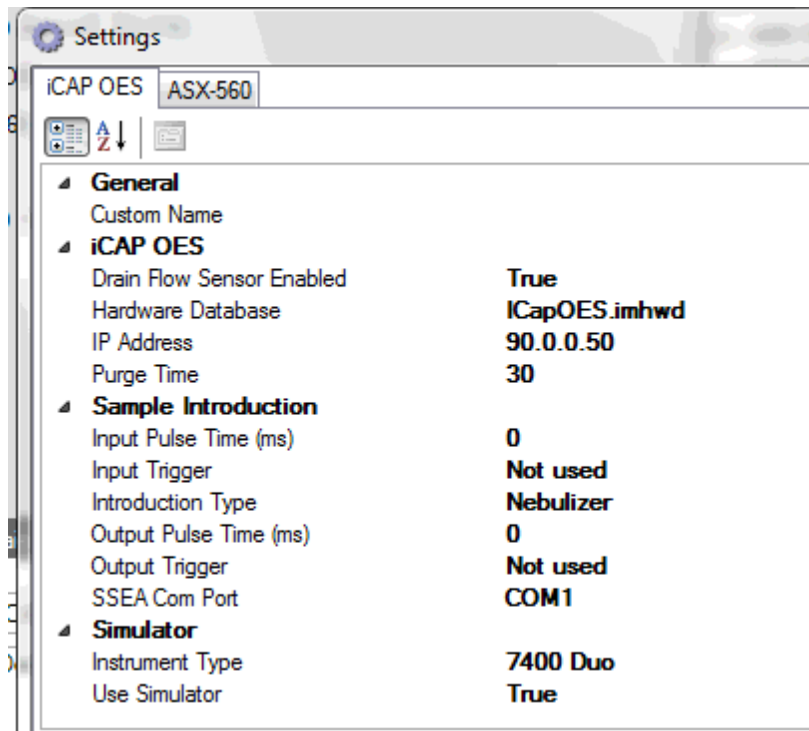


Figure 4: iCAP OES plugin settings

- 7) If you have additional plugins in that configuration, they will appear as an additional tab as can be seen with the ASX-560 plugin in Figure 5.

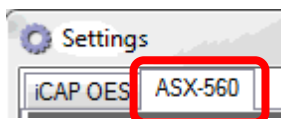


Figure 5: Plugin tab

Click on the additional plugin tab(s), if applicable to your configuration, and note their Setting(s) as demonstrated in Figure 6.

Note: Some plugins, such as ESI, will have multiple menus you will need to capture settings from.

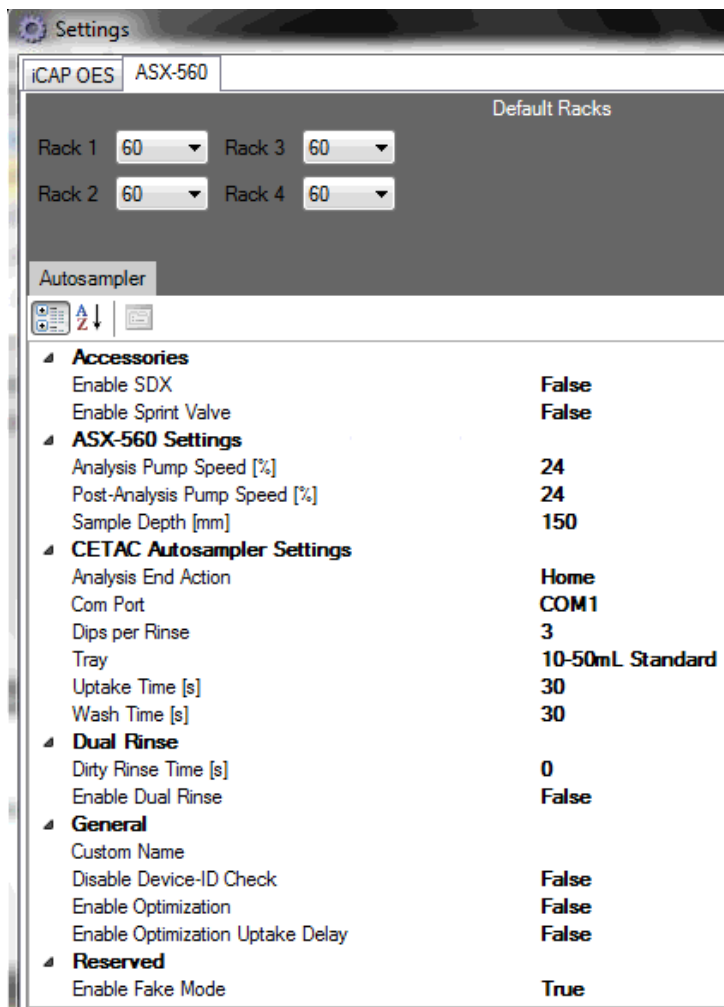


Figure 6: ASX-560 plugin settings

Note: If you use multiple configurations in your day-to-day analysis, you will need to note the Settings of any additional plugins contained in those configurations as well.




New Instrument License Seat

When Qtegra is licensed on a PC, the license key is bound to a hardware ID created from the computer component ID's. The hardware ID is generated from CPU, HDD and motherboard ID's. If one of these components has been changed or you seek to install Qtegra on a new PC, a new seat of the license for instrument control must be requested. This can be done by contacting: Qtegra.license.support@thermofisher.com. In the email you will need to provide your Qtegra license key and state that you would like another seat to be added to the license for instrument control.

Setting Up the New Computer

Viewing File Extensions

By default the “Program Data” folder is hidden. Please do the following to make it viewable:


- 1) Click on File Explorer  in the Windows taskbar. If File Explorer is not present there, click on Search  and then type in “File Explorer”. 
- 2) Click on the View tab




- 3) Check the Hidden items checkbox in the Show/hide section of the View ribbon



- 4) The ProgramData folder will now be visible when navigating to the C: drive

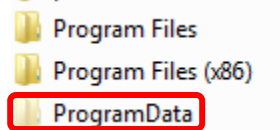


Network Setup

- 1) Click Search  and then type in “control panel” in the search box




- 2) An entry entitled **Control Panel** will appear. Click on this entry.

Note: You can right-click and select **Pin to Start** or **Pin to taskbar** to provide easy access to the Control Panel in the future.

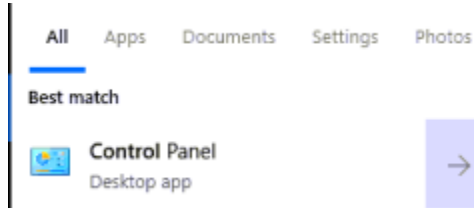


- 3) Click on **Network and Internet**

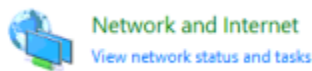

- 4) Click on **Network and Sharing Center**




- 5) Click on **Change adapter settings**

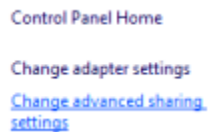


- 6) <Right-click> on the adapter your instrument is connected to and select **Properties**


Note: The name of the network connection may differ on your computer to the one referenced in this example

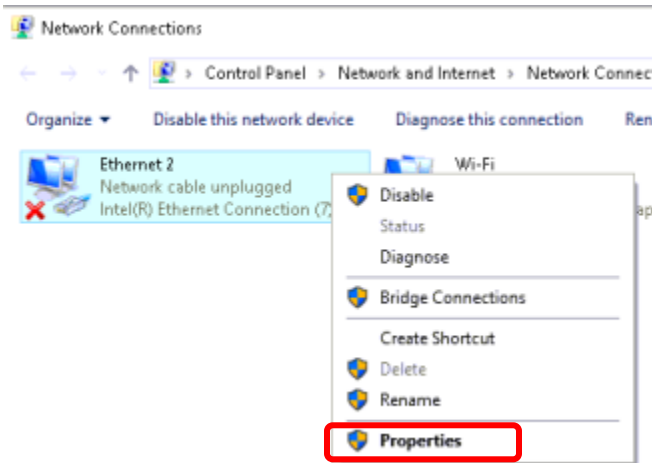


- 7) Deselect all of the checkboxes except for **Internet Protocol Version 4 (TCP/IPv4)**

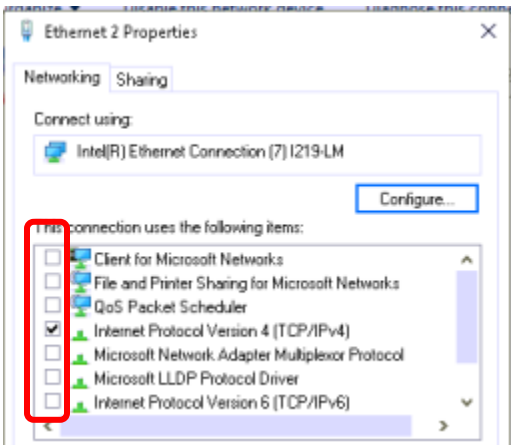

- 8) Select **Internet Protocol Version 4 (TCP/IPv4)** and click on **Properties**

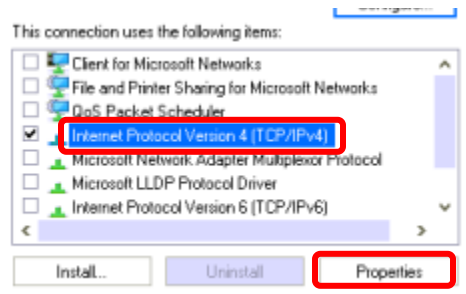

- 9) Click on the radio button **Use the following IP address** and enter into the **IP address** field: “90.0.0.1” and the **Subnet mask** field: “255.255.255.0” and then click **OK**

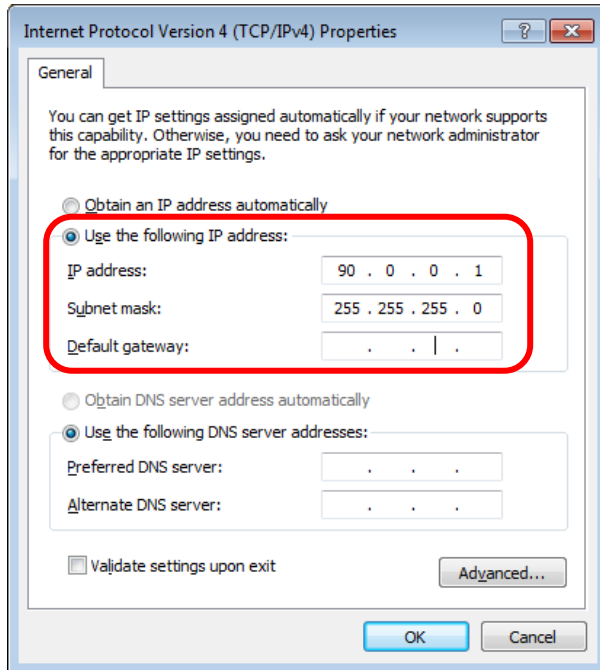


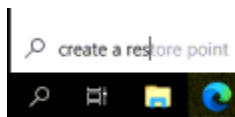
Windows Defender and Firewall

If the computer will not be networked, it is advisable to permanently turn OFF Windows Firewall and Windows Defender by following the attached tech note entitled “**Disabling Windows Defender and Firewall in Windows 10**”


If you the computer will be networked for example for access to the world wide web and firewall and/or antivirus will be enabled as a safeguard measure, exclusions will need to made in Security applications such as Firewall, Host Intrusion, and/or Antivirus. Please reference the attached tech note entitled “**Setting Up Exclusions for Qtegra for ICP-OES in Security Applications in Windows 10**” for more information on doing this.

Create a System Restore Point

- 10) Click Search  and then type in “Create a restore point” in the search box




11) An entry entitled “Create a restore point” will appear. Click on this entry.

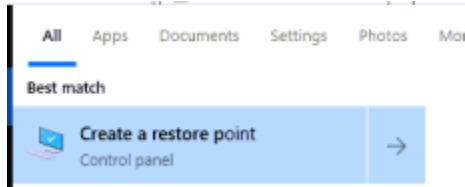


12) The System Properties popup window will appear. Click




Note: If System Protection is turned off, you will need to turn it on by clicking

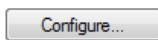

and then selecting the radio button: **Turn on system protection**

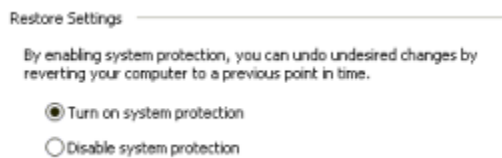

13) Name the restore point and click

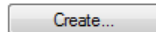

Choose a name that is relevant to the changes which will be made after the restore point has been created. For instance, since you are installing the Qtegra software, you might want to choose a name which references the Qtegra software version you will install.

Installing Qtegra on New Computer

Install Qtegra using the “*Qtegra for ICP-OES and ICP-MS Installation Guide*”, see PDF attachment in this document.

Installing Backed Up Files

Copy the files contained in the folders referenced in [Table 1](#) below from the backed up Qtegra location to the same folders on the new PC


Qtegra for ICP-OES Directory Reference Table	
File Type	File Path
Labbooks	
Labbook	C:\ProgramData\Thermo\Qtegra_Application Data\Workspace\LabBooks
Template	C:\ProgramData\Thermo\Qtegra_Application Data\Workspace\Templates

Reports	C:\ProgramData\Thermo\Qtegra_Application Data\Reporting
Export Schemes	C:\ProgramData\Thermo\Qtegra_Application Data\ExportSchemes

Table 1: Qtegra for ICP-OES Directory Reference Table

Note: If you are using an ESI FAST system and have created a custom FAST method, you will also want to copy custom methods from the backed up location: ***Qtegra_Application Data\ESI\Methods*** to ***C:\ProgramData\Thermo\Qtegra_Application Data\ESI\Methods*** on the new computer.

Create Configurations

Open  Configurator and create similar configurations to those which were recorded in the Capturing Qtegra Configurations section of this tech note.

Restore Plugin Settings

Input any changes to the default settings for the different plugins being used according to what was recorded in the **Capturing Qtegra Configurations** section of this document.