

Shutdown Procedure – iCAP 6000 & 7000 ICP OES

1. Make sure plasma is off and instrument is in Idle state.
2. Close Qtegra dashboard.
3. Turn off chiller with front control on/off push button. Then turn off chiller main power switch on rear above power cord.
4. Wait 15 minutes.
5. Turn off gas supplies.
6. Release tension on pump tubing.
7. Turn off main iCAP power switch on left side.



8. Turn off computer and accessories power; Auto-sampler, Fast Valve, Humidifier, etc.

Startup Procedure

1. Turn on iCAP main power switch.
2. Turn on gas supplies.
3. Turn on computer and accessories.
4. Wait 1 hour before turning on chiller to ignite plasma.