

Audit Manager Release Notes

Table of Contents

About This Document	3
Trademarks	3
Audit Manager v1.8	4
Release Date: Mar-2024	4
Supported Operating Systems	4
Supported Databases	4
New Features	5
Resolved Issues	5
Known Issues	5
Audit Manager v1.7 Hotfix 1	5
Release Date: April 2023	5
Supported Operating Systems	5
Supported Databases	6
New Features	6
Resolved Issues	6
Known Issues	6
Audit Manager v1.7	7
Release Date: December 2022	7
Supported Operating Systems	7
Supported Databases	7
New Features	7
Resolved Issues	7
Known Issues	8
Audit Manager v1.6 Hotfix 1	8
Release Date: August 2021	8
Supported Operating Systems	8
Supported Databases	8
New Features	9
Resolved Issues	9
Known Issues	9
Audit Manager v1.6	9
Release Date: April 2021	9
Supported Operating Systems	10
Supported Databases	10
New Features	10
Resolved Issues	10

Known Issues	10
Audit Manager v1.5	11
Release Date: December-2020	11
Supported Operating Systems	11
Supported Databases	11
New Features	11
Resolved Issues	12
Known Issues	12
Audit Manager v1.4 Hot Fix 1	12
Release Date: November-2020	12
Supported Operating Systems	12
Supported Databases	12
New Features	13
Resolved Issues	13
Known Issues	13
Audit Manager v1.4	13
Release Date: September-2020	13
Supported Operating Systems	13
Supported Databases	14
New Features	14
Resolved Issues	14
Known Issues	14
Audit Manager v1.3	15
Release Date: July-2020	15
Supported Operating Systems	15
Supported Databases	15
New Features	16
Resolved Issues	16
Known Issues	16
Audit Manager v1.2	16
Release Date: March-2020	16
Supported Operating Systems	16
Supported Databases	17
New Features	17
Resolved Issues	17
Known Issues	17
Audit Manager v1.1	18
Release Date: July-2019	18
Supported Operating Systems	18
Supported Databases	18
New Features	19
Resolved Issues	19
Known Issues	19

Audit Manager v1.0 Hot Fix 2	19
Release Date: 11-Apr-2018	19
Supported Operating Systems	19
Supported Databases	19
New Features	20
Resolved Issues	20
Known Issues	20
Audit Manager v1.0 Hot Fix 1	21
Release Date: 2-Mar-2018	21
Supported Operating Systems	21
New Features	21
Resolved Issues	21
Known Issues	21
Audit Manager v1.0	22
Release Date: 12-Jan-2018	22
Supported Operating Systems	22
Supported Databases	22
New Features	22
Resolved Issues	23
Known Issues	23
How to Contact Us	23

About This Document

This document contains a revision history of Audit Manager, including new features that may not be included in the User's Guide, resolved issues, and known issues.

Trademarks

Microsoft, Internet Explorer, and Windows are either trademarks or registered trademarks of Microsoft Corporation in the United States and/or other countries.

ToolBook is either a trademark or registered trademark of SumTotal in the United States and/or other countries.

Semprex is either a trademark or registered trademark of Semprex Corporation in the United States and/or other countries.

PerkinElmer is either a trademark or registered trademark of PerkinElmer, Inc. in the United States and/or other countries.

Sadtler and SearchMaster are trademarks of Bio-Rad Laboratories, Inc. in the United States and/or other

countries.

Bruker and OPUS are either trademarks or registered trademarks of Bruker Optics, Inc. in the United States and/or other countries.

All other trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries.

Audit Manager v1.8

Release Date: Mar-2024

Supported Operating Systems

- Server
 - Windows Server 2012 R2
 - Windows Server 2016 w/ Desktop Experience
 - Windows Server 2019 w/ Desktop Experience
- Workstation
 - Windows 10 64-bit Version 1607 and newer
 - Windows 11 64-bit

Supported Databases

- Maria DB
 - 10.3
 - 10.4
 - 10.5
- SQL Server
 - 2014
 - 2016
 - 2017
 - 2019
- Oracle
 - 19c

New Features

None

Resolved Issues

Only the First Failure to Log In is Recorded

When a user fails to log in, only the first failure is recorded. Subsequent failures are not recorded. All login attempts are now recorded in the audit log.

Known Issues

Selecting Refresh initiates Clear Filter

Workaround: Create a Preference with your favorite settings. Then when the filters are cleared, simply select your favorite preference to easily restore your favorites

Duplicate events can be created during import of ETX files

When events are imported from Windows Event Viewer (ETX files) multiple times, the import process can create duplicate events. No information is lost. Only occurs when user is upgraded from a release before 2017 and elects to import the events from older system.

Audit Manager warning message - report size is too large.

When creating an Audit Manager Report, a warning may appear stating that the report size is too large. The current design limits the user to 65535 bytes.

Audit Manager v1.7 Hotfix 1

Release Date: April 2023

Supported Operating Systems

- Server
 - Windows Server 2012 R2
 - Windows Server 2016 w/ Desktop Experience
 - Windows Server 2019 w/ Desktop Experience
- Workstation
 - Windows 10 64-bit Version 1607 and newer
 - Windows 11 64-bit

Supported Databases

- Maria DB
 - 10.3
 - 10.4
 - 10.5
- SQL Server
 - 2014
 - 2016
 - 2017
 - 2019
- Oracle
 - 19c

New Features

Added / updated events for OMNIC Paradigm

Added and updated event IDs used by OMNIC Paradigm.

Resolved Issues

Unable to retrieve user's display name

Under certain Active Directory configurations, Audit Manager is unable to properly retrieve a user's display name.

Known Issues

Selecting Refresh initiates Clear Filter

Workaround: Create a Preference with your favorite settings. Then when the filters are cleared, simply select your favorite preference to easily restore your favorites

Duplicate events can be created during import of ETX files

When events are imported from Windows Event Viewer (ETX files) multiple times, the import process can create duplicate events. No information is lost. Only occurs when user is upgraded from a release before 2017 and elects to import the events from older system.

Audit Manager warning message - report size is too large.

When creating an Audit Manager Report, a warning may appear stating that the report size is too large. The current design limits the user to 65535 bytes.

Audit Manager v1.7

Release Date: December 2022

Supported Operating Systems

- Server
 - Windows Server 2012 R2
 - Windows Server 2016 w/ Desktop Experience
 - Windows Server 2019 w/ Desktop Experience
- Workstation
 - Windows 10 64-bit Version 1607 and newer

Supported Databases

- Maria DB
 - 10.3
 - 10.4
 - 10.5
- SQL Server
 - 2014
 - 2016
 - 2017
 - 2019
- Oracle
 - 19c

New Features

Enabled creation of dump files upon application crash

Set up the system to automatically create dump files in the %ProgramData%\Thermo Scientific\Audit Manager\Dump Files folder if any of the applications crash.

Resolved Issues

Audit Manager warning message – report size is too large.

When creating an Audit Manager Report, a warning may appear stating that the report size is too large. The previous design limited the user to 65535 bytes.

Known Issues

Selecting Refresh initiates Clear Filter

Workaround: Create a Preference with your favorite settings. Then when the filters are cleared, simply select your favorite preference to easily restore your favorites

Duplicate events can be created during import of ETX files

When events are imported from Windows Event Viewer (ETX files) multiple times, the import process can create duplicate events. No information is lost. Only occurs when user is upgraded from a release before 2017 and elects to import the events from older system.

Audit Manager v1.6 Hotfix 1

Release Date: August 2021

Supported Operating Systems

- Server
 - Windows Server 2012 R2
 - Windows Server 2016 w/ Desktop Experience
 - Windows Server 2019 w/ Desktop Experience
- Workstation
 - Windows 10 64-bit Version 1607 and newer

Supported Databases

- Maria DB
 - 10.2
 - 10.3
 - 10.4
 - 10.5
- SQL Server
 - 2012
 - 2014
 - 2016
 - 2017
 - 2019
- Oracle

- 12c Release 1
- 12c Release 2
- 19c

New Features

None

Resolved Issues

Thermo Scientific Audit Log Service stops running

Certain combinations of software packages could cause an exception in Thermo Scientific Audit Log Service that would cause it to crash. Most commonly observed on systems with multiple client packages.

Fixed Audit Manager security Access control "Allow configuring database connection"

The Access Control "Allow configuring database connection" did not prevent the denied user from running the Configuration Utility (for database connections). This is now fixed.

Known Issues

Selecting Refresh initiates Clear Filter

Workaround: Create a Preference with your favorite settings. Then when the filters are cleared, simply select your favorite preference to easily restore your favorites

Duplicate events can be created during import of ETX files

When events are imported from Windows Event Viewer (ETX files) multiple times, the import process can create duplicate events. No information is lost. Only occurs when user is upgraded from a release before 2017 and elects to import the events from older system.

Audit Manager warning message - report size is too large.

When creating an Audit Manager Report, a warning may appear stating that the report size is too large. The current design limits the user to 65535 bytes.

Audit Manager v1.6

Release Date: April 2021

Supported Operating Systems

- Server
 - Windows Server 2012 R2
 - Windows Server 2016 w/ Desktop Experience
 - Windows Server 2019 w/ Desktop Experience
- Workstation
 - Windows 10 64-bit Version 1607 and newer

Supported Databases

- Maria DB
 - 10.2
 - 10.3
 - 10.4
- SQL Server
 - 2012
 - 2014
 - 2016
 - 2017
 - 2019
- Oracle
 - 12c Release 1
 - 12c Release 2
 - 18c

New Features

Added / updated events for OMNIC Paradigm

Added and updated event IDs used by OMNIC Paradigm.

Resolved Issues

None

Known Issues

None

Audit Manager v1.5

Release Date: December-2020

Supported Operating Systems

- Server
 - Windows Server 2012 R2
 - Windows Server 2016 w/ Desktop Experience
 - Windows Server 2019 w/ Desktop Experience
- Workstation
 - Windows 10 64-bit Version 1607 and newer

Supported Databases

- Maria DB
 - 10.0
 - 10.1
 - 10.2
 - 10.3
 - 10.4
- SQL Server
 - 2012
 - 2014
 - 2016
 - 2017
 - 2019
- Oracle
 - 12c Release 1
 - 12c Release 2
 - 18c

New Features

None

Resolved Issues

Security server name not read properly on OMNIC Paradigm systems

On Summit instruments and OMNIC Paradigm workstation installations, the security server name could fail to be picked up properly by Audit Manager resulting in an inability to review the audit log.

Known Issues

None

Audit Manager v1.4 Hot Fix 1

Release Date: November-2020

Supported Operating Systems

- Server
 - Windows Server 2012 R2
 - Windows Server 2016 w/ Desktop Experience
 - Windows Server 2019 w/ Desktop Experience
- Workstation
 - Windows 10 64-bit Version 1607 and newer

Supported Databases

- Maria DB
 - 10.0
 - 10.1
 - 10.2
 - 10.3
 - 10.4
- SQL Server
 - 2012
 - 2014
 - 2016
 - 2017
 - 2019

- Oracle
 - 12c Release 1
 - 12c Release 2
 - 18c

New Features

None

Resolved Issues

Upgrading from 1.3 to 1.4 failed to install all necessary files

Upgrading from Audit Log Service 1.3 to Audit Log Service 1.4 correctly attempted to downgrade several third-party DLLs; Windows Installer prohibited this move, resulting in the files not being present on the computer.

Known Issues

Security server name not read properly on OMNIC Paradigm systems

On Summit instruments and OMNIC Paradigm workstation installations, the security server name could fail to be picked up properly by Audit Manager resulting in an inability to review the audit log. Running ChangeServer and clicking OK is the workaround for this.

Audit Manager v1.4

Release Date: September-2020

Supported Operating Systems

- Server
 - Windows Server 2008 R2
 - Windows Server 2012
 - Windows Server 2012 R2
 - Windows Server 2016 w/ Desktop Experience
 - Windows Server 2019 w/ Desktop Experience
- Workstation
 - Windows 7 32-Bit
 - Windows 7 64-bit

- Windows 10 64-bit Version 1607 and newer

Supported Databases

- Maria DB
 - 10.0
 - 10.1
 - 10.2
 - 10.3
 - 10.4
- SQL Server
 - 2012
 - 2014
 - 2016
 - 2017
 - 2019
- Oracle
 - 11g Release 2
 - 12c Release 1
 - 12c Release 2
 - 18c

New Features

None

Resolved Issues

Deleting a measurement now logs an audit event and requires a reason

Previously, deleting a measurement in OMNIC Paradigm software did not require the user to enter a reason. Now, a reason is required, and the reason is logged as an event in Audit Manager.

Audit Manager no longer displays Unknown Account

If the event could not detect a Windows domain account, the account was displayed as unknown account. Now, the audit event captures the user currently signed into Windows.

Known Issues

Security server name not read properly on OMNIC Paradigm systems

On Summit instruments and OMNIC Paradigm workstation installations, the security server name could fail to

be picked up properly by Audit Manager resulting in an inability to review the audit log. Running ChangeServer and clicking OK is the workaround for this.

Audit Manager v1.3

Release Date: July-2020

Supported Operating Systems

- Server
 - Windows Server 2008 R2
 - Windows Server 2012
 - Windows Server 2012 R2
 - Windows Server 2016 w/ Desktop Experience
 - Windows Server 2019 w/ Desktop Experience
- Workstation
 - Windows 7 32-Bit
 - Windows 7 64-bit
 - Windows 10 64-bit Version 1607 and newer

Supported Databases

- Maria DB
 - 10.0
 - 10.1
 - 10.2
 - 10.3
 - 10.4
- SQL Server
 - 2012
 - 2014
 - 2016
 - 2017
 - 2019
- Oracle

- 11g Release 2
- 12c Release 1
- 12c Release 2
- 18c

New Features

None

Resolved Issues

Incorrect characters in some event descriptions

When using mixed languages between the application and the server, multi-byte characters may not appear correctly. The malformed characters will not be recoverable.

Known Issues

Translations are not complete

All user interface text and some event descriptions will be in English only.

Security server name not read properly on OMNIC Paradigm systems

On Summit instruments and OMNIC Paradigm workstation installations, the security server name could fail to be picked up properly by Audit Manager resulting in an inability to review the audit log. Running ChangeServer and clicking OK is the workaround for this.

Audit Manager v1.2

Release Date: March-2020

Supported Operating Systems

- Server
 - Windows Server 2008 R2
 - Windows Server 2012
 - Windows Server 2012 R2
 - Windows Server 2016 w/ Desktop Experience
 - Windows Server 2019 w/ Desktop Experience
- Workstation

- Windows 7 32-Bit
- Windows 7 64-bit
- Windows 10 64-bit Version 1607 and newer

Supported Databases

- Maria DB
 - 10.0
 - 10.1
 - 10.2
- SQL Server
 - 2012
 - 2014
 - 2016
- Oracle
 - 11g Release 2
 - 12c Release 1
 - 12c Release 2

New Features

Added Support for New Features in OMNIC Paradigm Software

Database Maintenance – Backup and Restore

Options – Changes on display for line weight, colors, and digits

Report Options

Resolved Issues

None

Known Issues

Translations are not complete

All user interface text and some event descriptions will be in English only.

Incorrect characters in some event descriptions

When using mixed languages between the application and the server, multi-byte characters may not appear correctly. The malformed characters will not be recoverable.

Security server name not read properly on OMNIC Paradigm systems

On Summit instruments and OMNIC Paradigm workstation installations, the security server name could fail to be picked up properly by Audit Manager resulting in an inability to review the audit log. Running ChangeServer and clicking OK is the workaround for this.

Audit Manager v1.1

Release Date: July-2019

Supported Operating Systems

- Server
 - Windows Server 2008 R2
 - Windows Server 2012
 - Windows Server 2012 R2
 - Windows Server 2016 w/ Desktop Experience
- Workstation
 - Windows 7 32-Bit
 - Windows 7 64-bit
 - Windows 10 64-bit Version 1607 and newer

Supported Databases

- Maria DB
 - 10.0
 - 10.1
 - 10.2
- SQL Server
 - 2012
 - 2014
 - 2016
- Oracle
 - 11g Release 2
 - 12c Release 1
 - 12c Release 2

New Features

Added support for OMNIC Paradigm

Support for the OMNIC Paradigm software

Resolved Issues

None

Known Issues

Translations are not complete

All user interface text and some event descriptions will be in English only.

Incorrect characters in some event descriptions

When using mixed languages between the application and the server, multi-byte characters may not appear correctly. The malformed characters will not be recoverable.

Audit Manager v1.0 Hot Fix 2

Release Date: 11-Apr-2018

Supported Operating Systems

- Server
 - Windows Server 2008 R2
 - Windows Server 2012
 - Windows Server 2012 R2
 - Windows Server 2016 w/ Desktop Experience
- Workstation
 - Windows 7 32-Bit
 - Windows 7 64-bit
 - Windows 8.1 64-bit
 - Windows 10 64-bit

Supported Databases

- Maria DB

- 10.0
- 10.1
- 10.2
- SQL Server
 - 2012
 - 2014
 - 2016
- Oracle
 - 11g Release 2
 - 12c Release 1
 - 12c Release 2

New Features

None

Resolved Issues

Certain network configurations prohibited retrieving the user's full name during service startup.

This was found to be a timing issue with the service startup when on a domain computer but using a local service account.

Certain network configurations prohibited retrieving the group membership for an account during database configuration.

The group membership check is no longer necessary and has been removed to avoid requiring additional permissions during database configuration.

Known Issues

Translations are not complete

All user interface text and some event descriptions will be in English only.

Incorrect characters in some event descriptions

When using mixed languages between the application and the server, multi-byte characters may not appear correctly. The malformed characters will not be recoverable.

Audit Manager v1.0 Hot Fix 1

Release Date: 2-Mar-2018

Supported Operating Systems

- Server
 - Windows Server 2008 R2
 - Windows Server 2012
 - Windows Server 2012 R2
 - Windows Server 2016 w/ Desktop Experience
- Workstation
 - Windows 7 32-Bit
 - Windows 7 64-bit
 - Windows 8.1 64-bit
 - Windows 10 64-bit

New Features

None

Resolved Issues

Certain Active Directory security configurations prohibited authenticating the service account during database configuration.

This was found to be an issue with the authentication options specified when communicating with the domain controller.

NTFS folder permissions could prevent secure data storage from saving data

The default folder permissions for user folders on newer Windows installations could cause the Secure Data Storage service to be unable to place data files in the secure location.

Known Issues

Translations are not complete

All user interface text and some event descriptions will be in English only.

Incorrect characters in some event descriptions

When using mixed languages between the application and the server, multi-byte characters may not appear correctly. The malformed characters will not be recoverable.

Audit Manager v1.0

Release Date: 12-Jan-2018

Supported Operating Systems

- Server
 - Windows Server 2008 R2
 - Windows Server 2012
 - Windows Server 2012 R2
 - Windows Server 2016 w/ Desktop Experience
- Workstation
 - Windows 7 32-Bit
 - Windows 7 64-bit
 - Windows 8.1 64-bit
 - Windows 10 64-bit

Supported Databases

- Maria DB
 - 10.0
 - 10.1
 - 10.2
- SQL Server
 - 2012
 - 2014
 - 2016
- Oracle
 - 11g Release 2
 - 12c Release 1
 - 12c Release 2

New Features

Initial Release, all features are new

Resolved Issues

Initial Release, all features are new

Known Issues

Translations are not complete

All user interface text and some event descriptions will be in English only.

Incorrect characters in some event descriptions

When using mixed languages between the application and the server, multi-byte characters may not appear correctly. The malformed characters will not be recoverable.

How to Contact Us

Current contact information is located at <https://www.thermofisher.com>

Select the "Contact Us" icon at the top of the screen