


KNOWLEDGE SHARING

Question/Issue/Symptom/Problem:

How to Replace Legacy ULT Upright Freezer Battery

Resolution/Answer:

Follow Procedure Below:

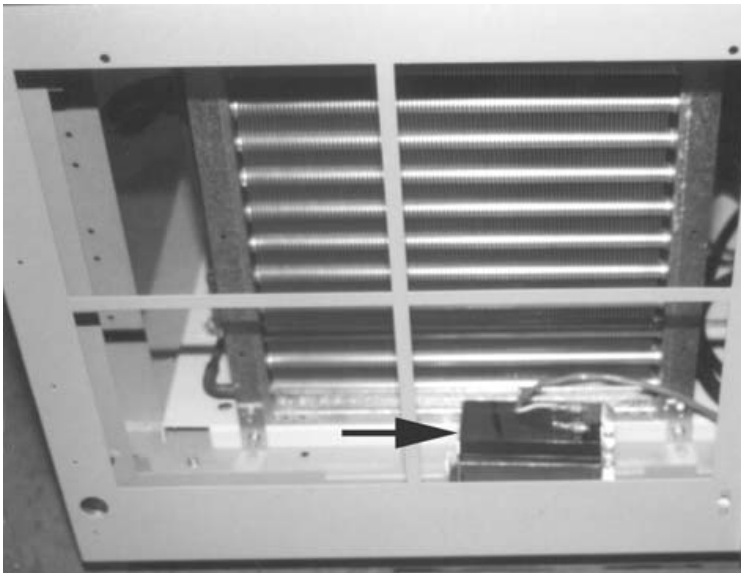

1. Turn the unit Off, unplug the freezer from the wall receptacle.
2. Locate the alarm battery, mounted on the floor behind the condenser/air filter grill. Please see example photo above after front bottom grill panel is removed.
3. If battery wires are not labeled, then label the positive and negative wire and disconnect the two leads from the gel cell battery.
4. Remove the screws and metal retaining bracket; remove battery.
5. Install new battery in the reverse process. Allow about a week to ten days for battery to fully charge.

Title:

Legacy ULT Upright Freezer Battery Replacement Instructions

Audience:

Public

Author:

Ron Prickett
