

# Service

# Spare Parts



Material	Equipment-number
75002430	41547414
75002431	41540951
75002432	41540958
75002435	41545641
75002436	41546059
75002449	41553372
75002450	41540959
75002461	41540964

Material	Equipment-number
75002410	41540923
75002411	41540940
75002415	41540942
75002416	41540950
75002465	41545656
75002491	41540981
75002494	41540982
75002495	41540979

The original version of this manual is valid. Thermo Scientific cannot guarantee the accuracy or completeness of subsequent copies.

## 6.0 SPARE PARTS LIST





# Microliter ventilated spareparts

**Heraeus:** Pico 17/21

**Sorvall:** Legend Micro 17/21

**Thermo Scientific:** MicroCL 17/21

**Fisher:** AccuSpin Micro 17


# Centrifuges Overview – Zentrifugen Übersicht

---

**Heraeus:** Pico 17/21  
**Sorvall:** Legend Micro 17/21  
**Thermo Scientific:** Micro CL 17/21  
**Fisher:** AccuSpin Micro 17



- casing and electronics –  
Gehäuse und Elektronik

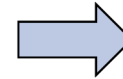



# Spareparts Microliter centrifuge

Pos.	Part Nr.:	Description:	Beschreibung:	Heraeus/ Thermo 120V	Heraeus/ Thermo 230V	Sorvall/ Fisher 120V	Sorvall/ Fisher 230V
100	20058271	Housing Sorvall	Gehäuse Sorvall			X	X
100	20058095	Housing Heraeus	Gehäuse Heraeus	X	X		
102	20057553	Holder tool	Werkzeughalter	X	X	X	X
103	20360104	Screw driver 5mm black	Sechskant 5mm	X	X	X	X
104	20058230	Gasket	Deckeldichtung	X	X	X	X
105	70060419	Lid chassis	Deckelblech	X	X	X	X
106	20058296	Lid cover Sorvall	Deckeloberschale Sorvall			X	X
106	20058096	Lid cover Heraeus	Deckeloberschale Heraeus	X	X		
107	20050543	Window	Sichtfenster	X	X	X	X
108	20058092	Lid axis	Deckelachse	X	X	X	X
109	20350211	Collar	Gleitbuchse	X	X	X	X
110	20059010	Nut	Deckelmutter	X	X	X	X
111	70902024	Foam silencer	Schaumblock	X	X	X	X
113	20058028	Grill 3x3mm	Gitternetz fein 3x3mm	X	X	X	X
114	20058319	Foam rear	Dämmeinsatz hinten	X	X	X	X
115	20058093	Spring	Deckelfeder	X	X	X	X
125	20058715	Key foil „Pico 17“	Tasterfolie „Pico 17“	X	X		
125	20058716	Key foil „Pico 21“	Tasterfolie „Pico 21“	X	X		
125	20058719	Key foil „Legend Micro 17“	Tasterfolie „Legend Micro 17“			X	X
125	20058720	Key foil „Legend Micro 21“	Tasterfolie „Legend Micro 21“			X	X
125	20058723	Key foil „Micro CL 17“	Tasterfolie „Micro CL 17“	X	X		
125	20058724	Key foil „Micro CL 21“	Tasterfolie „Micro CL 21“	X	X		
125	20058727	Key foil „Pico 17“	Tasterfolie „Pico 17“			X	X

# Spareparts Microliter centrifuge

Pos.	Part Nr.:	Description:	Beschreibung:	Heraeus/ Thermo 120V	Heraeus/ Thermo 230V	Sorvall/ Fisher 120V	Sorvall/ Fisher 230V
130	20150328	PCB display	Anzeigeplatte	X	X	X	X
131	50104102	Flat ribbon cable	Flachbandkabel	X	X	X	X
135	20901161	Latch	Deckelschloss	X	X	X	X
140	70902030	Motor 3PH	Motor 3PH	X	X	X	X
141	20150317	PCB tachboard	Drehzahlerfassung Platine	X	X	X	X
150	20150396	PCB main assy	Hauptplatine		X		X
150	20150397	PCB main assy	Hauptplatine	X		X	
151	<b>see next page</b>						
152	50141342	Splash guard	Spritzschutz	X	X	X	X
153	20160751	Resistor	Vorwiderstand		X		X
154	20230218	Fuse 4 AMP slow blown	Sicherung 4A Träge		X		X
154	20230228	Fuse 6,3 AMP slow blown	Sicherung 6,3A Träge	X		X	
155	70902023	Service kit AV mounts	Service kit Gummipuffer	X	X	X	X
160	70060482	Bottom plate Sorvall	Bodenplatte Sorvall			X	X
160	70060418	Bottom plate Heraeus	Bodenplatte Heraeus	X	X		
161	20058120	Foot	Stellfüsse	X	X	X	X
162	20058029	Foam air channel	Schaumstoff Luftführung	X	X	X	X
165	20180628	Socket	Geräteanschluss	X	X	X	X
180	20190246	Mains cable for power supply 230V	Geräteanschlussleitung 230V		X		X
180	20190247	Mains cable for power supply 120V	Geräteanschlussleitung 120V	X		X	

