

Instruction Sheet		
GP3 Centrifuges	Clearing Password	February 3, 2015

To reset the pass code assigned to one of the program memories, proceed as follows:

- Switch off the centrifuge
- Press program key 1 and 2 and switch on the centrifuge
- When asked "Enter service menu?", press START
- Press ENTER three times until "Clr prog passcode?" appears in the display
- Press 1 on numerical keypad to change a NO into a YES
- Press ENTER
- "Master passcode" appears on the display
- Enter the master passcode of this centrifuge (0000 if not changed) press ENTER
- "Clear code program n°" appears on the screen
- Enter program number of which to clear the passcode
- Press ENTER
- "Program passcode cleared" appears on the display for three seconds
- "Clear code program n°" appears on the screen
- Continue with other program number or press STOP two times to leave service menu