



To add the macro to the toolbar perform the following:


- In OMNIC - Go to Edit: Edit Toolbar.
- Select Add Item in the upper right.

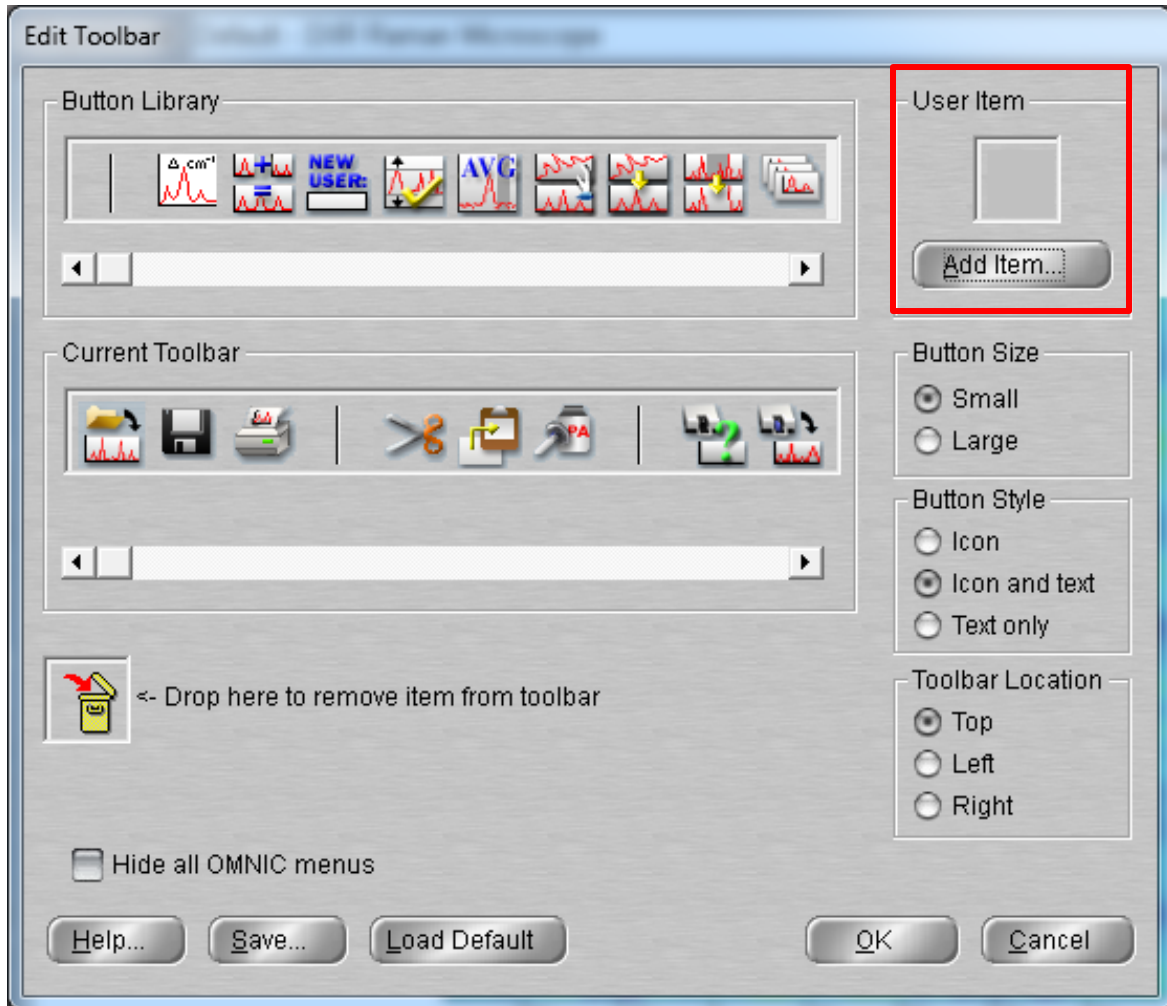



- Select Run Basic Macro and then click Browse.

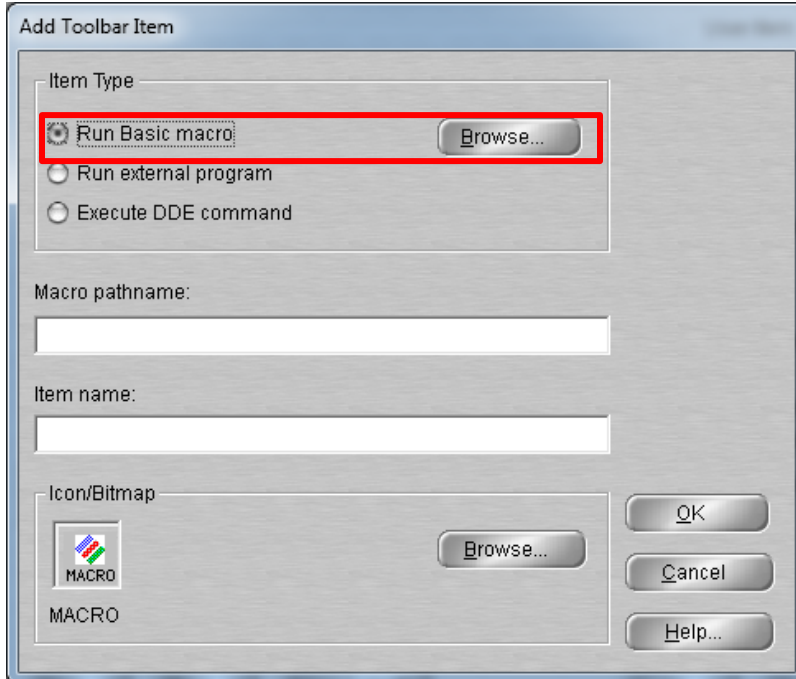

- Browse to the macro and select it. Default location is C:\My Documents\Omic\Macro\
- Modify Item name and change the icon if desired and Click OK.

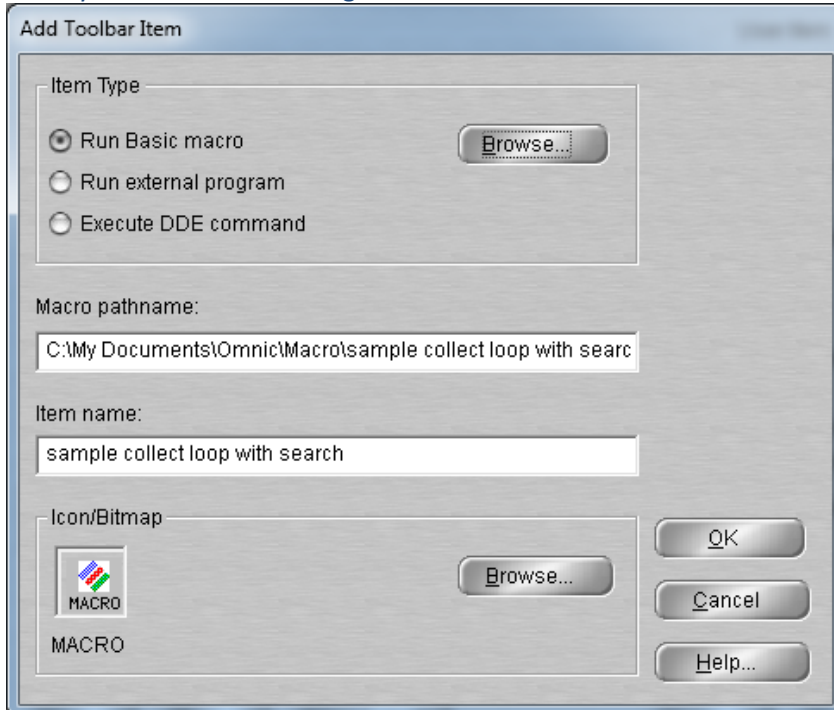

- Drag and drop the Macro icon from the User Item to the Current Toolbar location you want it displayed.

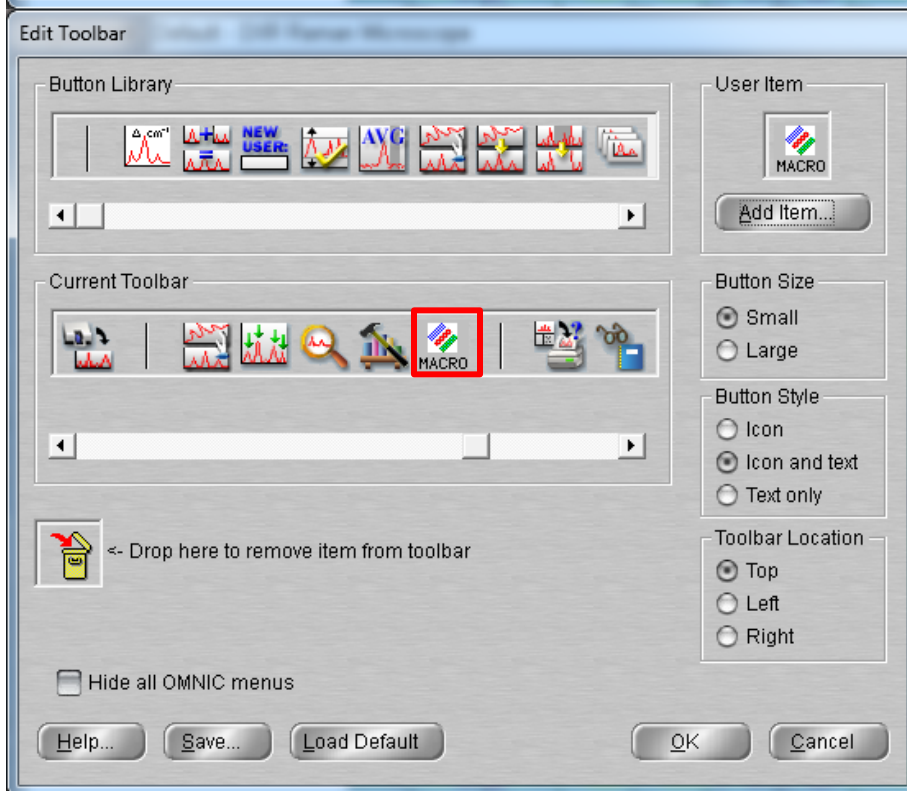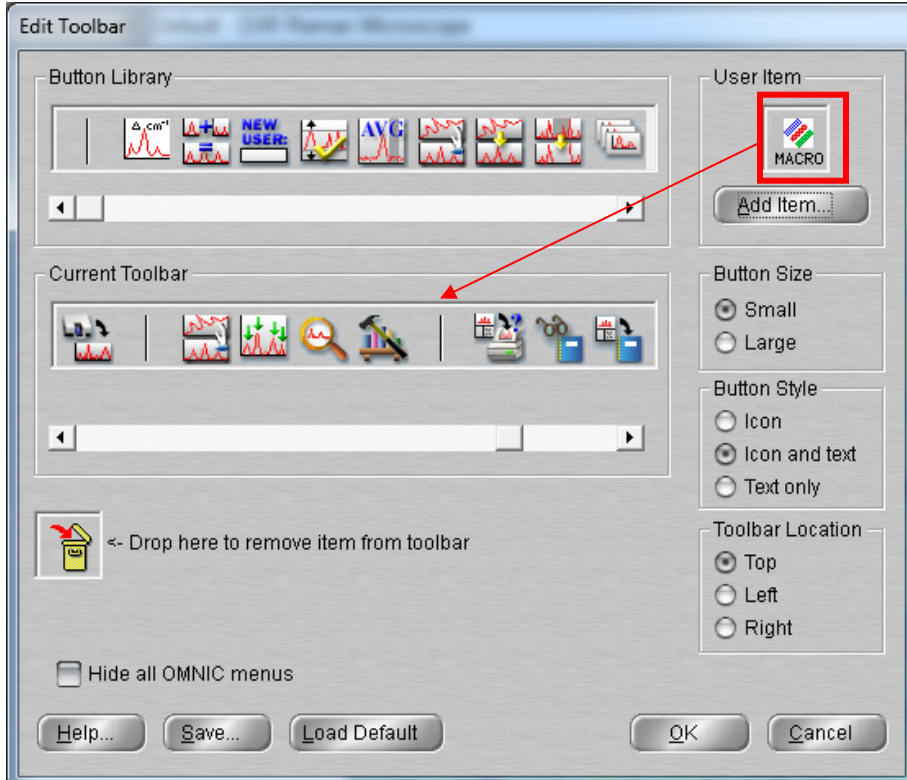

- Click OK
- Save the configuration using “Save Configuration As” from the file menu. If you do not save the configuration, the icon will not be there the next time OMNIC is opened.
- Click on the icon to run the macro when needed.