IsoTemp™ General Purpose Laboratory Refrigerators and Freezers

Installation and Operation

331950H03 • Revision B • 7/22/21

IMPORTANT Read this instruction manual. Failure to follow the instructions in this manual can result in damage to the unit, injury to operating personnel and poor equipment performance.

CAUTION All internal adjustments and maintenance must be performed by qualified service personnel.

Material in this manual is for informational purposes only. The contents and the product it describes are subject to change without notice. The Fisher Scientific makes no representations or warranties with respect to this manual. In no event shall Fisher Scientific be held liable for any damages, direct or incidental, arising from or related to the use of this manual.

© 2021 Fisher Scientific Inc. All rights reserved.

Contents

Models	1
Safety Information	2
Alert Signals	2
Intended Use	3
General Purpose Refrigerator & Freezer	4
Product Specification	4
Unpacking and Inspection	5
Unpacking	5
Visible Loss or Damage	5
Concealed Loss or Damage	5
Operating Conditions	6
Installation	7
Selecting a Location	7
Leveling the Unit	7
Electrical Connection	7
Installing the Door Handle for FBV18CPSA Model ...	7
Be Advised	9
Operation	10
Start-Up Procedure	11
Control panel FBV18CPSA	11
Control panel FBV20RPSA and FBV20FPSA	11
Restart Procedure	12
How to Save Energy	12

Safety Tips	12
Maintenance	13
Cleaning the Unit	13
Cleaning Interior/Exterior/Door Gaskets	13
Cleaning the Condenser	13
Transportation and Storage	13
Troubleshooting	14
Wiring Diagram	16
End of Life Care	18
Warranty	19
Regulatory Compliance	20
Product Safety	20
Electromagnetic Compatibility	20
Contact Information	22

Models

The table below shows the units covered in this operation and installation manual by model number.

Table 1. Applicable Models

Refrigerator - Freezer	Convertible Refrigerator / Freezer
FBV18CPSA	FBV20RPSA / FBV20FPSA

Safety Information

It is your responsibility to understand the proper function and operational characteristics of the unit. This instruction manual should be thoroughly read and all operators must be trained before attempting to place these units in service. Awareness of the stated cautions and warnings, and compliance with recommended operating parameters together with maintenance requirements are important for safe and satisfactory operation. The units should be used for its intended application, alterations or modifications will void the Warranty.

These products are not intended, nor can they be used, as a sterile or patient connected device. In addition, these units are not designed for use in Class I, II or III locations as defined by the US National Electrical Code, unless otherwise noted.

Alert Signals

Safety Alert: Important operating instructions. To reduce the risk of injury or poor performance of the unit. Read the user manual before putting the equipment into operation.

WARNING: Indicates an immediately hazardous situation, which if not avoided, will result in death or serious injury.

CAUTION: Indicates an immediately hazardous situation, which if not avoided, may result in minor to moderate injury.

WARNING: Avoid contact with cold freezer surfaces potential for cold burns or skin sticking to cold surfaces.

WARNING: Use of this symbol involves risk of electric shock.

Do not modify or use power supplies other than OEM equipment. Connection of the power supply may require a properly grounded receptacle. Potential for electrical shock or equipment damage exists if precautions are not followed.

Note: Notes alert you to pertinent facts and conditions.

DANGER RISK OF CHILD ENTRAPMENT
Before you throw away your old refrigerator or freezer:

- Take off doors.
- Leave the shelves in the place so that children may not easily climb inside.

WARNING: If the equipment is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired.

WARNING: Units are charged with hydrocarbon refrigerant (R600a). Only qualified service personnel should service this unit.

WARNING: Risk of fire or explosion. Flammable refrigerant used. To be repaired only by trained service personnel. Do not puncture refrigerant tubing.

WARNING: No equipment that uses an open flame should be placed inside the refrigerator. This will harm the unit, hamper functionality and compromise your safety.

WARNING: Risk of fire or explosion. Flammable refrigerant used. Do not use mechanical devices to defrost the unit.

WARNING: This product can expose you to chemicals including lead, which is known to the State of California to cause cancer, and bisphenol A (BPA) and lead which are known to the State of California to cause birth defects or other reproductive harm. For more information go to www.P65Warnings.ca.gov.

Intended Use

The General Purpose refrigerator and freezer described in this manual is for professional use only.

These products are intended for use in research for the storage of samples or inventory in the following temperature ranges:

Refrigerator +33.8°F to +53.6°F (+1°C to +12°C)

Freezer -11.2°F to +10.4°F (-24°C to -12°C)

These units are not considered as medical devices and have therefore not been registered with a medical device regulatory body (e.g. FDA): that is, these have not been evaluated for the storage of samples for diagnostic use or for samples to be re-introduced to the body.

Note: Do not store corrosive materials in these units. Any damage which occurs due to storage of corrosives will not be covered under warranty claims.

Follow site protocols to identify materials of corrosive nature.

General Purpose Refrigerator & Freezer

These are general purpose units available in all refrigerator and freezer:

Model: FBV18CPSA - 18 Cu.ft

30" Wide, Refrigerator and Freezer

Features:

- Color: White
- Defrost Type: Auto
- Interior Light: Yes
- Control Type: Electronic
- Defrost Drain: Yes
- Refrigerator: 2 Adjustable shelves, 2 Bulk basket storage
- Freezer: 1 Adjustable shelf
- Leveling Legs: 2

Model: FBV20RPSA - 21 Cu.ft

33" Wide, Convertible Refrigerator / Freezer

Features:

- Color: White
- Defrost Type: Auto
- Interior Light: Yes
- Control Type: Electronic
- Defrost Drain: Yes
- Cabinet: 4 Adjustable shelves, 1 Bulk basket storage
- Leveling Legs: 2
- Convertible: Yes
- Default setting: Refrigerator

Model: FBV20FPSA - 21 Cu.ft

33" Wide, Convertible Refrigerator / Freezer

Features:

- Color: White
- Defrost Type: Auto
- Interior Light: Yes
- Control Type: Electronic
- Defrost Drain: Yes
- Cabinet: 4 Adjustable shelves, 1 Bulk basket storage
- Leveling Legs: 2
- Convertible: Yes
- Default setting: Freezer

Product Specification

	FBV18CPSA	FBV20RPSA / FBV20FPSA
Exterior Dimensions H x W x D inches (cm)	66.61" x 29.53" x 32.79" 169 x 75 x 83.28 cm	76.77" x 32.76" x 28.94" 194.9 x 83.21 x 73.5 cm
Net Weight lbs. (kg)	175 lbs. (79 kg)	196 lbs. (89 kg)
Shipping Weight lbs. (kg)	224 lbs. (101.6 kg)	244 lbs. (110.6 kg)
Refrigerator Temperature Range °F (°C)	33.8°F to 53.6°F +1°C to +12°C	33.8°F to 53.6°F +1°C to +12°C
Freezer Temperature Range °F (°C)	-11.2°F to 10.4°F -24°C to -12°C	-11.2°F to 10.4°F -24°C to -12°C
Electrical Characteristics Volts/Hz, Amps	115 V / 60 Hz, 1.1 A	115 V / 60 Hz, 2.2 A
Maximum allowed load on the shelves lbs (kg)	55 lbs. (25 kg)	55 lbs. (25 kg)

Note: Amps listed are at normal run mode, starting amps may be higher.

Unpacking and Inspection

CAUTION: DO NOT REMOVE, under any circumstance, the grounding prongs from the 3-prong power cord supplied with all units.

CAUTION: DO NOT USE electrical extension cords that may result in voltage loss and possible hazardous operation.

Unpacking

This should be inspected upon delivery. When received, carefully examine for any shipping damage before unpacking. If damage is discovered, the delivering carrier should both specify and sign for the damage on your copy of the delivery receipt.

Visible Loss or Damage

Note any external evidence of loss or damage on the freight bill, or express receipt, and have it signed by the carrier's agent. Failure to adequately describe such external evidence of loss or damage may result in the carrier's refusing to honor your damage claim. The form required to file such a claim will be supplied by the carrier.

IMPORTANT: Failure to request an inspection of damage within a few days after receipt of shipment absolves the carrier from any liability for damage. You must call for a damage inspection promptly.

Concealed Loss or Damage

Concealed loss or damage refers to loss or damage, which does not become apparent until the merchandise has been unpacked and inspected. If either occur, make a written request for the carrier's agent within 15 days of the delivery date; then file a claim with the carrier since the damage is the carrier's responsibility.

If you follow the above instructions carefully, we will guarantee our support of your claim to be compensated for loss from concealed damage.

DO NOT, FOR ANY REASON RETURN THIS UNIT WITHOUT FIRST OBTAINING AUTHORIZATION.

Location of parts

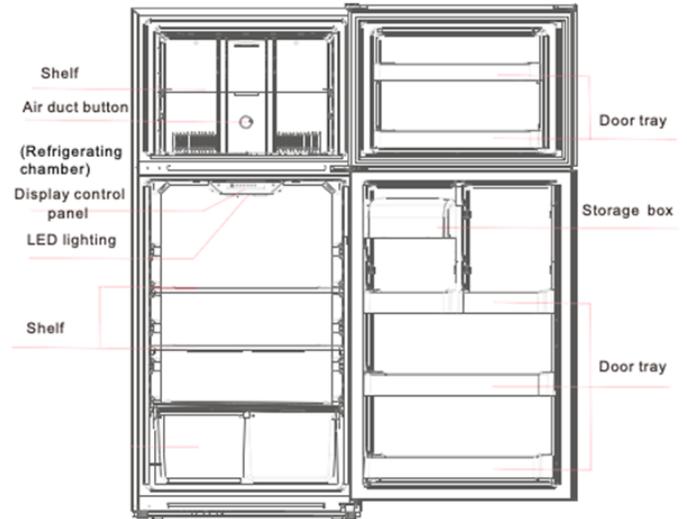
Figure 1. FBV18CPSA Model

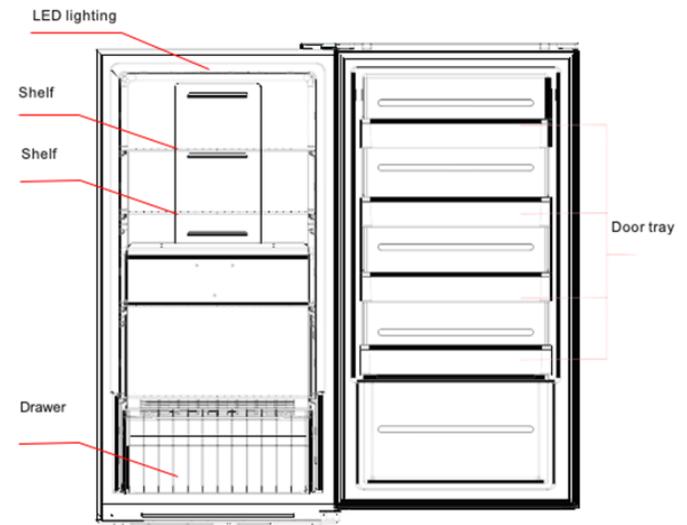
Figure 2. FBV20RPSA and FBV20FPSA Model

Note: Figure 1 & 2 are only for reference. The actual configuration will depend on the physical product or statement by the distributor.

Operating Conditions

The units described in this manual are classified for use as stationary equipment in a Pollution Degree 2 and Over voltage Category II environment.

These units are designed to operate under the following environmental conditions:

- Indoor use
- Altitude up to 2,000 m (6512 feet)
- Maximum relative humidity 65%, non-condensing
- Temperatures: 59°F to 90°F (15°C to 32°C)
- Main supply voltage fluctuations should not exceed by $\pm 10\%$ of the nominal voltage.

Installation

Selecting a Location

Choose a location for the FBV20RPSA, FBV20FPSA and FBV18CPSA that provides a clearance of 11" (30 cm) at the top, 4" (11 cm) at the rear and 4" (11 cm) on each side.

Appropriate electrical power must be available. Attach the refrigerator and freezer to the facility's electrical supply as directed by the National Electrical Code.

Installation Instructions

Remove all packing materials, including the bottom cushions, foam pads, and tape inside the unit. Tear off all protective film on the doors and on the refrigerator and freezer body.

- Place the unit in a well ventilated place, on flat sturdy ground.
- Avoid locating the unit in moist or humid areas. Excess moisture in the air may cause frost to form quickly on the evaporator and cabinet, requiring frequent defrosting.
- Locate the unit away from direct sunlight and other heat sources. Direct sunlight may affect the performance of the unit.

Let the freezer stand upright for approximately two hours before connecting it to power. This reduces the possibility of a malfunction in the cooling system from improper handling during transportation.

While moving the unit, DO NOT tilt it beyond 45° from upright.

Leveling the Unit

It is very important for the refrigerator and freezer to be level. If the refrigerator and freezer is not leveled during installation, the doors may not close or seal properly, causing cooling, frost or moisture problems.

To level your refrigerator and freezer, rotate the leveling feet clockwise (to raise) or counter-clockwise (to lower).

Note: Have someone push against the top of the refrigerator and freezer to help take some weight off the leveling foot, making it easier to adjust.

Electrical Connection

Determine the total amount of current presently being used by other apparatus connected to the circuit that will be used by this refrigerator and freezer. It is critical that this added current demand and other equipment on this circuit not exceed the rating of the fuse or circuit breaker in use.

The frequency and nominal voltage requirements for the unit are specified on the data plate. Supply this unit with an electrical source that meets these requirements only. Low line voltage is often the cause of service complaints. With the unit running, check that the line voltage is within $\pm 10\%$ of that specified on the data plate.

CAUTION: Be sure the voltage supplied to the refrigerator and freezer is equal to that specified on the data plate.

WARNING: For personal safety, this unit must be properly grounded before use.

General Purpose units must be connected to a grounded outlet matching the nameplate and / or the information furnished in this manual. If you are not sure about the outlet, you should contact a qualified electrician for assistance.

The Unit should always be connected to a dedicated power source.

WARNING: DO NOT REMOVE, under any circumstance, the grounding prongs from the 3-prong power cord supplied with all units.

WARNING: DO NOT USE electrical extension cords, it may result in voltage loss and possible hazardous operation.

Installing the Door Handle for FBV18CPSA Model

Note: The figures shown in this user manual are only for reference. The actual product may differ slightly. The screws and wrench are in the plastic bag.

A Phillips head screwdriver is the only equipment needed to install the door handle. Follow these steps for a proper installation.

1. Use your Phillips head screwdriver to fix the bolts (1) into the bolt holes (2) on the refrigerator door.
2. Put the handle on the door through the bolt (1), making sure the handle contacts the surface of the door.
3. Use the handle screws (3) to fix the handle to the door. Make sure the handle is attached firmly.
4. Put the cap (4) on the handle screw hole.
5. Follow the same procedure for installing the door handle on the freezer door.

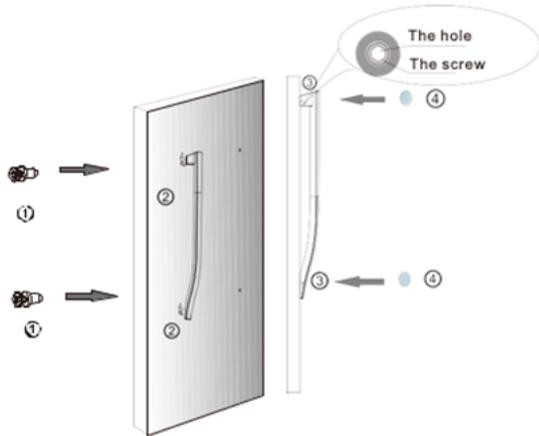
Figure 3. FBV18CPSA door handle installation


Reversing the Direction of the Door Swing

This unit has the ability to open from either the right or the left side. Reversing the door swing can be done yourself using the following tools:

- Phillips head screwdriver
- Putty knife or flat head screwdriver
- 5/16" socket and ratchet
- Masking tape

Follow these steps to reverse the direction of the door swing:


1. Unplug the refrigerator, and remove all the contents. Close the door and tape the door to the cabinet.
2. Remove the left cover plate and the top right hinge cover. Unscrew the screws, and remove the top right hinge. Save the screws, as you will need them to reattach the door.

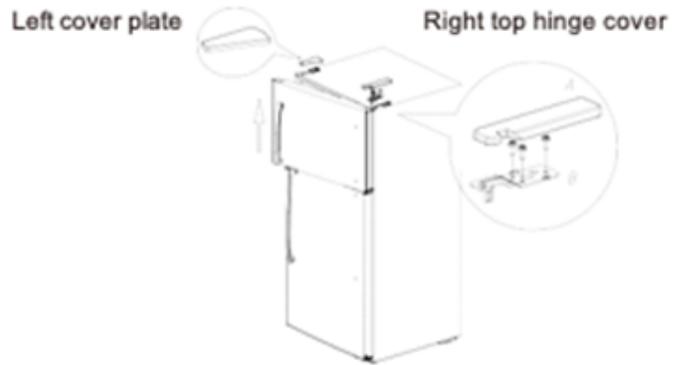3. Lift the freezer door up and away from the cabinet.
4. Unscrew the screws of the middle hinge, remove the hinge, and then lift the refrigerator door up and away from the cabinet. Save the screws, as you will need them to reattach the door.


5. Unscrew the screws of the bottom hinge, then remove the bottom hinge from the bottom right side of the refrigerator.
6. Unscrew the hinge shaft from right side of the bottom hinge, and move it to the left side.
7. Secure the bottom hinge to the left side of the refrigerator, using the screws you previously removed.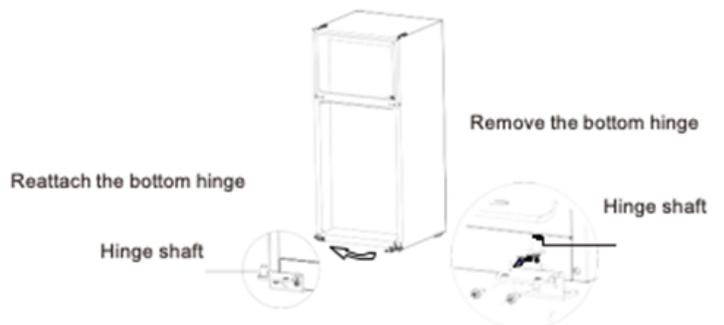


8. Dismantle the sleeves from the top right of the doors and the hole cap from the top left side of the doors. Assemble them on the opposite side.

9. Remove the refrigerator door stopper, and install it on the other side of the door.
10. Remove the freezer door self-locking block and stopper. Save them to reuse later.
11. Take out the self-locking block and stopper from the accessory bag in the refrigerator and install on the left side of the door.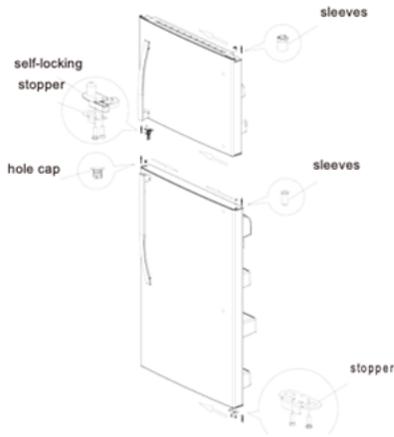

12. Set the refrigerator door down onto the bottom hinge (vertically), then take out the middle hinge for the other side from the accessory bag, and assemble the middle hinge.

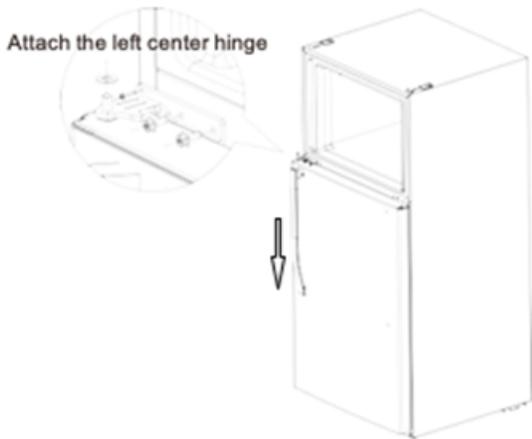13. Set the freezer door down onto the middle hinge (vertically), then take out the top hinge and cover, attach the top hinge, hinge cover, and hole cover.

Be Advised

WARNING: Storage by user of any materials in the product that may cause a deterioration of the product shall be deemed to constitute abnormal and improper usage of the product for purposes of this warranty.

WARNING: Before connecting the final power supply, check the electrical characteristics of the unit nameplate to see that it is in agreement with the power supplied. In addition, power should be wired to the unit according to the electrical schematic and all applicable codes. Only qualified electricians should work on the electrical portion of any unit installation.

Operation

CAUTION: Do not use in the presence of flammable or combustible materials or explosive gases. Fire or explosion could result, causing death or severe injury.

Follow site protocols for loading and unloading samples. Overloading the unit forces the compressor to run longer.

Wipe containers dry before placing items in the unit. This cuts down on frost build-up inside the unit.

The unit storage shelves should not be lined with aluminum foil, wax paper or paper toweling. Liners interfere with cold air circulation, making the unit less efficient.

Organize and label to reduce door openings and extended searches. Remove as many items as needed at one time, and close the door as soon as possible.


Before Using the Unit

1. Keep the unit upright for half an hour before connecting to the power supply.
2. Connect the unit to a power supply, and let the unit run for 2-3 hours.

Note: If the ambient temperature of the room is higher than normal (starting the machine in the summertime, for example), it may take up to 4 hours to start using the unit.

Start-Up Procedure

Control panel FBV18CPSA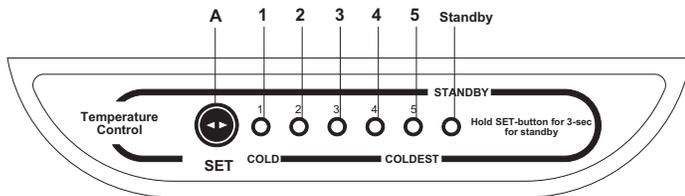

Once the unit is turned on, the display screen will display for three seconds. The unit will run on Setting 3 automatically.

Temperature Setting

The temperature settings run from cold (1) to coldest (5), each time the Temperature Setting button (A) is pressed. After the appropriate setting is selected, the unit will run according to that setting.


Note: The unit will start running on the set value 25 seconds after choosing the particular setting.

Standby Mode


- Press the Temperature Setting (A) button for three seconds to activate standby mode. The Standby indicator light will illuminate.
- Press the Temperature Settings (A) button for three seconds to resume normal refrigerator operation. The Standby indicator light will go out.

Setting the Freezer Temperature

- Turn the freezer temperature dial clockwise to reach a colder temperature.
- Turn the freezer temperature dial counterclockwise to reach a less cold temperature.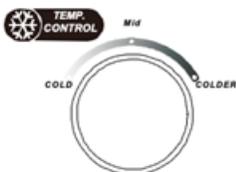

Control panel FBV20RPSA and FBV20FPSA

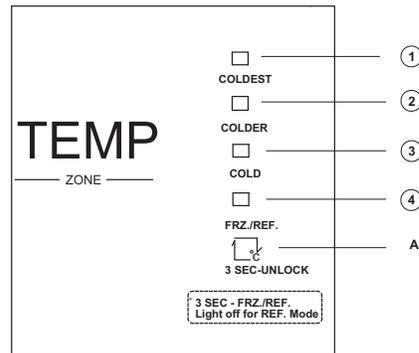1. Coldest
 2. Colder
 3. Cold
 4. Freezer and Refrigerator
- A. Temperature setting
- Once powered on, the display screen will fully display for three seconds, and then the normal operation functions will display.

Lock/Unlock

- When the temperature control panel is unused for 15 seconds, the panel will automatically lock.
- Press the Temperature Setting key (A) for three seconds to unlock the temperature control panel, if locked. All of the LED lights will blink for two seconds when the panel is unlocked.

Temperature Setting

- Press the Temperature Setting key (A) each time to change the temperature setting.
- The settings will cycle from cold to coldest, in order.
- After 15 seconds, the temperature will change to the appropriate setting.

Switching between Freezer and Refrigerator modes

This is a convertible unit which can function as refrigerator or freezer. Allow 6 to 8 hours for change of mode to come to effect. Remove samples before initiating the change of mode to avoid loss of samples and add them after 8 hours of initiating the change of mode. Follow site protocols for loading and unloading samples.

1. Press the Temperature Setting key (A) for three seconds, the Freezer and Refrigerator (4) light will illuminate, and the unit will run as a freezer.
2. Press the Temperature Setting key (A) for another three seconds, the Freezer and Refrigerator (4) light will turn off, and the unit will run as a refrigerator.
3. Do not change the mode (Refrigerator to Freezer or vice versa) or temperature setting for at least 1 hour after unit resumes to run from any power interruption including movement of unit which requires unplugging it from power supply. This would allow unit to stabilize and run properly. Also, if there is any power interruption within 1 hour of change in mode, verify the mode after power resumes to ensure.

Restart Procedure

If unit is unplugged or turned off, allow 5 minutes before restarting or plugging it back in.

How to Save Energy

- Be sure to follow location suggestions as mentioned in the **Installation**.
- Wipe moisture from glassware or other materials before placing them in a unit.
- Don't overcrowd the unit. Too many items can increase electrical energy demand in order to keep everything cool.
- Close the door as soon as possible in hot, humid weather.
- Make certain that the door is closed tightly.
- Keep containers covered, when possible, to reduce moisture buildup.
- Set operating temperature no colder than necessary for the items being refrigerated.

Safety Tips

CAUTION: After a unit is in operation, do not touch the cold surfaces, particularly when hands are damp. Skin may adhere to the cold surfaces.

WARNING: We recommend handling samples by wearing gloves to avoid frost bite.

WARNING: Never disconnect your unit by pulling the power cord. Always grip the plug securely and pull straight out from the outlet.

WARNING: Do not use a power cord that shows cracks or abrasions. Have a qualified electrician repair or replace damaged cords immediately.

Maintenance

Unit should be cleaned at least once a year following instructions below.

CAUTION: When servicing the unit, disconnect from the electrical power source.

CAUTION: Refer servicing to qualified personnel.

Note: Follow site protocols for cleaning the unit.

Cleaning the Unit

Before beginning cleaning unit follow these instruction listed below:

- Disconnect power cord from its outlet.
- The unit designs permit easy and rapid cleaning and should not take more than a few minutes. Remember to wear protective gloves to prevent frost bite, especially when removing items from freezer units.
- Do not use abrasive scouring powders, waxes, solvents, furniture polish, undiluted detergents or cleansers containing petroleum products on the surfaces of units.

Cleaning Interior/Exterior/Door Gaskets

A solution of mild soap/detergent and lukewarm water can be used for cleaning the interior, exterior and door gaskets with a soft, clean cloth. Rinse with clean water and dry thoroughly before reconnecting and turning on the unit.

The cabinet interior should be cleaned frequently. Any spilled liquid should be wiped off immediately since stains resulting from some spills could be permanent if not quickly removed. The most convenient time to clean the interior is after defrosting.

CAUTION: Solutions used to clean coils or neutralize bacteria growth must not be corrosive to metals (enamel-coated steel) and materials used in the maintenance of this equipment— damage can result: use a soft cloth and warm water to clean.

Cleaning the Condenser

With forced-fan vented units, remove the screws that mount the grill to the unit. Pull temperature control knobs straight out. This will expose condenser for cleaning.

These surfaces may be warm to the touch. The condenser should be cleaned before becoming clogged with dirt/dust. Construction or other dirt causing environments may significantly increase the required frequency. The condenser should be cleaned at least once a year.

CAUTION: Never clean around the condensers with your fingers. Some surfaces are sharp.

Transportation and Storage

Use all required local facility procedures and PPE while transporting and storing the product.

Troubleshooting

CAUTION: DO NOT try to repair your freezer yourself. Doing so invalidates the warranty.

For model FBV18CPSA

Warning Lights

The following warning lights appearing on your display indicate corresponding faults within the refrigerator and freezer. The refrigerating and freezing functions may still operate when these indicators appear, though you should contact a qualified technician if these faults appear to ensure normal refrigerating and freezing functions.

Indicator Lights

In the case of a program failure, the following LED indicator lights will blink:

Error Code FBV20RPSA, FBV20FPSA	Error Code FBV18CPSA	Error Description
Indicator light 3 and 4	Indicator light 1 and 2 are illuminated	Refrigerator temperature sensor circuit
Indicator light 2 and 4	Indicator light 1 and 3 are illuminated	Freezer defrost sensor
Indicator light 1 and 4	Indicator light 1 and 4 are illuminated	Ambient temperature sensor

General

This section is a guide to troubleshoot general operational problems.

Problem	Probable Solution
The unit does not operate.	Make sure the unit is connected to the power supply. Ensure the voltage is not too low. Check the circuit has not been tripped.
There is an odor.	Odorous samples should be tightly closed. Clean the inside of the appliance.
The compressor will not stop running.	Long operation of the appliance is normal during the summer/in warmer weather when the ambient temperature is high. Too much samples have been added to the appliance at one time. Ensure samples are cooled before adding to the appliance. The door is opened too frequently.
The light does not turn on.	Make sure the unit is connected to the power supply. Replace the light bulb.
The doors will not close properly.	The unit is blocked by samples. The unit is not properly levelled.
There is a loud noise coming from the unit.	Check whether the floor is level, and whether the unit is placed stably.
The door does not seal tightly.	Remove any foreign matter from the door seal. Heat the door seal, and then cool it for restoration.
The water pan overflows.	There is too much samples in the unit, The doors are not closed properly, resulting in frosting due to the entry of air and increased water due to defrosting.
The unit housing is hot.	Heat dissipation of the built-in compressor, via the housing, which is normal. The ambient temperature is too high. There is too much sample stored in the unit, or the compressor has shut down, providing for insufficient heat dissipation.
There is condensation on the surface of the unit.	When the ambient temperature is high, condensation on the outside of the appliance is normal. Wipe the condensate clean with a towel.
There is an abnormal noise coming from the unit	The compressor may produce a loud buzzing sound during operation, particularly upon start up or stop. This is normal. Refrigerant flowing inside of the unit may produce a creaking sound, which is normal.

Wiring Diagram

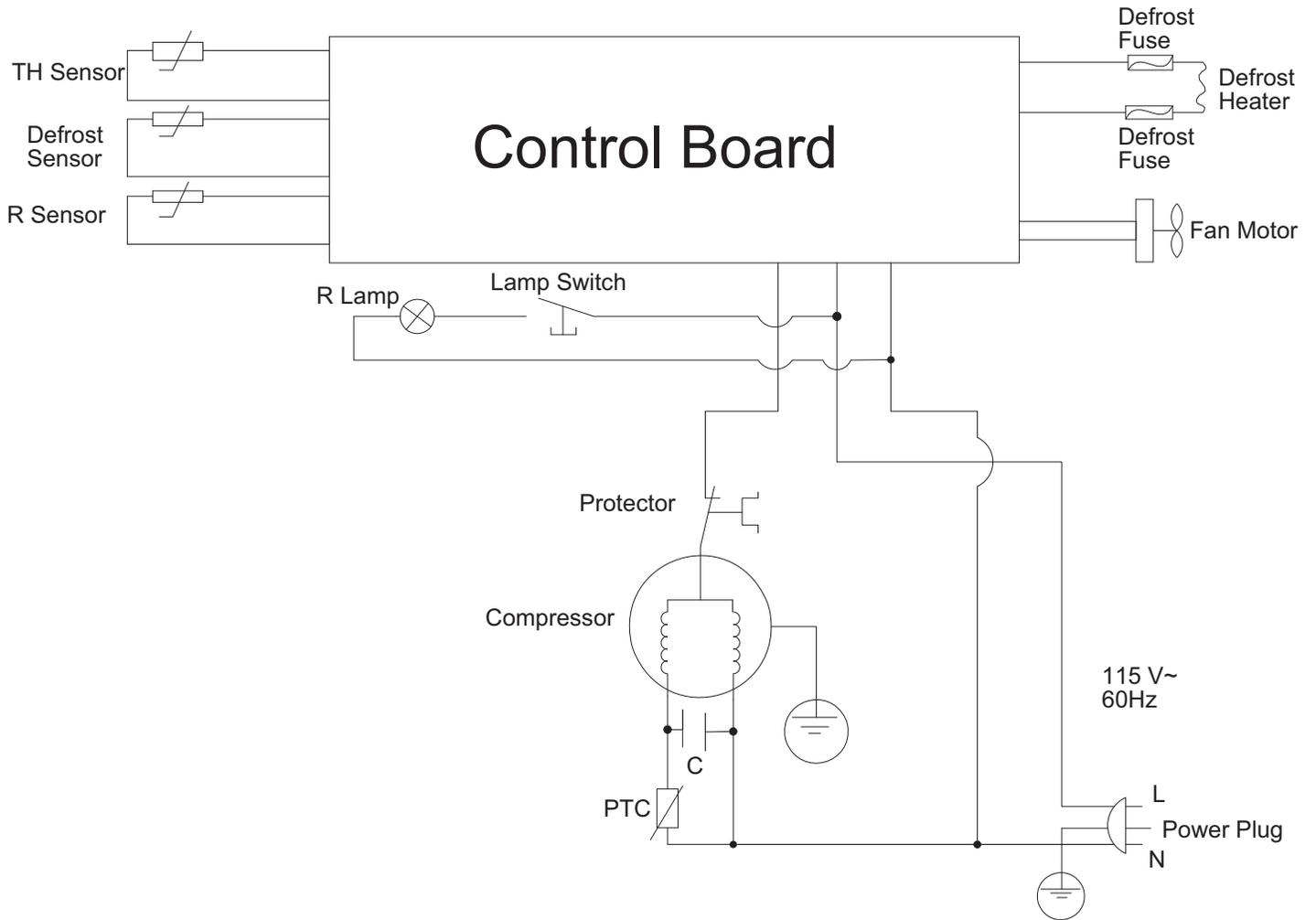
Figure 4. Wiring Diagram for FBV18CPSA model

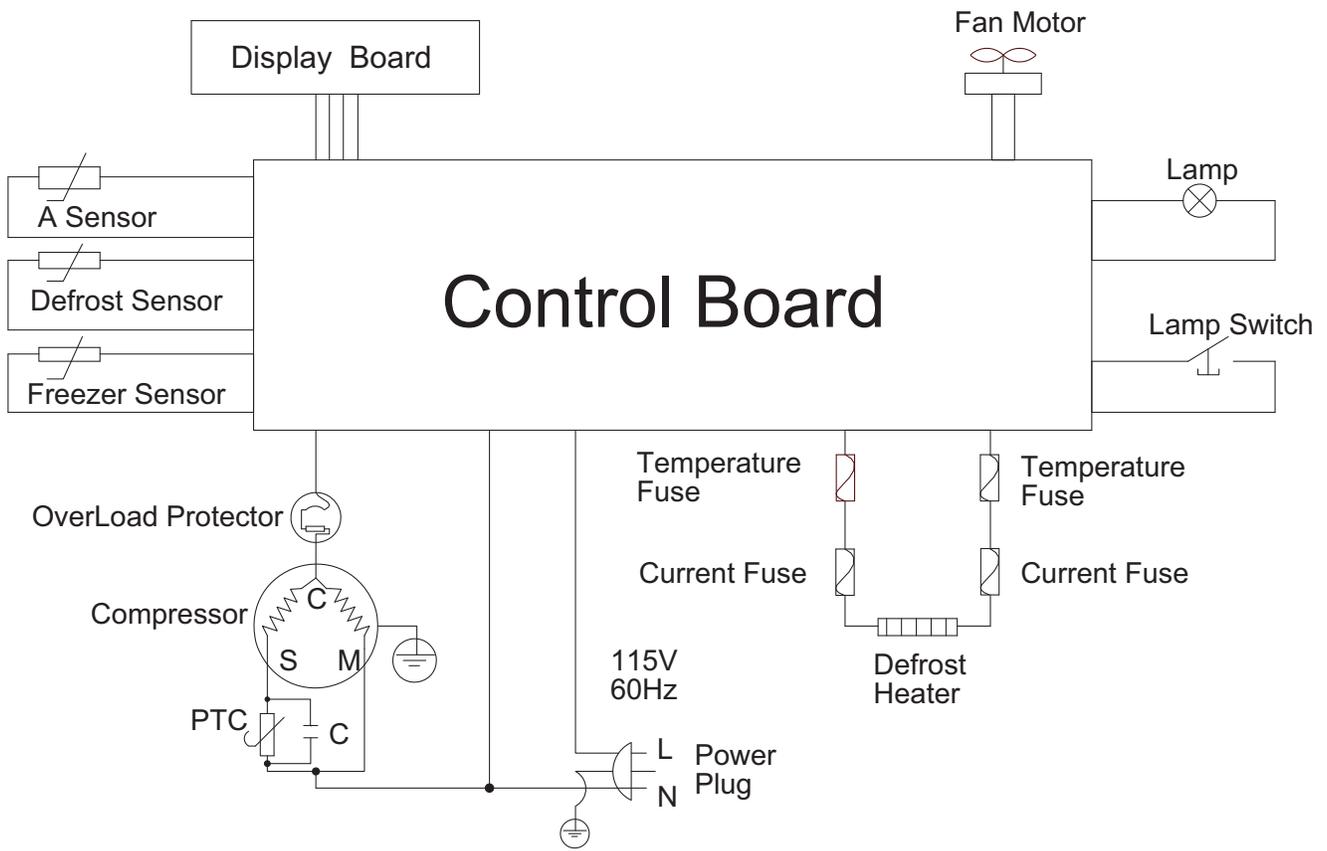
Figure 5. Wiring Diagram for FBV20RPSA and FBV20FPSA model

End of Life Care

Be sure to follow local regulations when disposing of an old unit. Some suggestions are listed below:

1. Remove items and defrost unit. Be sure to clean up any biological safety hazards.
2. Remove the cabinet door to help prevent entrapment inside of a unit.
3. Have a certified technician remove the refrigerant and compressor, then drain the compressor and oil from the system. Disposal of the unit should be done by following local regulations.

Warranty

Laboratory instruments and equipment manufactured by Fisher Scientific Company L.L.C. – Laboratory Equipment Division (hereinafter called “the Company”) are warranted only as stated below.

Subject to the exceptions and upon the conditions specified below, the Company agrees, at its election, to correct by repair, by replacement, or by credit to the purchaser, any defect of materials or workmanship which develops within one year (13 months for refrigerator and freezer products) from the date of purchase by the original purchaser by the Company or by an authorized dealer of the Company provided that investigation or factory inspection by the Company discloses that such defect developed under normal and proper use

The exceptions and conditions mentioned above are the following:

- a. The Company makes no warranty concerning components or accessories not manufactured by it, such as tubes, batteries, etc. However, in the event of the failure of any component or accessory not manufactured by the Company, the Company will give reasonable assistance to the purchaser in obtaining from the respective manufacturer whatever adjustment is reasonable in the light of the manufacturer’s own warranty.
- b. The Company shall be released from all obligations under its warranty in the event repairs or modifications are made by persons other than its own service personnel or authorized dealer personnel unless such repairs by others are made with the written consent of the Company.
- c. **THE COMPANY MAKES NO WARRANTY OF MERCHANTABILITY, FITNESS FOR ANY PARTICULAR PURPOSE, OR ANY OTHER WARRANTY, EXPRESS OR IMPLIED, EITHER IN FACT OR BY OPERATION OF LAW,...STATUTORY OR OTHERWISE.**
- d. The above warranty and the above obligations to repair, replace, or credit are complete and exclusive and the Company expressly disclaims liability for lost profits or for special, indirect, incidental, consequential, or exemplary damages of any nature whether attributable to contract, warranty, negligence, strict liability, or otherwise even if the Company has been advised of the possibility of such damages.

- e. Representations and warranties made by any person, including dealers and representatives of the Company, which are inconsistent or in conflict with the foregoing warranty shall not be binding upon the Company unless reduced to writing and signed by an officer of the Company.

Important

For your future reference and when contacting the factory, please have the following information readily available:

Model Number: _____
Serial Number: _____
Date Purchased: _____

The above information can be found on the dataplate attached to the equipment. If available, please provide the date purchased, the source of purchase (manufacturer or specific agent/rep organization), and purchase order number.

If you need assistance:

Fisher Scientific products are backed by a global technical support team ready to support your applications. We offer cold storage accessories, including remote alarms, temperature recorders, and validation services.

Visit fishersci.com or call:

SALES DIVISION	
Phone	1-828-658-2711
	1-800-252-7100
FAX	1-828-645-3368
LABORATORY PARTS and SERVICE	
Phone	1-800-438-4851
FAX	1-828-658-2576
TECHNICAL SUPPORT	
Phone	1-800-438-4851

Regulatory Compliance

Product Safety

Product Testing

This product family has been tested to applicable product safety standards by Underwriters Laboratories (UL), which is a Nationally Recognized Test Laboratory (NRTL). The product bears the following mark

Hydrocarbon Refrigerants

According to U.S. Code of Federal Regulation 40 Part 82, this freezer employs the natural hydrocarbon refrigerant R600a. Because of the nature of hydrocarbon refrigerants, for mechanical repair, such as recharge or compressor replacement, should only be carried out by a certified refrigeration technician. The safety of this equipment is listed by Underwriter Laboratory (UL) under Standard UL471, section SB – “natural refrigerant”.

Electromagnetic Compatibility

FCC Statement (USA)

Any changes or modifications not expressly approved by the party responsible for compliance could void the user’s authority to operate the equipment.

Note: The equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

Canadian ISED IC Notice

This ISM digital apparatus complies with Canadian ICES-001, Class A

This page is
intentionally
left blank.

This page is
intentionally
left blank.

This page is
intentionally
left blank.

Contact Information

Find out more at fishersci.com

In the United States:

Order online: fishersci.com

Fax an order: 1-800-926-1166

Call customer service: 1-800-766-7000

In Canada:

Order online: fishersci.ca

Fax an order: 1-800-463-2996

Call customer service: 1-800-234-7437

© 2021 Fisher Scientific Inc. All rights reserved. Trademarks used are owned as indicated at fishersci.com/trademarks. 331950H03 0721

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