

2.5 Lid Unlock



1. Switch the centrifuge on through the main on the left side of the unit.
2. When the centrifuge is switched on but stationary: LID led on at the opening.
3. The mechanic unlocks and consequent opening of the lid is obtained pulling the lever on the top right side of the machine towards the centrifuge front.

Warning: before proceeding to the installation operations, it is advisable to disconnect the centrifuge from the mains, by moving the POWER switch on “off” position.


4. After opening the lid, remove any possible packaging around the driving spindle. Unscrew the equipment-blocking nut from the engine shaft. Carefully clean the inside of the centrifugation chamber removing any packing residues. Due to the air - whirling, solid particles accidentally left in the centrifugation chamber would create an excessive wear of the chamber itself and of the outer Equipment rotating bottoms.

2.6 Emergency lid unlock



In the event of mains non-availability or power failure, the lid can be opened using the special tools supplied with the centrifuge.

This operation should always be effected with the centrifuge mains plug disconnected.

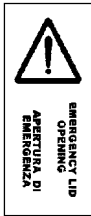

1. Unscrew the nylon screw in the hole on the right side of the centrifuge and here below identified.
2. Insert horizontally the unlocking tool supplied with the centrifuge in the hole (fig. a).
3. Lift the unlocking tool and at the same time pull the lever (fig. b).
4. At the end screw again the nylon screw.

Warning - Should the emergency occur due to electricity failure, during centrifugation, wait at least 40 minutes for the end of the Rotating Equipment rotation. Upon opening, lift the lid by hand and observe the rotor (be careful in this operation): if it is still rotating, close the lid and wait 10 minutes more before repeating the operation.

