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### 3.5 Check of Acceleration Rate

Whenever the rotor #3765 and #8171 (only LABO-FUGE 400 / 400R) is set, the microprocessor will check the acceleration rate to protect the drive system or other types from damage in the event of false speed or RCF programming (see 11.1 E-18).

**NOTE!** With this rotor setting slow acceleration will not be accepted (E-18).

### 3.6 Emergency Lid Release

#### ATTENTION - DANGER!

It is very dangerous to open the lid while the rotor is spinning!

In the event of a failure of the electrical supply, the automatic lid release is rendered inoperative.

To remove the samples when the rotor has come to **standstill**, a manually operated release is provided.

Performance as follows:

- look through the lid window to ensure rotor is not spinning
- **turn off the power, remove mains plug**
- remove plastic stopper in the center underneath the front casing using a screwdriver, spatula or knife
- pull the stopper and its attached rip cord downwards (see Fig. 2) to unlock the lid mechanically

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### 3.7 Overtemperature Protection

The LABOFUGE 400R is equipped with overtemperature protection. If the temperature in the display exceeds 50 °C, the microprocessor will indicate an overtemperature alarm (E-07, see 11.1).