



Always set the mains switch to the “O” position before performing this manual procedure, even in the case of a mains power outage. ▲

Should the emergency occur due to electricity failure during centrifugation, wait at least 10 minutes for the rotating equipment to stop spinning. In spite of the absence of noise, the rotor may still be rotating when you begin to open the lid manually.

- Insert the unlocking tool (or a flat screwdriver) horizontally into the hole on the right hand side of the instrument by piercing the protection sticker
- By using the tools, turn the gear inward until the lid latch is disengaged.
- Upon opening, lift the lid by hand and observe the rotor (be careful with this operation): if it is still rotating, close the lid and wait.



The protection sticker must be replaced after it has been pierced (item 85241801)

