

# KNOWLEDGE SHARING

## Question/Issue/Symptom/Problem:

How to change alarm setpoints on TSX High Performance Refrigerators and Freezers.

## Resolution/Answer:




Follow Steps below:



1. Turn the key to the power on position.
2. Enter the programming mode by pressing the Scan button and holding for 5 seconds. The display will then flash "Prg" to indicate that you have entered the programming mode (see table below).
3. Release Scan button. The first number is the control setpoint.
4. Press scan button once to move through menu. The next setting is cold alarm setpoint. Adjust up and down arrow to change.
5. Press scan button again to move on to warm alarm setpoint. Use up and down arrow to adjust.
6. Push scan button one more time to exit program menu.

*Note: If keypad is untouched for 30 seconds while in program (PRG) mode, the PRG mode shall be exited.*

**Table 9. Setpoint Programming Functions**

Screen/Button		Function	Summary
1	Initial Screen	Unit Setpoint	The temperature that the unit is set to run
2	Press Scan (  )	Cold Alarm Temperature	The temperature that will actuate the cold alarm (if active)
3	Press Scan (  )	Warm Alarm Temperature	The temperature that will actuate the warm alarm (if active)
4	Press Scan (  )	Exit Program Mode	Returns to normal operating screen

**Title:**

Instructions: TSX High Performance Refrigerators and Freezers - How to Change Alarm Setpoints

**Audience:**

Public

**Author:**

Alex Franklin

---