


Instruction Sheet


Control Rate Freezer Reports

Description here: How to run a report on the New Control Rate Freezer


New CRF Models: TSCM
Revision Date: July 23, 2021
Created By: Rick Pressley

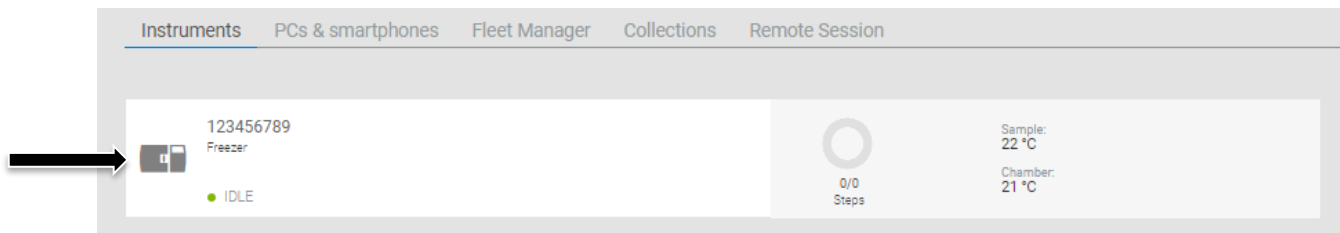

Choose the Instrument you want to Run the report on and Click the Instrument.

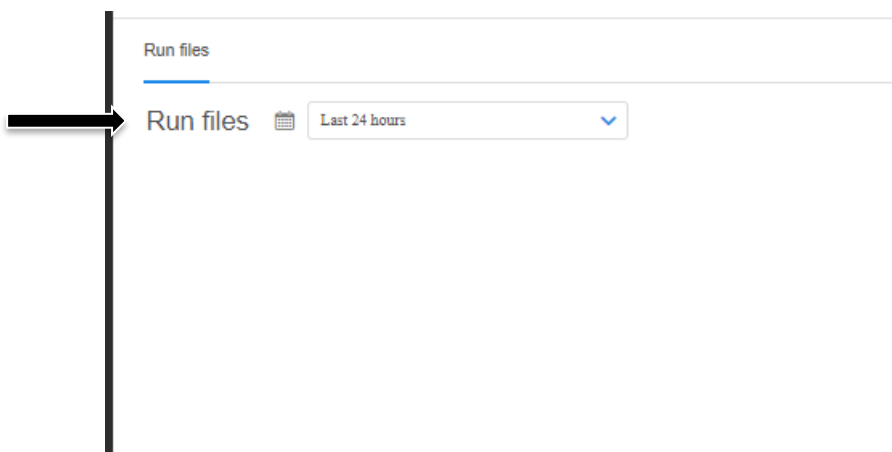


This will Bring up the run page. Choose the Date range for the report.

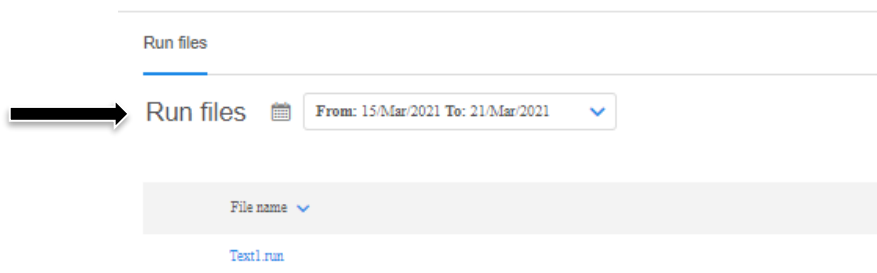


Then Choose the Date of the report.


Run files

Run files  From: 15/3/2021 To: 21/3/2021 

Compare Rates 

File name	File creation date	Download
Text1.run	Mar 18, 2021 - 11:03 AM	 Download

Click on Download for the report you want to View

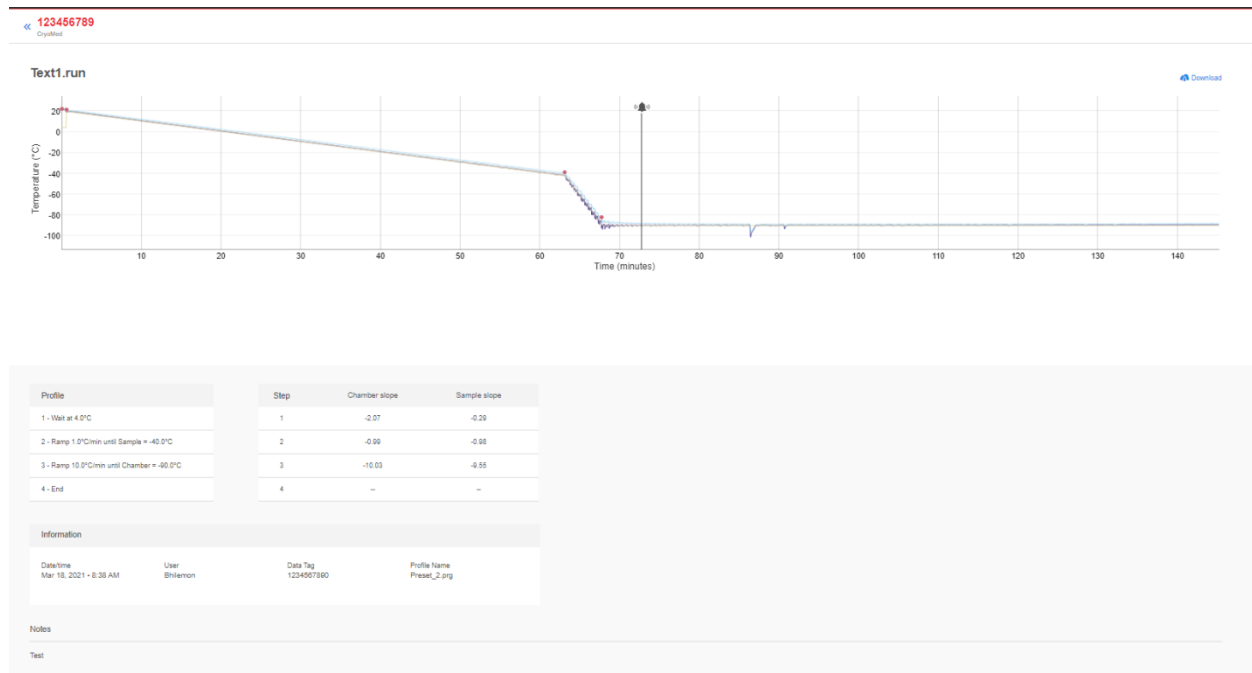

After downloading open the report.