

Instruction Sheet.

Heracell Telemetry Settings


Description here: How to set up the Telemetry Cable Set for Heracell Incubator

TSX, STP, PEEK, Heracell, Vios
Revision Date: April 25, 2023
Created By Rick Pressley

328928H03 (DLCPTDATA103) 150i /240i




The BAUD Rate needs to be changed on the Incubator and the DLC to 9600
Follow this path to the Interface settings. Use the plus and minus button to change the Baud rate and then press enter.

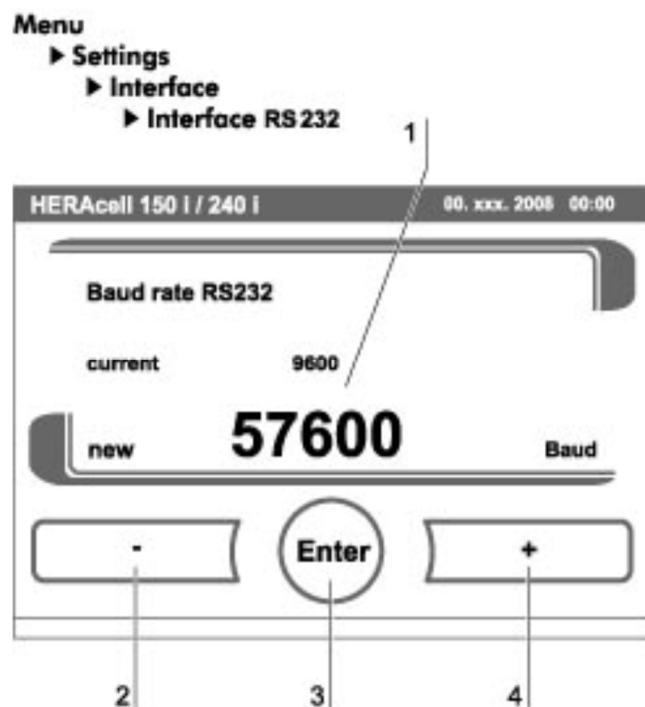


Change Baud rate to 9600

From the menu on the left go to sensors,
Add the sensor type, the USB that it is in, the baud rate, and the voltage of the unit, then submit.

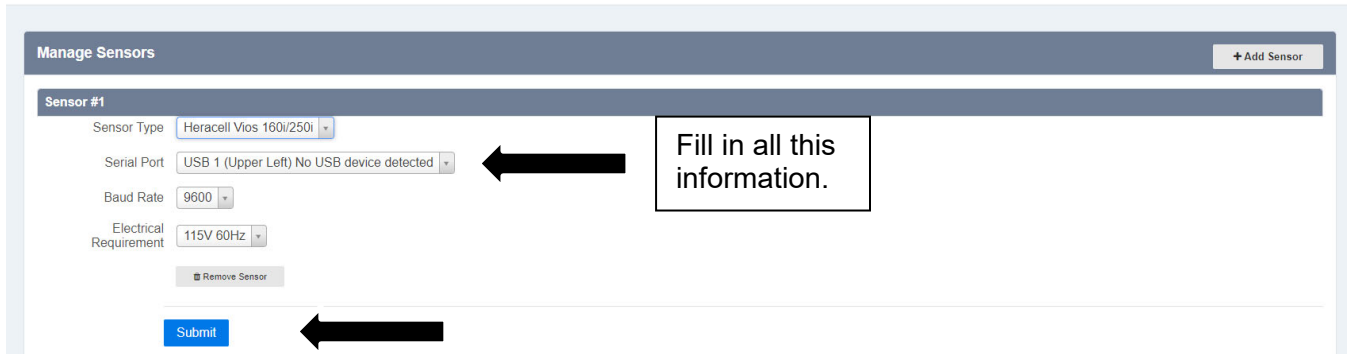

Then add your account email and the serial number of the device, then click deploy.

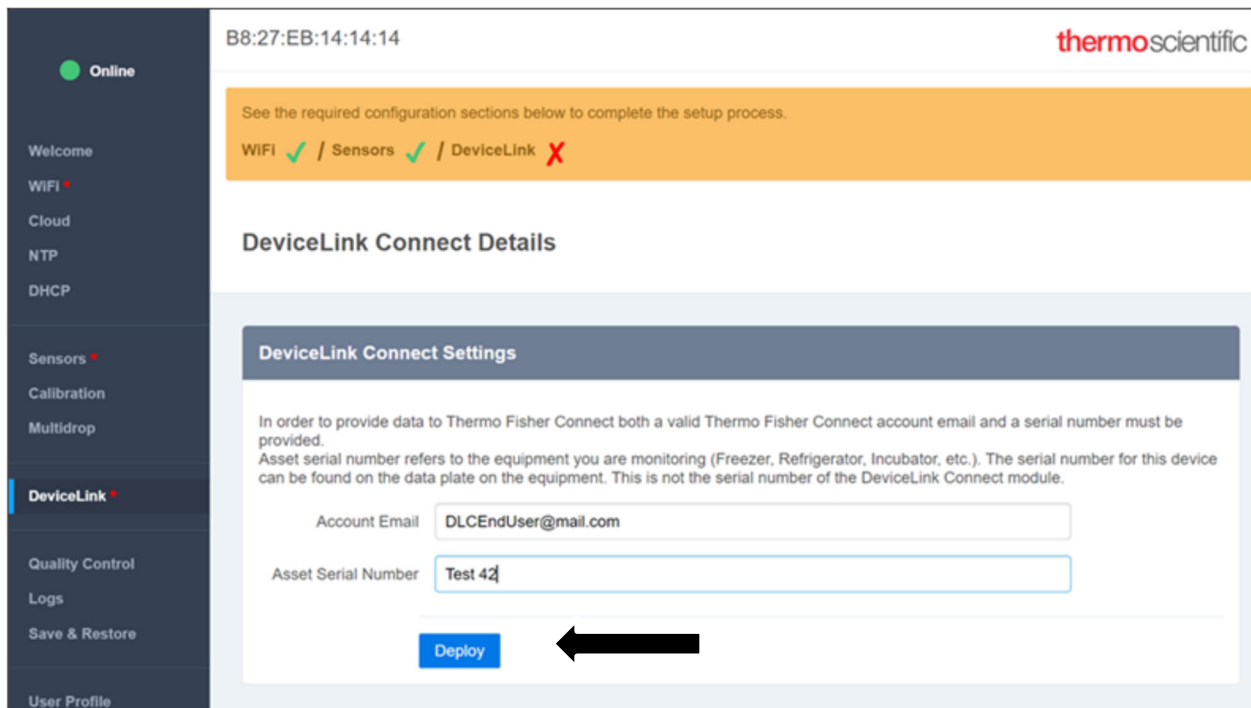

Note: Any time something is changed in the settings (WIFI, Sensors) you must deploy for the changes to be saved.