

B 39432 H01

-140/150°
CFC

REVCO INSPECTION AND ENGINEERING APPROVAL MUST BE OBTAINED PRIOR TO INITIAL PRODUCTION RUN.

IF THE SPECIFICATIONS OR TOLERANCES NOTED ON THE DRAWING CAN NOT BE MET, IMMEDIATELY NOTIFY REVCO ENGINEERING DEPARTMENT IN WRITING.

THIS DRAWING IS FURNISHED WITH THE UNDERSTANDING THAT THE ESSENCE THEREOF WILL NOT BE REPRODUCED IN WHOLE OR IN PART, WITHOUT WRITTEN AUTHORIZATION BY REVCO. ALL DESIGNS ORIGINATED BY REVCO ARE THE PROPERTY OF SAID COMPANY AND WILL BE PROTECTED BY PATENTS.


GS REVCO/LINDBERG
ASHEVILLE N.C.

TITLE WIRING DIAGRAM

SCALE NONE DIST. F B 39432 H01

08/28/95	05/18/93	DATE	TOLERANCES UNLESS OTHERWISE SPECIFIED	
REV. TERM STRIP #3; ADDED STRIP 1&2 BY STAR/LIN CAP'S	AMBIENT J7 HARNESSE ADDED	REVISONS	DECIMALS .000 ± .010 .00 ± .03 ANGULAR ±	
			DO NOT SCALE DRAWING	
			DRAWN	MGM 01/26/95
			CHECKED	BME 06/19/95
16942	17132	CHANGE	APPROVED	BME 06/19/95
B	A	LETTER	RELEASED	16955

MGM *SMH* JER

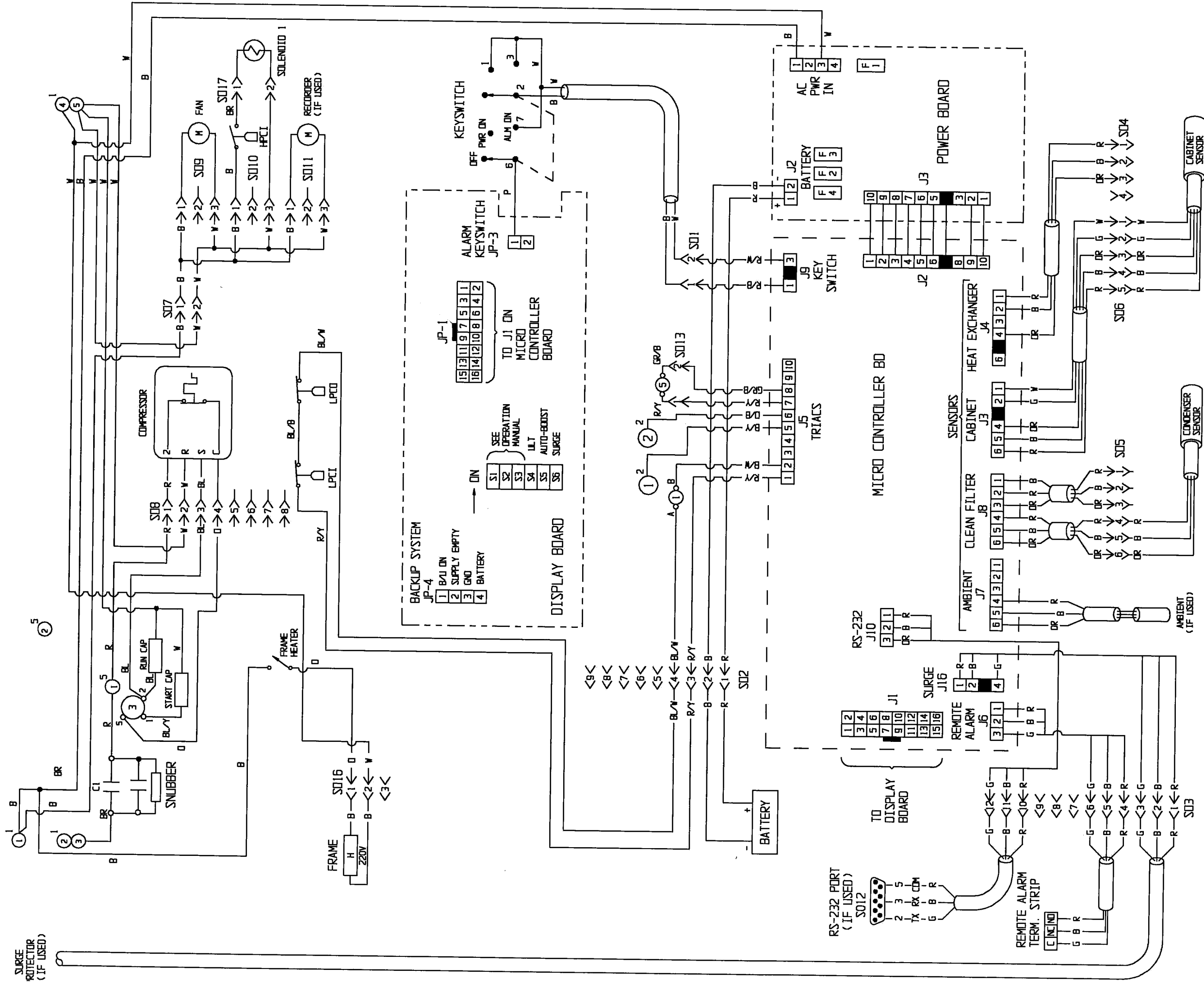

WIRES NOT CONNECTED
 WIRES CONNECTED

① TERM. STRIP # OUTSIDE IS STRIP #, INSIDE IS LUG #
 Ⓞ RELAY COIL # OUTSIDE IS TERM #, INSIDE IS RELAY #
 ○ LARGER # IS RELAY #
 ○ SMALLER # IS TERM. #

NO PRESS. SW. CLOSSES ON DROP
 NC PRESS. SW. OPENS ON RISE

SOLENOID

MOTOR
 HEATER

LEGEND:
 R - RED
 BL - BLUE
 GR - GREY
 Y - YELLOW
 O - ORANGE
 T - TAN
 B - BLACK
 BR - BROWN
 G - GREEN
 W - WHITE
 P - PURPLE
 / - SIGNIFIES WIRE TRACE

B 39432 HO1B
 WIRING DIAGRAM
 CFC -140/-150°