

# Instruction Sheet

## *How to Replace a Battery in Chest Freezer(s)*


How to Replace an Alarm Battery in Chest Freezer(s)

Chest Freezer (s)
<b>Revision Date: February 5, 2018</b>
Created by: Wayne Anders

## All Models Except 3 cu ft

The following instructions describe the battery replacement procedure for specific models.


1. Open the front door by grasping the handle and pulling.
2. Locate the battery power switch (Figure 5-1). Turn the battery power switch to the Off position (O).
3. Remove the four screws holding the recorder bezel to gain access to the battery. If you do not have a recorder, remove the blank panel above the filter panel by pulling straight out on the panel, and then remove the four screws holding the dummy panel in place where the recorder would have been installed.
4. Remove the three nuts securing the battery bracket. See Figure 5-1.
5. Remove the bracket and old battery. Discard properly. Install the new battery and secure.
6. Reconnect the battery (red to positive and black to negative).
7. Replace the recorder bezel.
8. Turn the battery power switch to Standby mode.
9. Close lower panel door.

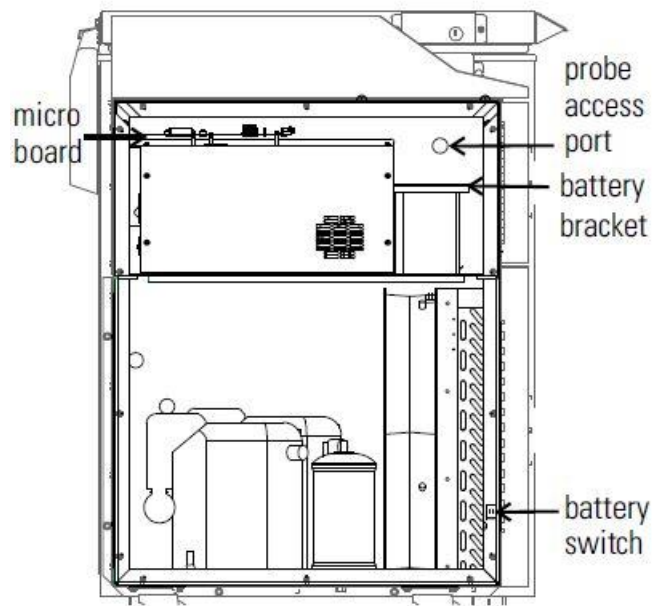

**Figure 5-1.** Battery and Switch location

### 3 cu ft Models

1. Open the front door by grasping the handle and pulling.
2. Locate the battery power switch (Figure 5-2). Turn the battery power switch to the Off position (O).
3. Remove the two nuts securing the battery bracket. See Figure 5-2.
5. Remove the bracket and old battery. Discard properly. Install the new battery and secure.
6. Reconnect the battery (red to positive and black to negative).
7. Replace the recorder bezel.
8. Turn the battery power switch to Standby mode ( ).
9. Close lower panel door.

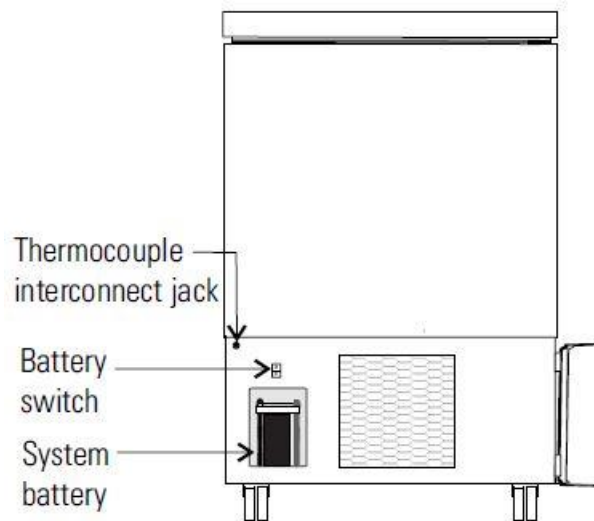

**Figure 5-2.** Battery and Switch location