


KNOWLEDGE SHARING

Question/Issue/Symptom/Problem:

How to check event log on PEEK Series ULT Freezers.

Resolution/Answer:

Follow Steps below:

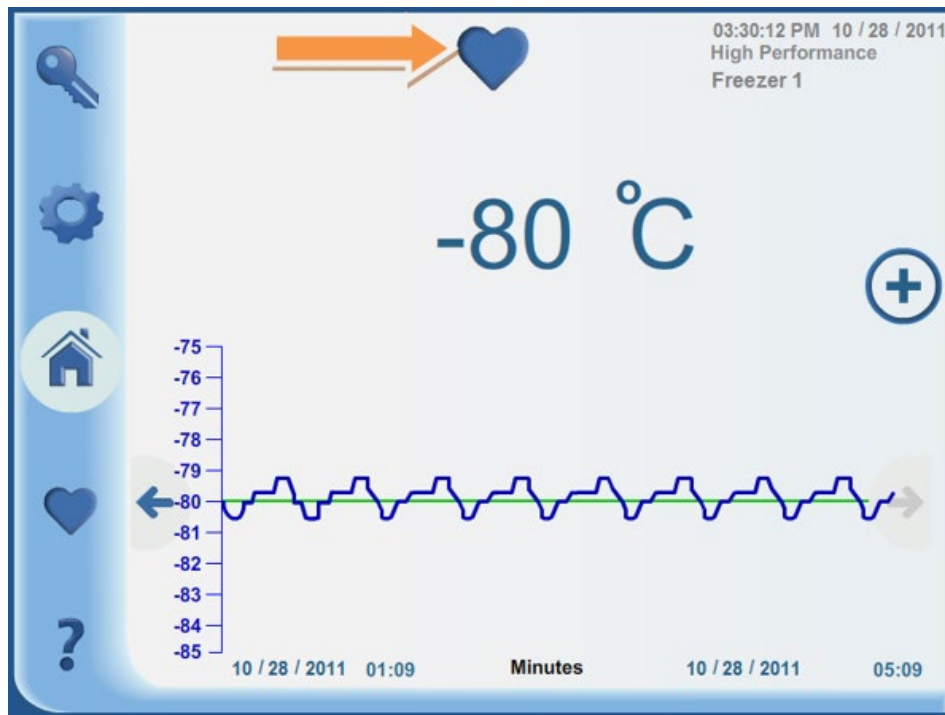

1. Press blue, yellow, or red heart on home screen to gain access the Health System Status Overview.



2. Select event notepad with pen icon to bring up extensive list of events per date and time.

Date	Time	Event
05/17/2011	05:12:01	Door Close
06/22/2011	03:20:05	System battery PM Reset.
06/23/2011	05:22:01	UI/Main Lost Communications
08/23/2011	03:14:05	System battery PM Expired.
08/25/2011	02:17:01	Power Failure Active
08/30/2011	11:12:01	Door Close
09/06/2011	11:15:01	Door Open
09/08/2011	09:22:01	Set point Settings Change
09/09/2011	05:14:01	Warm Alarm Setting Change
10/02/2011	06:32:01	UI/Main Lost Communications

3. The event log table will pull on screen. All items can be viewed, or icon filter can be used to narrow certain events. Example: Press door icon which will then filter event table to just door opening activity.

Title:

Instructions: PEEK Series ULT Freezers - How to Check Event Log

Audience:

Public

Author:

Alex Franklin
