

OMNIC Service Pack Release Notes

About This Document	11
OMNIC Service Pack (Nov-2021)	11
<i>Updated Products</i>	<i>11</i>
<i>Added Products.....</i>	<i>11</i>
<i>Deleted Products.....</i>	<i>11</i>
OMNIC Service Pack (23-July-2021).....	11
<i>Updated Products</i>	<i>11</i>
<i>Added Products.....</i>	<i>11</i>
<i>Deleted Products.....</i>	<i>11</i>
OMNIC Service Pack (28-May-2021).....	11
<i>Updated Products</i>	<i>11</i>
<i>Added Products.....</i>	<i>12</i>
<i>Deleted Products.....</i>	<i>12</i>
OMNIC Service Pack (10-May-2021).....	12
<i>Updated Products</i>	<i>12</i>
<i>Added Products.....</i>	<i>12</i>
<i>Deleted Products.....</i>	<i>12</i>
OMNIC Service Pack (26-Mar-2021)	12
<i>Updated Products</i>	<i>12</i>
<i>Added Products.....</i>	<i>12</i>
<i>Deleted Products.....</i>	<i>12</i>
OMNIC Service Pack (26-Feb-2021)	12
<i>Updated Products</i>	<i>12</i>
<i>Added Products.....</i>	<i>13</i>
<i>Deleted Products.....</i>	<i>13</i>
OMNIC Service Pack (12-Oct-2020)	13
<i>Updated Products</i>	<i>13</i>
<i>Added Products.....</i>	<i>13</i>
<i>Deleted Products.....</i>	<i>13</i>
OMNIC Service Pack (31-Aug-2020)	13
<i>Updated Products</i>	<i>13</i>

Added Products 13

Deleted Products 13

OMNIC Service Pack (17-Jul-2020) 13

Updated Products 13

Added Products 14

Deleted Products 14

OMNIC Service Pack (13-Jul-2020) 14

Updated Products 14

Added Products 14

Deleted Products 14

OMNIC Service Pack (29-May-2020) 14

Updated Products 14

Added Products 14

Deleted Products 14

OMNIC Service Pack (11-May-2020) 14

Updated Products 14

Added Products 14

Deleted Products 15

OMNIC Service Pack (10-Apr-2020) 15

Updated Products 15

Added Products 15

Deleted Products 15

OMNIC Service Pack (05-Feb-2020) 15

Updated Products 15

Added Products 15

Deleted Products 15

OMNIC Service Pack (22-Nov-2019) 15

Updated Products 15

Added Products 15

Deleted Products 15

OMNIC Service Pack (01-Nov-2019) 16

Updated Products 16

Added Products 16

Deleted Products 16

OMNIC Service Pack (18-Oct-2019) 16

Updated Products 16

Added Products..... 16

Deleted Products..... 16

OMNIC Service Pack (30-Sep-2019) 16

Updated Products 16

Added Products..... 16

Deleted Products..... 17

OMNIC Service Pack (03-Sep-2019) 17

Updated Products 17

Added Products..... 17

Deleted Products..... 17

OMNIC Service Pack (26-Aug-2019) 17

Updated Products 17

Added Products..... 17

Deleted Products..... 17

OMNIC Service Pack (26-Jul-2019) 17

Updated Products 17

Added Products..... 18

Deleted Products..... 18

OMNIC Service Pack (12-Jul-2019) 18

Updated Products 18

Added Products..... 18

Deleted Products..... 18

OMNIC Service Pack (14-Jun-2019) 18

Updated Products 18

Added Products..... 19

Deleted Products..... 19

OMNIC Service Pack (15-Apr-2019) 19

Updated Products 19

Added Products..... 19

Deleted Products..... 19

OMNIC Service Pack (26-Feb-2019) 19

Updated Products 19

Added Products 20

Deleted Products 20

OMNIC Service Pack (12-Feb-2019) 20

Updated Products 20

Added Products 20

Deleted Products 20

OMNIC Service Pack (18-Jan-2019) 20

Updated Products 20

Added Products 20

Deleted Products 20

OMNIC Service Pack (30-Nov-2018) 20

Updated Products 20

Added Products 21

Deleted Products 21

OMNIC Service Pack (21-Nov-2018) 21

Updated Products 21

Added Products 21

Deleted Products 21

OMNIC Service Pack (06-Nov-2018) 21

Updated Products 21

Added Products 21

Deleted Products 21

OMNIC Service Pack (11-Oct-2018) 21

Updated Products 21

Added Products 22

Deleted Products 22

OMNIC Service Pack (19-Sep-2018) 22

Updated Products 22

Added Products 22

Deleted Products 22

OMNIC Service Pack (14-Sep-2018) 22

Updated Products 22

Added Products 23

Deleted Products 23

OMNIC Service Pack (24-Jul-2018) 23

Updated Products 23

Added Products..... 23

Deleted Products..... 23

OMNIC Service Pack (Jun-2018) 23

Updated Products 23

Added Products..... 23

Deleted Products..... 23

OMNIC Service Pack (11-Apr-2018) 23

Updated Products 23

Added Products..... 24

Deleted Products..... 24

OMNIC Service Pack (21-Feb-2018) 24

Updated Products 24

Added Products..... 24

Deleted Products..... 24

OMNIC Service Pack (2-Feb-2018)..... 24

Updated Products 24

Added Products..... 24

Deleted Products..... 24

OMNIC Service Pack (12-Jan-2018) 24

Updated Products 24

Added Products..... 25

Deleted Products..... 25

OMNIC Service Pack (8-Nov-2017) 25

Updated Products 25

Added Products..... 25

Deleted Products..... 25

OMNIC Service Pack (3-Nov-2017) 25

Updated Products 25

Added Products..... 26

Deleted Products..... 26

OMNIC Service Pack (31-Aug-2017) 26

Updated Products 26

Added Products 26

Deleted Products 26

OMNIC Service Pack (04-Aug-2017) 26

Updated Products 26

Added Products 27

Deleted Products 27

OMNIC Service Pack (29-Jun-2017) 27

Updated Products 27

Added Products 27

Deleted Products 27

OMNIC Service Pack (20-Jun-2017) 27

Updated Products 27

Added Products 27

Deleted Products 27

OMNIC Service Pack (14-Jun-2017) 27

Updated Products 27

Added Products 28

Deleted Products 29

OMNIC Service Pack (20-Apr-2017) 29

Updated Products 29

Added Products 29

Deleted Products 29

OMNIC Service Pack (10-Mar-2017) 29

Updated Products 29

Added Products 30

Deleted Products 30

OMNIC Service Pack (13-Jan-2017) 30

Updated Products 30

Added Products 30

Deleted Products 30

OMNIC Service Pack (21-Dec-2016) 30

Updated Products 30

Added Products 30

Deleted Products 30

OMNIC Service Pack (16-Dec-2016)..... 30

Updated Products 30

Added Products..... 31

Deleted Products..... 31

OMNIC Service Pack (09-Nov-2016) 31

Updated Products 31

Added Products..... 31

Deleted Products..... 31

OMNIC Service Pack (12-Oct-2016) 31

Updated Products 31

Added Products..... 31

Deleted Products..... 31

OMNIC Service Pack (04-Oct-2016) 32

Updated Products 32

Added Products..... 32

Deleted Products..... 32

OMNIC Service Pack (12-Aug-2016) 32

Updated Products 32

Added Products..... 32

Deleted Products..... 32

OMNIC Service Pack (27-Jun-2016) 32

Updated Products 32

Added Products..... 32

Deleted Products..... 33

OMNIC Service Pack (22-Jun-2016) 33

Updated Products 33

Added Products..... 33

Deleted Products..... 33

OMNIC Service Pack (21-Jun-2016) 33

Updated Products 33

Added Products..... 33

Deleted Products..... 33

OMNIC Service Pack (31-May-2016)..... 33

Updated Products 33

<i>Added Products</i>	33
<i>Deleted Products</i>	33
OMNIC Service Pack (06-May-2016)	34
<i>Updated Products</i>	34
<i>Added Products</i>	34
<i>Deleted Products</i>	34
OMNIC Service Pack (23-Feb-2016)	34
<i>Updated Products</i>	34
<i>Added Products</i>	34
<i>Deleted Products</i>	34
OMNIC Service Pack (18-Feb-2016)	34
<i>Updated Products</i>	34
<i>Added Products</i>	35
<i>Deleted Products</i>	35
OMNIC Service Pack (04-Feb-2016)	35
<i>Updated Products</i>	35
<i>Added Products</i>	35
<i>Deleted Products</i>	35
OMNIC Service Pack (12-Jan-2016)	36
<i>Updated Products</i>	36
<i>Added Products</i>	36
<i>Deleted Products</i>	36
OMNIC Service Pack (23-Nov-2015)	36
<i>Updated Products</i>	36
<i>Added Products</i>	37
<i>Deleted Products</i>	37
OMNIC Service Pack (28-Oct-2015)	37
<i>Updated Products</i>	37
<i>Added Products</i>	37
<i>Deleted Products</i>	38
OMNIC Service Pack (08-May-2015)	38
<i>Updated Products</i>	38
<i>Added Products</i>	38
<i>Deleted Products</i>	38

OMNIC Service Pack (22-Jan-2015) 38

Updated Products 38

Added Products..... 38

Deleted Products..... 38

OMNIC Service Pack (10-Nov-2014) 38

Updated Products 38

Added Products..... 39

Deleted Products..... 39

OMNIC Service Pack (02-Apr-2014) 39

Updated Products 39

Added Products..... 39

Deleted Products..... 39

OMNIC Service Pack (28-Feb-2014) 39

Updated Products 39

Added Products..... 39

Deleted Products..... 39

OMNIC Service Pack (09-Sep-2013) 40

Updated Products 40

Added Products..... 40

Deleted Products..... 40

OMNIC Service Pack (22-Jul-2013) 40

Updated Products 40

Added Products..... 40

Deleted Products..... 40

OMNIC Service Pack (15-Jun-2013) 40

Updated Products 40

Added Products..... 40

Deleted Products..... 40

OMNIC Service Pack (01-May-2013)..... 41

Updated Products 41

Added Products..... 41

Deleted Products..... 41

OMNIC Service Pack (08-Mar-2013) 41

Updated Products 41

Added Products..... 41
Deleted Products..... 41
OMNIC Service Pack (04-Oct-2012) 41
Updated Products 41
Added Products..... 42
Deleted Products..... 42
OMNIC Service Pack (09-Aug-2012) 42
Updated Products 42
Added Products..... 42
Deleted Products..... 42
How to Contact Us 42
Trademarks 43

About This Document

This document contains a revision history of the OMNIC Suite OMNIC Service Pack, including updated products, added products and Deleted Products. Change details are available within the Release Notes associated with each product.

OMNIC Service Pack (Nov-2021)

Updated Products

Service Pack Release Notes	This document
OMNIC v9.12 HF6	See OMNIC Release Notes
OMNIC 9 FTIR Universal Driver Files v9.12 HF6	See OMNIC Release Notes
OMNIC for Dispersive Raman v9.12 HF6	See OMNIC for Dispersive Raman Release Notes
OMNIC Release Notes v9.12 HF6	The OMNIC Release Notes
Raman AutoSampler v4.1 HF2	Removed support for Macros Basic

Added Products

None

Deleted Products

None

OMNIC Service Pack (23-July-2021)

Updated Products

Service Pack Release Notes	This document
OMNIC 9.12 HF5	See OMNIC Release Notes
OMNIC for Dispersive Raman 9.12 HF5	See OMNIC for Dispersive Raman Release Notes
OMNIC Release Notes v9.12 HF5	The OMNIC Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (28-May-2021)

Updated Products

Service Pack Release Notes	This document
OMNIC 9.12 HF4	See OMNIC Release Notes
OMNIC for Dispersive Raman 9.12 HF4	See OMNIC for Dispersive Raman Release Notes
OMNIC Release Notes v9.12 HF4	The OMNIC Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (10-May-2021)

Updated Products

Service Pack Release Notes	This document
OMNIC 9.12 HF3	See OMNIC Release Notes
AtIµs 9.12.HF3	See AtIµs Release Notes
OMNIC Picta 1.8	See OMNIC Picta Release Notes
ValPro Qualification JP Methods for FTIR	See ValPro Qualification JP Methods for FTIR Release Notes
ValPro Qualification PV Methods for FTIR	See ValPro Qualification PV Methods for FTIR Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (26-Mar-2021)

Updated Products

Service Pack Release Notes	This document
OMNIC 9.12 HF2	See OMNIC Release Notes
OMNIC for Dispersive Raman 9.12 HF2	See OMNIC for Dispersive Raman Release Notes
Array Automation 9.12 HF2	See Array Automation Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (26-Feb-2021)

Updated Products

Service Pack Release Notes	This document
OMNIC 9.12 HF1	See OMNIC Release Notes

OMNIC for Dispersive Raman 9.12 HF1	See OMNIC for Dispersive Raman Release Notes
Array Automation 9.12 HF1	See Array Automation Release Notes
Atlus 9.12 HF1	See Atlus Release Notes
TQ Analyst 9.12 HF1	See TQ Analyst Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (12-Oct-2020)

Updated Products

Service Pack Release Notes	This document
OMNIC 9.12	See OMNIC Release Notes
OMNIC for Dispersive Raman 9.12	See OMNIC for Dispersive Raman Release Notes
TQ Analyst 9.12	See TQ Analyst Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (31-Aug-2020)

Updated Products

Service Pack Release Notes	This document
OMNIC 9.11 HF7	See OMNIC Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (17-Jul-2020)

Updated Products

Service Pack Release Notes	This document
Picta 1.7 HF6	See OMNIC Picta Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (13-Jul-2020)

Updated Products

Service Pack Release Notes	This document
Security Suite 2.3	See Security Suite Release Notes
Audit Manager 1.3	See Audit Manager Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (29-May-2020)

Updated Products

Service Pack Release Notes	This document
Picta 1.7 HF5	See OMNIC Picta Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (11-May-2020)

Updated Products

Service Pack Release Notes	This document
OMNIC 9.11 HF6	See OMNIC Release Notes
Atlus 9.11 HF6	See Atlus Release Notes
Security Suite Plug-ins v4.10 HF1	Added logic to prevent upgrading OMNICxi from 1.x to 2.x

Added Products

None

Deleted Products

None

OMNIC Service Pack (10-Apr-2020)

Updated Products

Service Pack Release Notes	This document
OMNIC Picta 1.7 HF4	See OMNIC Picta Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (05-Feb-2020)

Updated Products

Service Pack Release Notes	This document
Atlus 9.11 HF5	See Atlus Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (22-Nov-2019)

Updated Products

Service Pack Release Notes	This document
OMNIC Picta 1.7 HF2	See OMNIC Picta Release Notes
TQ Analyst 9.11	See TQ Analyst Release Notes
OMNIC 9.11 HF4	See OMNIC Release Notes
Atlus 9.11 HF4	See Atlus Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (01-Nov-2019)

Updated Products

Service Pack Release Notes	This document
OMNIC 9.11 HF3	See OMNIC Release Notes
At μ s 9.11 HF3	See At μ s Release Notes
ValPro Qualification PHEUR Methods for FTIR v1.7	See ValPro Qualification PHEUR Methods for FTIR Release Notes
ValPro Qualification PV Methods for FTIR v1.7	See ValPro Qualification PV Methods for FTIR Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (18-Oct-2019)

Updated Products

OMNIC 9 FTIR Universal Driver Files v9.11 HF2	See OMNIC Release Notes
OMNIC v9.11 HF2	The OMNIC Release Notes
Service Pack Release Notes	This document

Added Products

None

Deleted Products

None

OMNIC Service Pack (30-Sep-2019)

Updated Products

At μ s v9.11 HF1	See At μ s Release Notes
OMNIC v9.11 HF1	See OMNIC Release Notes
OMNIC for Dispersive Raman v9.11 HF1	See OMNIC for Dispersive Raman Release Notes
OMNIC 9 FTIR Universal Driver Files v9.11 HF1	See OMNIC Release Notes
OMNIC Release Notes v9.11 HF1	The OMNIC Release Notes
Service Pack Release Notes	This document
ValPro Sys Qual FTIR Raman v2.9 HF1	See ValPro Qualification Release Notes
Visual C++ 2015-2019 x86 Redistributable v14.22	Updated to latest version from Microsoft

Added Products

None

Deleted Products

None

OMNIC Service Pack (03-Sep-2019)

Updated Products

OMNICxi Optical Component Files	See OMNICxi Optical Component Files Release Notes
OMNICxi v1.7	See OMNICxi Release Notes
Service Pack Release Notes	This document

Added Products

None

Deleted Products

None

OMNIC Service Pack (26-Aug-2019)

Updated Products

OMNIC v9.11	See OMNIC Release Notes
OMNIC for Dispersive Raman v9.11	See OMNIC for Dispersive Raman Release Notes
OMNIC Release Notes v9.11	The OMNIC Release Notes
Service Pack Release Notes	This document
Setup v4.9 HF3	Installation now sets default registry items correctly
Thermo Software IQ v3.2 HF4	See Thermo Software IQ Release Notes
TQ Analyst Data Security v9.8 HF2	See TQ Analyst Release Notes
TQ Analyst Release Notes v9.8 HF2	TQ Analyst Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (26-Jul-2019)

Updated Products

Service Pack Release Notes	This document
Thermo Software IQ v3.2 HF3	See Thermo Software IQ Release Notes
ValPro Qualification PHEUR Methods for FTIR v1.6	See ValPro Qualification PHEUR Methods for FTIR Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (12-Jul-2019)

Updated Products

Audit Log Service v1.1	See Audit Manager Release Notes
Audit Manager v1.1	See Audit Manager Release Notes
Security Suite Plug-ins v4.9 HF2	Improved diagnostic logging
Security Suite Supported Products	Added: <ul style="list-style-type: none"> • OMNIC Paradigm • TQ Analyst for OMNIC Paradigm
Service Pack Release Notes	This document
Setup v4.9 HF2	Improved diagnostic logging
Security Suite v2.2	See Security Suite Release Notes
Thermo Software IQ v3.2 HF2	Updated timestamp format to be consistent with other products
TQ Analyst v9.8 HF1	See TQ Analysts Release Notes
TQ Analyst Data Security v9.8 HF1	See TQ Analysts Release Notes
TQ Analyst - EZ Edition v9.8 HF1	See TQ Analysts Release Notes
TQ Analyst Release Notes v9.8 HF1	TQ Analysts Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (14-Jun-2019)

Updated Products

Service Pack Release Notes	This document
ValPro Qualification CP Methods for FTIR v1.4	See ValPro Qualification CP Methods for FTIR Release Notes
ValPro Qualification JP Methods for FTIR v1.4	See ValPro Qualification JP Methods for FTIR Release Notes
ValPro Qualification PHEUR Methods for FTIR v1.5	See ValPro Qualification PHEUR Methods for FTIR Release Notes
ValPro Qualification PV Methods for FTIR v1.4	See ValPro Qualification PV Methods for FTIR Release Notes
ValPro Qualification USP Methods for FTIR v1.4	See ValPro Qualification USP Methods for FTIR

	Release Notes
--	---------------

Added Products

None

Deleted Products

None

OMNIC Service Pack (15-Apr-2019)

Updated Products

OMNIC v9.9 HF3	See OMNIC Release Notes
OMNIC Data Security v9.9 HF3	See OMNIC Release Notes
Wafer Analyzer SW for OMNIC 9 v9.9 HF3	See OMNIC Release Notes
OMNIC Release Notes v9.9 HF3	The OMNIC Release Notes
Security Suite Plug-ins v4.9 HF1	Updated to work with shipping version of Setup Utility
Security Suite Supported Products	See OMNIC Release Notes
Service Pack Release Notes	This document
ValPro Qualification PHEUR Methods for FTIR v1.4	See ValPro Qualification PHEUR Methods for FTIR Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (26-Feb-2019)

Updated Products

Service Pack Release Notes	This document
ValPro Sys Qual FTIR Raman Methods v2.9 HF1	See ValPro Qualification Methods Release Notes
ValPro Qualification CP Methods for FTIR v1.3	See ValPro Qualification CP Methods for FTIR Release Notes
ValPro Qualification JP Methods for FTIR v1.3	See ValPro Qualification JP Methods for FTIR Release Notes
ValPro Qualification PHEUR Methods for FTIR v1.3	See ValPro Qualification PHEUR Methods for FTIR Release Notes
ValPro Qualification PV Methods for FTIR v1.3	See ValPro Qualification PV Methods for FTIR Release Notes
ValPro Qualification USP Methods for FTIR v1.3	See ValPro Qualification USP Methods for FTIR Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (12-Feb-2019)

Updated Products

Service Pack Release Notes	This document
OMNIC v9.9 HF2	See OMNIC Release Notes
OMNIC for Dispersive Raman v9.9 HF2	See OMNIC for Dispersive Raman Release Notes
OMNIC for Particle Analysis v9.9 HF2	See OMNIC Particle Analysis Release Notes
OMNIC Picta v1.7 HF1	See OMNIC Picta Release Notes
Atl μ s v9.9 HF2	See Atl μ s Release Note
SST PEM & TRS Experiments v9.9	See OMNIC SST Release Notes
Visualization Tools v10.0 HF3	See Support for saving maps with security installed

Added Products

None

Deleted Products

None

OMNIC Service Pack (18-Jan-2019)

Updated Products

Service Pack Release Notes	This document
OMNIC for Raman iS50 v9.9	See OMNIC for Raman iS50 Release Notes
OMNIC for Raman NXR v9.9	See OMNIC for Raman NXR Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (30-Nov-2018)

Updated Products

Service Pack Release Notes	This document
iS20 Documentation Update v1.1	Updates

Demo Instrument Config (for 9.9)	Added support for iS20 simulation
----------------------------------	-----------------------------------

Added Products

None

Deleted Products

None

OMNIC Service Pack (21-Nov-2018)

Updated Products

OMNICxi v1.6 HF1	See OMNICxi Release Notes
Service Pack Release Notes	This document

Added Products

OMNIC Cloud Connector v1.0	See OMNIC Cloud Connector Release Notes
----------------------------	---

Deleted Products

None

OMNIC Service Pack (06-Nov-2018)

Updated Products

OMNIC v9.9 HF1	See OMNIC Release Notes
OMNIC 9 FTIR Universal Driver Files v9.9 HF1	See OMNIC Release Notes
OMNIC Release Notes	Updated OMNIC Release Notes
Service Pack Release Notes	This document
Setup v4.9	Improved .NET Framework installation

Added Products

Videology Tools v1.0.151	Restored – See OMNIC Release Notes
--------------------------	------------------------------------

Deleted Products

None

OMNIC Service Pack (11-Oct-2018)

Updated Products

Atlμs v9.9 HF1	See Atlμs Release Notes
OMNIC Picta v1.7	See Picta Release Notes
Security Suite Plug-ins v4.8 HF3	Corrected issue with creating distributed media Added check to catch improper media configuration
Security Suite Supported Products	Added missing files causing failure to do distributed installs

Service Pack Release Notes	This document
Visualization Tools v10.0 HF2	Changes to support Atlµs and Picta fixes

Added Products

Particle Analysis Software v9.9	Initial release (in OMNIC for Dispersive Raman only)
---------------------------------	--

Deleted Products

None

OMNIC Service Pack (19-Sep-2018)

Updated Products

Nicolet iS5 Documentation v9.9	Brought up to current branding standard
Service Pack Release Notes	This document
Setup v4.8	Allow user to opt out of potential server reboot during distributed installation

Added Products

None

Deleted Products

None

OMNIC Service Pack (14-Sep-2018)

Updated Products

Array Automation v9.9	See Array Automation Release Notes
Atlµs v9.9	See Atlµs Release Notes
Gas App OM for Antaris IGS and iG50 SP v9.9	See OMNIC Release Notes
MicroView v9.9	See Atlµs Release Notes
OMNIC 9 FTIR Universal Driver Files v9.9	See OMNIC Release Notes
OMNIC Data Security v9.9	See OMNIC Release Notes
OMNIC for Dispersive Raman v9.9	See OMNIC for Dispersive Raman Release Notes
OMNIC MC v9.9	See OMNIC Release Notes
OMNIC Release Notes	Updated OMNIC Release Notes
OMNIC v9.9	See OMNIC Release Notes
Series Data Collect v9.9	See OMNIC Release Notes
Service Pack Plug-in v4.8	Removed ability to scan the service pack itself
Service Pack Release Notes	This document
Thermo Software IQ v3.2	Alphabetized the product list. Added universal pass/fail note at the end of the report
Wafer Analyzer SW for OMNIC 9 v9.9	See OMNIC Release Notes

Added Products

Gas Analysis for OMNIC v9.9	Initial release
-----------------------------	-----------------

Deleted Products

None

OMNIC Service Pack (24-Jul-2018)

Updated Products

None

Added Products

None

Deleted Products

qPro Developer v2.0	Product decision to not include on Service Pack
qPro Operator v2.0	Product decision to not include on Service Pack
qPro Server v2.0	Product decision to not include on Service Pack

OMNIC Service Pack (Jun-2018)

Updated Products

OMNIC Picta v1.6 HF1	See OMNIC Picta Release Notes
Visualization Tools v10.0 HF1	See OMNIC Picta Release Notes
Service Pack Release Notes	This document

Added Products

None

Deleted Products

None

OMNIC Service Pack (11-Apr-2018)

Updated Products

OMNIC 9.8 HF2 – Data Security	See OMNIC Release Notes
OMNIC 9.8 HF2 – Wafer Analyzer	See OMNIC Release Notes
OMNIC for Dispersive Raman 9.8 HF2	See OMNIC for Dispersive Raman Release Notes
OMNIC Release Notes	Updated OMNIC Release Notes
Security Suite v2.1 HF2	See Security Suite Release Notes
Security Suite Plug-ins	See Security Suite Release Notes
Audit Log Service v1.0 HF2	See Audit Manager Release Notes
Audit Manager v1.0 HF2	See Audit Manager Release Notes

Service Pack Release Notes	This document
----------------------------	---------------

Added Products

None

Deleted Products

None

OMNIC Service Pack (21-Feb-2018)

Updated Products

Audit Log Service v1.0 HF1	See Audit Manager Release Notes
Audit Manager v1.0	See Audit Manager Release Notes
Security Suite Documentation v2.1 HF1	Updated for this release
Security Suite Plug-ins v2.1	See Security Suite Release Notes
Security Suite v2.1 HF1	See Security Suite Release Notes
Service Pack Release Notes	This document

Added Products

None

Deleted Products

None

OMNIC Service Pack (2-Feb-2018)

Updated Products

OMNICxi Optical Components v1.1	See OMNICxi Optical Components Files Release Notes
Service Pack Release Notes	This document

Added Products

qPro Developer v2.0	See qPro Developer Release Notes
qPro Operator v2.0	See qPro Operator Release Notes
qPro Server v2.0	See qPro Developer and Operator Release Notes

Deleted Products

None

OMNIC Service Pack (12-Jan-2018)

Updated Products

Security Suite Documentation v2.1	Updated for this release
Security Suite Language Pack v2.1	See Security Suite Release Notes

Security Suite Plug-ins v2.1	See Security Suite Release Notes
Security Suite Supported Products	See Security Suite Release Notes
Security Suite v2.1	See Security Suite Release Notes
Service Pack Release Notes	This document
Thermo Software IQ v3.1 HF1	Updated how file sizes were calculated

Added Products

Audit Log Service v1.0	See Audit Manager Release Notes
Audit Manager v1.0	See Audit Manager Release Notes
Security Suite SP Batch ID v1.00	Needed to create proper media structure

Deleted Products

None

OMNIC Service Pack (8-Nov-2017)

Updated Products

OMNIC Picta Data Security v1.6 HF1	Fixed upgrade over 1.6
Service Pack Release Notes	This document

Added Products

None

Deleted Products

None

OMNIC Service Pack (3-Nov-2017)

Updated Products

Array Automation v9.8 HF1	See Array Automation Release Notes
Atlμs v9.8 HF1	Fixed upgrade from OMNIC Picta 1.3 and older Also see Atlμs Release Notes
Diagnostics for Picta v1.6 HF1	Fixed upgrade from OMNIC Picta 1.3 and older
Installer for Thermo Products	Fixed upgrade from OMNIC Picta 1.3 and older
Nicolet iN10 Documentation v9.4	Fixed upgrade from OMNIC Picta 1.3 and older
Nicolet iS50 Raman for OMNIC 9 v9.8 HF1	See OMNIC for Raman iS50 Release Notes
NXR FT-Raman Accessory Module Software v9.8 HF1	See OMNIC for Raman NXR Release Notes
OMNIC 9 FTIR Universal Driver Files v9.8	Fixed upgrade from OMNIC Picta 1.3 and older
OMNIC 9 Service Pack Plug-in v4.6 HF1	Fixed upgrade from OMNIC Picta 1.3 and older
OMNIC Data Security v9.8 HF1	See OMNIC Release Notes
OMNIC for Dispersive Raman v9.8 HF1	See OMNIC for Dispersive Raman Release Notes
OMNIC MC v9.8 HF1	Fixed upgrade from OMNIC Picta 1.3 and older
OMNIC Picta Data Security v1.6 HF1	Fixed upgrade from OMNIC Picta 1.3 and older
OMNIC Picta v1.6 HF1	Fixed upgrade from OMNIC Picta 1.3 and older

OMNIC Release Notes v9.8 HF1	Updated for this release
OMNIC v9.8 HF1	See OMNIC Release Notes
Security Suite Documentation v2.0 HF2	Fixed removing old version on upgrade
Security Suite Plug-ins v2.0 HF3	See Security Suite Release Notes
Security Suite v2.0 HF3	See Security Suite Release Notes
Series Data Collect v9.8 HF1	See OMNIC Release Notes
SST PEM & TRS Experiments v9.8 HF1	See OMNIC SST Release Notes
TQ Analyst - EZ Edition v9.8	Fixed upgrade from OMNIC Picta 1.3 and older
TQ Analyst v9.8	Fixed upgrade from OMNIC Picta 1.3 and older

Added Products

Service Pack Release Notes	This document's initial release
----------------------------	---------------------------------

Deleted Products

None

OMNIC Service Pack (31-Aug-2017)

Updated Products

Security Suite v2.0 HF2	See Security Suite Release Notes
-------------------------	----------------------------------

Added Products

None

Deleted Products

None

OMNIC Service Pack (04-Aug-2017)

Updated Products

9.8 HF1 ATlμS	Update of Atlμs Release Notes only
9.8 TQ ANALYST PROFESSIONAL EDITION	See TQ Analyst Release Notes
9.8 TQ ANALYST – EZ EDITION	See TQ Analyst Release Notes
9.8 HF1 OMNIC MC	See Atlμs Release Notes
1.6 OMNICxi FOR DISPERSIVE RAMAN	See Security Suite Release Notes
1.0 TGA-IR DOCUMENTATION	Moved from standalone DVD to Program DVD
1.6 INTERNATIONALIZATION FOR OMNICxi	Changed to support OMNICxi 1.6
1.6 OMNICxi DS	Changed to support OMNICxi 1.6
9.8 TQ ANALYSTS DATA SECURITY	Changed to match current Security 2.0
SECURITY SUITE SUPPORTED PRODUCTS	Changed to support OMNICxi & TQ
2.0 SECURITY SUITE DOCUMENTATION	Changed to match current Security 2.0
9.8 TQ ANALYSTS RELEASE NOTES	See TQ Analyst Release Notes
2.0 SECURITY SUITE SERVICE PACK	See Security Suite Release Notes

Added Products

None

Deleted Products

None

OMNIC Service Pack (29-Jun-2017)

Updated Products

9.8 HF1 ATμS	See Atμs Release Notes
--------------	------------------------

Added Products

None

Deleted Products

None

OMNIC Service Pack (20-Jun-2017)

Updated Products

2.0 HF1 SECURITY SUITE PLUG-INS	See Security Suite Release Notes
---------------------------------	----------------------------------

Added Products

None

Deleted Products

None

OMNIC Service Pack (14-Jun-2017)

Updated Products

9.8 GAS APP OMNIC FOR ANTARIS IGS	Rebuilt for compatibility with OMNIC 9.8
9.8 OMNIC LIBRARY CONVERTER	Rebuilt for compatibility with OMNIC 9.8
9.8 OMNIC INTERNATIONALIZATION	Rebuilt for compatibility with OMNIC 9.8
9.8 CUSTOM OMNIC	See OMNIC Release Notes
10.0 OMNIC MACROS BASIC UTILITY	No code changes, changed package structure
3.1 THERMO SOFTWARE IQ	Better # Version digits, Auto-scan & ToolBook support
9.8 ARRAY AUTOMATION	See Array Automation Release Notes
9.8 SERIES DATA COLLECT	Rebuilt for compatibility with OMNIC 9.8
9.8 MICROVIEW	Rebuilt for compatibility with OMNIC 9.8
9.8 SST PEM & TRS EXPERIMENTS	See OMNIC SST Release Notes
9.8 ATμS	See Atμs Release Notes

9.8 OMNIC DATA SECURITY	See OMNIC Release Notes
9.8 NICOLET iS50 RAMAN FOR OMNIC 9	Rebuilt for compatibility with OMNIC 9.8
9.8 TQ ANALYST PROFESSIONAL EDITION	See TQ Analyst Release Notes
9.8 TQ ANALYST - EZ EDITION	See TQ Analyst Release Notes
9.8 OMNIC MC	Rebuilt for compatibility with OMNIC 9.8
10.0 VISUALIZATION TOOLS	Rebuilt for compatibility with OMNIC 9.8
9.8 OMNIC FOR DISPERSIVE RAMAN	See OMNIC for Dispersive Raman Release Notes
9.8 NICOLET iS5 DRIVER AND HELP	Rebuilt for compatibility with OMNIC 9.8
9.8 NICOLET iS5N DRIVER AND HELP	Rebuilt for compatibility with OMNIC 9.8
9.8 NICOLET iS10 DRIVER AND HELP	Rebuilt for compatibility with OMNIC 9.8
9.8 NICOLET iS50 DRIVER AND HELP	Rebuilt for compatibility with OMNIC 9.8
9.8 NICOLET iS50 NIR FOR OMNIC 9	Rebuilt for compatibility with OMNIC 9.8
9.8 NIC 330/370 USB DRIVER AND HELP	Rebuilt for compatibility with OMNIC 9.8
9.8 NICOLET 380 DRIVER AND HELP	Rebuilt for compatibility with OMNIC 9.8
9.8 NICOLET 4700 USB DRIVER-HELP	Rebuilt for compatibility with OMNIC 9.8
9.8 NICOLET 5700 USB DRIVER-HELP	Rebuilt for compatibility with OMNIC 9.8
9.8 NICOLET 6700 DRIVER AND HELP	Rebuilt for compatibility with OMNIC 9.8
9.8 NIC 8700 USB DRIVER AND HELP	Rebuilt for compatibility with OMNIC 9.8
9.8 NICOLET IG50 DRIVER AND HELP	Rebuilt for compatibility with OMNIC 9.8
9.8 RAMAN UNIVERSAL DRIVER FILES	Rebuilt for compatibility with OMNIC 9.8
9.8 WAFER ANALYZER SW FOR OMNIC 9	See OMNIC Release Notes
9.8 NXR FT-RAMAN ACCY MODULE SW	Rebuilt for compatibility with OMNIC 9.8
4.6 OMNIC 9 UPGRADE PLUG-IN	Updated to support Security Suite 2.0
9.8 OMNIC DEMO INSTRUMENT CONFIG	Rebuilt for compatibility with OMNIC 9.8
9.8 CHROMELEON GC-IR INTERFACE	Rebuilt for compatibility with OMNIC 9.8
9.8 SERVICE PACK-ADD-IN	Updated to support Security Suite 2.0
9.8 LINKAM STAGE SUPPORT FOR OMNIC	Rebuilt for compatibility with OMNIC 9.8
1.5 DXR RAMAN FAMILY DOCUMENTATION	Updated for compatibility with OFDR 9.8
2.2 SPECTA	See OMNIC Spectra Release Notes
2.8 VALPRO SOFTWARE	See ValPro Qualification Release Notes
2.8 VALPRO METHODS	See ValPro Qualification Methods Release Notes
4.6 THERMO SCIENTIFIC SETUP FILES	Rebuilt for compatibility with Security Suite 2.0
9.8 GAS APP FOR IGS AND IG50 SP	Rebuilt for compatibility with OMNIC 9.8
1.4 OMNICXI DATA SECURITY	Rebuilt for compatibility with Security Suite 2.0
9.8 CUSTOM OMNIC RELEASE NOTES	Release Notes for OMNIC 9.8
9.7 HF1 TQ ANALYST DATA SECURITY	Rebuilt for compatibility with Security Suite 2.0
2.2 OMNIC SPECTA DATA SECURITY	Rebuilt for compatibility with Security Suite 2.0
9.8 ANTARIS DRIVER	Rebuilt for compatibility with OMNIC 9.8
9.8 RESULT DRIVER	Rebuilt for compatibility with OMNIC 9.8
9.7 HF1 TQ ANALYST RELEASE NOTES	See TQ Analyst Release Notes
2.2 SPECTA INTERNATIONALIZATION	Rebuilt for compatibility with OMNIC Spectra 2.2
2.7 HF6 VALPRO SYS QUAL FTIR DOC	Updated ValPro Documentation

Added Products

Initial integration of OMNIC Picta into the Service Pack

1.6 OMNIC PICTA FOR OMNIC 9	See OMNIC PICTA Release Notes
-----------------------------	-------------------------------

9.8 NICOLET iN10 DRIVER AND HELP	Rebuilt for compatibility with PICTA 1.6
1.6 DIAGNOSTICS FOR OMNIC PICTA	Rebuilt for compatibility with PICTA 1.6
1.6 OMNIC PICTA INTERNATIONALIZATION	Rebuilt for compatibility with PICTA 1.6
1.6 OMNIC PICTA DATA SECURITY	Rebuilt for compatibility with Security 2.0
1.0 iN5 DTGS EXPERIMENT FILES	See OMNIC Release Notes
1.0 iN5 MCTA EXPERIMENT FILES	See OMNIC Release Notes
9.8 NICOLET iN5 MICROSCOPE DOCS	See OMNIC Release Notes
2.0 SECURITY SUITE	See Security Suite Release Notes
2.0 SECURITY SUITE LANGUAGE PACK	Updated for compatibility with Security Suite 2.0
2.0 SECURITY SUITE PLUG-INS	Updated for compatibility with Security Suite 2.0
2.0 SECURITY SUITE SUPPORTED PRODUCTS	Updated for compatibility with Security Suite 2.0
2.0 SECURITY SUITE DOCUMENTATION	Updated for compatibility with Security Suite 2.0
2.0 SECURITY SUITE SERVICE PACK JUNE 2017	Service Pack for Security Suite 2.0

Deleted Products

None

OMNIC Service Pack (20-Apr-2017)

Updated Products

1.5 HF1 OMNICXI FOR DISPERSIVE RAMAN SW

Added Products

None

Deleted Products

None

OMNIC Service Pack (10-Mar-2017)

Updated Products

1.8 THERMO SECURITY ADMINISTRATION	
9.7 HF3 MICROVIEW	
9.7 HF3 ATL μ S	
9.7 HF3 OMNIC DS	
9.7 HF3 OMNIC FOR DISPERSIVE RAMAN	
9.7 HF3 FTIR UNIVERSAL DRIVER FILES	
9.7 HF3 CHROMELEON GC-IR INTERFACE	
1.1 VIDEOLOGY USB CAMERA SUPPORT	
4.5 SERVICE PACK PLUG-IN	
3.05 SENTECH USB CAMERA SUPPORT	
4.5 THERMO SCIENTIFIC SETUP FILES	
9.7 HF3 CUSTOM OMNIC RELEASE NOTES	
9.5 HF1 TQ ANALYST DS	

9.7 HF3 ARRAY AUTOMATION	
4.2 HF1 RAMAN AUTOSAMPLER	

Added Products

None

Deleted Products

None

OMNIC Service Pack (13-Jan-2017)

Updated Products

2.7 HF3 VALPRO METHODS

Added Products

None

Deleted Products

None

OMNIC Service Pack (21-Dec-2016)

Updated Products

1.5 OMNICXI FOR DISPERSIVE RAMAN SW	
1.5 INTERNATIONALIZATION FOR OMNICXI	
9.7 HF1 OMNIC INTERNATIONALIZATION	
1.5 OMNICXI OPTICAL COMPONENTS	

Added Products

None

Deleted Products

None

OMNIC Service Pack (16-Dec-2016)

Updated Products

9.7 HF2 OMNIC 9 COMPLETE DRIVER DVD	
9.7 HF2 USB UNIV DRIVER FILES	
9.7 HF2 NICOLET iS50 DRIVER AND HELP	
9.7 HF2 NICOLET iG50 DRIVER AND HELP	
9.7 HF2 CUSTOM OMNIC RELEASE NOTES	
9.7 HF2 OMNIC EDUCATIONAL W/S 10 USER	

9.7 HF2 OMNIC FOR NIR	
9.7 HF2 CUSTOM OMNIC 9	
9.7 HF2 OMNIC FOR DISPERSIVE RAMAN	
THERMO SCIENTIFIC SETUP FILES	

Added Products

MACROS/PRO FOR OMNIC 9	
9.7 HF1 OMNIC INTERNATIONALIZATION	

Deleted Products

None	
------	--

OMNIC Service Pack (09-Nov-2016)

Updated Products

9.7 HF1 OMNIC INTERNATIONALIZATION	
9.7 HF1 CUSTOM OMNIC 9	
9.7 HF1 OMNIC FOR DISPERSIVE RAMAN	
9.7 HF1 AT μ S	
9.7 HF1 OMNIC FOR NIR	
9.7 HF1 CUSTOM OMNIC RELEASE NOTES	

Added Products

None

Deleted Products

None

OMNIC Service Pack (12-Oct-2016)

Updated Products

10.0 MACROS BASIC FOR OMNIC 9	
9.7 HF2 CUSTOM OMNIC 9	
9.7 TQ ANALYST PROFESSIONAL EDITION	
9.7 TQ ANALYST - EZ EDITION	
9.7 OMNIC FOR DISPERSIVE RAMAN	
9.7 OMNIC EDUCATIONAL W/S 10 USER	
9.7 OMNIC FOR NIR	

Added Products

None

Deleted Products

None

OMNIC Service Pack (04-Oct-2016)

Updated Products

9.7 NICOLET 4700 USB DRIVER-HELP	
9.7 NICOLET 5700 USB DRIVER-HELP	
3.02 HF3 SENTECH USB CAMERA SUPPORT	
9.7 OMNIC FOR DISPERSIVE RAMAN DRIVER	
9.7 GAS APP FOR IGS AND IG50 SP	
9.7 NICOLET IG50 DRIVER AND HELP	
11.5 TOOLBOOK REDIST	
9.7 CUSTOM OMNIC RELEASE NOTES	
9.7 NICOLET iS5N DRIVER AND HELP	

Added Products

None

Deleted Products

None

OMNIC Service Pack (12-Aug-2016)

Updated Products

9.6 OMNIC DS HF1	
9.5 TQ ANALYST PROFESSIONAL EDITION	
9.5 TQ ANALYST - EZ EDITION	
9.5 TQ ANALYST DS	
9.5 ACLS ANALYSIS FOR TQ ANALYST	

Added Products

None

Deleted Products

None

OMNIC Service Pack (27-Jun-2016)

Updated Products

2.7 VALPRO METHODS

Added Products

None

Deleted Products

None

OMNIC Service Pack (22-Jun-2016)

Updated Products

9.6 HF2 CUSTOM OMNIC 9	
9.6 HF2 CUSTOM OMNIC RELEASE NOTES	

Added Products

None

Deleted Products

None

OMNIC Service Pack (21-Jun-2016)

Updated Products

9.6 HF1 GAS APP OMNIC FOR ANTARIS IGS	
9.6 HF1 CUSTOM OMNIC 9	
9.6 HF1 SERIES DATA COLLECT	
9.6 HF1 OMNIC FOR DISPERSIVE RAMAN	
9.6 HF1 GAS APP FOR IGS AND IG50 SP	
9.6 HF1 CUSTOM OMNIC RELEASE NOTES	

Added Products

None

Deleted Products

None

OMNIC Service Pack (31-May-2016)

Updated Products

1.4 HF2 OMNICXI FOR DISPERSIVE RAMAN SW	
1.4 OMNICXI DS	

Added Products

None

Deleted Products

None

OMNIC Service Pack (06-May-2016)

Updated Products

9.6 CHROMELEON GC-IR INTERFACE	
9.6 LINKAM STAGE SUPPORT FOR OMNIC	
9.6 OMNIC FOR ANTARIS WORKSTATION	
9.6 NIC 330/370 USB DRIVER AND HELP	
9.6 NICOLET 4700 USB DRIVER-HELP	
9.6 NICOLET 5700 USB DRIVER-HELP	
2.2 IZ10 ACCESSORY HELP	
THERMO SCIENTIFIC SETUP FILES	
9.6 OMNIC FOR DISPERSIVE RAMAN DRIVER	
9.5 GAS APP FOR IGS AND IG50 SP	
9.6 NICOLET IG50 DRIVER AND HELP	
9.6 CUSTOM OMNIC RELEASE NOTES	
9.5 HF1 OMNIC DS	

Added Products

None

Deleted Products

None

OMNIC Service Pack (23-Feb-2016)

Updated Products

9.5 HF1 1.3 HF1 USB UNIV DRIVER FILES	
1.3 OMNICXI FOR DISPERSIVE RAMAN SOFTWARE	
1.3 HF1 OMNICXI 64 DRIVER	

Added Products

None

Deleted Products

None

OMNIC Service Pack (18-Feb-2016)

Updated Products

9.5 HF1 NICOLET IG50 DRIVER AND HELP	
9.5 NICOLET IG50 DOCUMENTATION	
9.5 HF1 CUSTOM OMNIC RELEASE NOTES	
9.5 HF1 NICOLET 6700 DRIVER AND HELP	
9.5 HF1 NICOLET iS10 DRIVER AND HELP	

9.5 HF1 NICOLET 380 DRIVER AND HELP	
9.5 HF1 NIC 8700 USB DRIVER AND HELP	
9.5 HF1 OMNIC 9 COMPLETE DRIVER DVD	
9.5 HF1 NICOLET iS50 DRIVER AND HELP	
9.5 HF1 NICOLET iS5 DRIVER AND HELP	
9.5 HF1 USB UNIVERSAL DRIVER FILES	
9.5 HF1 NIC 330/370 USB DRIVER AND HELP	
9.5 HF1 NICOLET 4700 USB DRIVER-HELP	
9.5 HF1 NICOLET 5700 USB DRIVER-HELP	
9.5 HF1 OMNIC FOR DISPERSIVE RAMAN DRIVER	
9.5 HF1 NICOLET iN10 DRIVER AND HELP	
9.5 HF1 OMNIC LIBRARY CONVERTER	
9.5 HF1 CUSTOM OMNIC 9	
VISUAL C++ 2012 UPDATE 4(X64) REDIST	
VISUAL C++ 2013 UPDATE 4 X64 REDIST	
VISUAL C++ 2015 UPDATE 4 X64 REDIST	
9.5 HF1 CUSTOM OMNIC RELEASE NOTES	
NICOLET iG50 DRIVER AND HELP - OMNIC 9	

Added Products

None

Deleted Products

None

OMNIC Service Pack (04-Feb-2016)

Updated Products

9.5 CUSTOM OMNIC 9	
9.5 OMNIC FOR DISPERSIVE RAMAN	
9.4 TQ ANALYST - EZ EDITION	
9.4 TQ ANALYST PROFESSIONAL EDITION	
9.5 OMNIC FOR ANTARIS WORKSTATION	
10.5 TOOLBOOK REDIST	
2.7 VALPRO METHODS	

Added Products

None

Deleted Products

None

OMNIC Service Pack (12-Jan-2016)

Updated Products

9.5 HF1 NICOLET iS50 DOCUMENTATION	
9.5 HF1 NICOLET iS50 DRIVER AND HELP	
9.5 HF1 USB UNIVERSAL DRIVER FILES	

Added Products

None

Deleted Products

None

OMNIC Service Pack (23-Nov-2015)

Updated Products

9.4 GAS APPLICATION OMNIC FOR ANTARIS IGS	
9.5 OMNIC LIBRARY CONVERTER	
9.5 OMNIC INTERNATIONALIZATION	
9.5 CUSTOM OMNIC 9	
3.1 THERMO IQ FOR OMNIC 9	
9.5 ARRAY AUTOMATION	
9.5 SERIES DATA COLLECT	
9.5 MICROVIEW	
9.5 SST PEM & TRS EXPERIMENTS SW	
9.5 AT _L μS	
9.5 OMNIC DS	
9.5 NICOLET iS50 RAMAN FOR OMNIC 9	
9.5 OMNIC MC	
9.5 VISUALIZATION TOOLS	
9.5 OMNIC FOR DISPERSIVE RAMAN	
9.5 OMNIC 9 COMPLETE DRIVER DVD	
9.5 NICOLET iS50 DRIVER AND HELP	
9.5 WAFER ANALYZER SW FOR OMNIC 9	
9.5 NXR FT-RAMAN ACCY MODULE SW	
9.5 USB UNIVERSAL DRIVER FILES	
9.5 OMNIC EDUCATIONAL W/S 10 USER	
9.5 OMNIC DEMO INSTRUMENT CONFIG	
1.1 VIDEOLOGY USB CAMERA SUPPORT	
9.5 LINKAM STAGE SUPPORT FOR OMNIC	
9.5 OMNIC FOR ANTARIS WORKSTATION	
9.5 GAS APP FOR ANTARIS IGS SP	
9.5 OMNIC FOR DISPERSIVE RAMAN DRIVER	

Added Products

Initial integration of OMNICxi into the Service Pack

1.3 OMNICXI FOR DISPERSIVE RAMAN SOFTWARE	
1.3 OMNICXI 64 DRIVER	
1.3 INTERNATIONALIZATION FOR OMNICXI	
1.3 DXR RAMAN FAMILY DOCUMENTATION	
9.4 VIS AND ANALYSIS TOOLS	
1.2 VISTOOLS RF	
VISUAL C++ 2012 UPDATE 4(X86) REDIST	
VISUAL C++ 2012 UPDATE 4(X64) REDIST	
9.2 HF2 OMNIC WORKSTATION FOR DXRXI	
OMNICXI RAMAN IMAGING SOFTWARE FOR DXRXI	
1.3 OMNICXI PRODUCT INFO	
VISUAL C++ 2013 UPDATE 4 X64 REDIST	
OMNIC 9 SERVICE PACK - 20-NOV-2015	
1.3 OMNICXI FOR DISPERSIVE RAMAN SOFTWARE	
1.3 DXR RAMAN FAMILY DOCUMENTATION	
1.3 INTERNATIONALIZATION FOR OMNICXI	
VISUAL C++ 2013 UPDATE 4 X64 REDIST	
9.4 GAS APPLICATION OMNIC FOR ANTARIS IGS	
VISUAL C++ 2013 UPDATE 4 X86 REDIST	
VISUAL C++ 2013 UPDATE 4 X64 REDIST	
1.3 OMNICXI PRODUCT INFO	
9.2 OMNIC DOCUMENTATION	
9.5 OMNIC LIBRARY CONVERTER	
9.1 OMNIC MACROS BASIC UTILITY	
OMNIC 9 ATLM μ S FOR DISPERSIVE RAMAN	
9.4 TQ ANALYST - EZ EDITION	
THERMO SOFTWARE IQ	
9.5 OMNIC FOR DISPERSIVE RAMAN	
1.2 DXR RAMAN FAMILY DOCUMENTATION	

Deleted Products

None

OMNIC Service Pack (28-Oct-2015)

Updated Products

9.3 UPGRADE XML	
9.3 SERVICE PACK UPDATE FILES	
9.1 OMNIC MACROS BASIC UTILITY	
1 SP5 OMNIC PICTA DS	

Added Products

None

Deleted Products

None

OMNIC Service Pack (08-May-2015)

Updated Products

9.3 NICOLET iN10 DOCUMENTATION	
1.0 HF4 NICOLET iS5 LS DOCUMENTATION	

Added Products

None

Deleted Products

None

OMNIC Service Pack (22-Jan-2015)

Updated Products

9.3 CUSTOM OMNIC 9	
9.2 HF5 OMNIC FOR DISPERSIVE RAMAN	

Added Products

None

Deleted Products

None

OMNIC Service Pack (10-Nov-2014)

Updated Products

2.1 OMNIC SPECTA	
9.2 HF5 OMNIC INTERNATIONALIZATION	
9.3 CUSTOM OMNIC 9	
9.2 HF5 ATLM μ S	
9.2 HF5 OMNIC MC	
9.2 HF5 VISUALIZATION TOOLS	
2 SP4 HF 5 OMNIC PICTA FOR OMNIC 9	
9.2 HF5 OMNIC FOR DISPERSIVE RAMAN	
9.2 HF5 NICOLET iN10 DRIVER AND HELP	
1 SP4 HF2 DIAGNOSTICS FOR OMNIC PICTA	
1 SP4 HF2 OMNIC PICTA INTERNATIONAL	
1 SP5 OMNIC PICTA DS	
3.02 HF2 SENTECH USB CAMERA SUPPORT	

1.1 VIDEOLOGY USB CAMERA SUPPORT	
2.0 IZ10 ACCESSORY HELP	

Added Products

None

Deleted Products

None

OMNIC Service Pack (02-Apr-2014)

Updated Products

9.1 NICOLET iS5 DOCUMENTATION	
9.2 HF4 USB UNIVERSAL DRIVER FILES	
9.2 iS5 DOCUMENTATION	

Added Products

None

Deleted Products

None

OMNIC Service Pack (28-Feb-2014)

Updated Products

9.2 HF2 CUSTOM OMNIC 9	
9.1 NICOLET iS50 RAMAN FOR OMNIC 9	
9.2 HF1 NIC NXR 9600 DRIVER AND HELP	
9.2 HF3 NICOLET iS50 DRIVER AND HELP	
9.1 NXR FT-RAMAN ACCY MODULE SW	
9.1 UPGRADE SETUP FILES	
9.2 HF1 USB UNIVERSAL DRIVER FILES	
3.02 SENTECH USB CAMERA SUPPORT	
9.2 HF1 NICOLET iS10 DRIVER AND HELP	
9.2 HF1 NICOLET iN10 DRIVER AND HELP	
9.3 HF3 NICOLET iS5 DRIVER AND HELP	

Added Products

None

Deleted Products

None

OMNIC Service Pack (09-Sep-2013)

Updated Products

9.2 HF1 TQ ANALYST PROF EDITION	
9.2 HF1 TQ ANALYST - EZ EDITION	

Added Products

None

Deleted Products

None

OMNIC Service Pack (22-Jul-2013)

Updated Products

9.2 HF1 NICOLET iS50 DOCUMENTATION

Added Products

None

Deleted Products

None

OMNIC Service Pack (15-Jun-2013)

Updated Products

9.2 NICOLET iS10 DOCUMENTATION	
3.0 OMNIC DS DOCUMENTATION	
M-P_NICOLET_iS10_DOC	
M-P_NICOLET_CENTAURUS_DOC	
M-P_TGA-IR_DOC	
M-P_SMART ORBIT_DOC	
9.2 NICOLET X700 DOCUMENTATION	

Added Products

None

Deleted Products

None

OMNIC Service Pack (01-May-2013)

Updated Products

9.1 AT μ S	
9.1 VISUALIZATION TOOLS	
1 SP4 OMNIC PICTA FOR OMNIC 9	
9.2 HF1 USB UNIVERSAL DRIVER FILES	

Added Products

None

Deleted Products

None

OMNIC Service Pack (08-Mar-2013)

Updated Products

9.2 HF1 CUSTOM OMNIC 9	
9.2 HF1 NICOLET 6700 DRIVER AND HELP	
9.2 HF1 NICOLET iS10 DRIVER AND HELP	
9.2 HF1 NICOLET 380 DRIVER AND HELP	
9.2 HF1 NIC NXR 9600 DRIVER AND HELP	
9.2 HF1 NIC 8700 USB DRIVER AND HELP	
9.2 HF1 OMNIC FOR DISPERSIVE RAMAN	
9.2 HF1 NICOLET iN10 DRIVER AND HELP	
9.2 HF1 NICOLET iS50 DRIVER AND HELP	
9.2 HF1 NICOLET iS5 DRIVER AND HELP	
9.2 HF1 NICOLET iS50 DOCUMENTATION	
9.2 HF1 USB UNIVERSAL DRIVER FILES	
9.2 HF1 OMNIC EDUCATIONAL W/S 10 USER	
9.2 HF1 OMNIC 9 FOR ANTARIS II	
9.2 HF1 OMNIC FOR ANTARIS WORKSTATION	
9.2 HF1 NIC 330/370 USB DRIVER AND HELP	

Added Products

None

Deleted Products

None

OMNIC Service Pack (04-Oct-2012)

Updated Products

9.1 SST PEM & TRS EXPERIMENTS SW	
----------------------------------	--

9.1.1 CUSTOM OMNIC 9	
9.1 USB UNIVERSAL DRIVER FILES	

Added Products

None

Deleted Products

None

OMNIC Service Pack (09-Aug-2012)

Updated Products

9.1 OMNIC DOCUMENTATION	
9 OMNIC LIBRARY CONVERTER	
9.1 MACROS BASIC FOR OMNIC 9	
9.1 OMNIC INTERNATIONALIZATION	
9.1 CUSTOM OMNIC 9	
2.2 GAML CONVERTERS PACKAGE	
9.1 OMNIC MACROS BASIC UTILITY	
2.2 THERMO IQ FOR OMNIC 9	
9.1 SERIES DATA COLLECT	
9.1 TQ ANALYST PROFESSIONAL EDITION	
9.1 VISUALIZATION TOOLS	
2.5 OMNIC DS DOCUMENTATION	
9.1 NICOLET iS50 DRIVER AND HELP	
9.1 NICOLET iS50 DOCUMENTATION	
9.1 UPGRADE SETUP FILES	
9.1 OMNIC 9 UPGRADE INDICATOR FILE	
NICOLET iS50 OMNIC NIR MODULE	

Added Products

None

Deleted Products

None

How to Contact Us

Current contact information is located at www.thermofisher.com

Select the Support > Contact Us

Trademarks

Microsoft and Windows are either trademarks or registered trademarks of Microsoft Corporation in the United States and/or other countries.