

# KNOWLEDGE SHARING

**Question/Issue/Symptom/Problem:**

How to Properly Turn Off Standard Performance Freezer

**Resolution/Answer:**

Follow Procedure Below:



1. Turn the power switch off on the back of the unit.
2. Once the switch is in the off position, the display will show "OFF" followed by "YES" and then "NO" in 2 second intervals.
3. The checkmark will also be illuminated while "YES" and "NO" are showing. Press the checkmark while "YES" is showing. With a flashing "YES", you are required to confirm by pressing the checkmark button a second time.
4. A count down will start on screen and then will go off. Power down will then be complete.

---

**Title:**

Standard Performance Freezer Power Down Procedure

**Audience:**

Public

**Author:**

Ron Prickett

---