


# Instruction Sheet

## *How to Change Battery on TSX Series ULT Freezers*




TSX, RLE, TLE, HLE, 89000
Revision Date: September 24, 2018

*How to Change battery on TSX Series ULT Freezers*

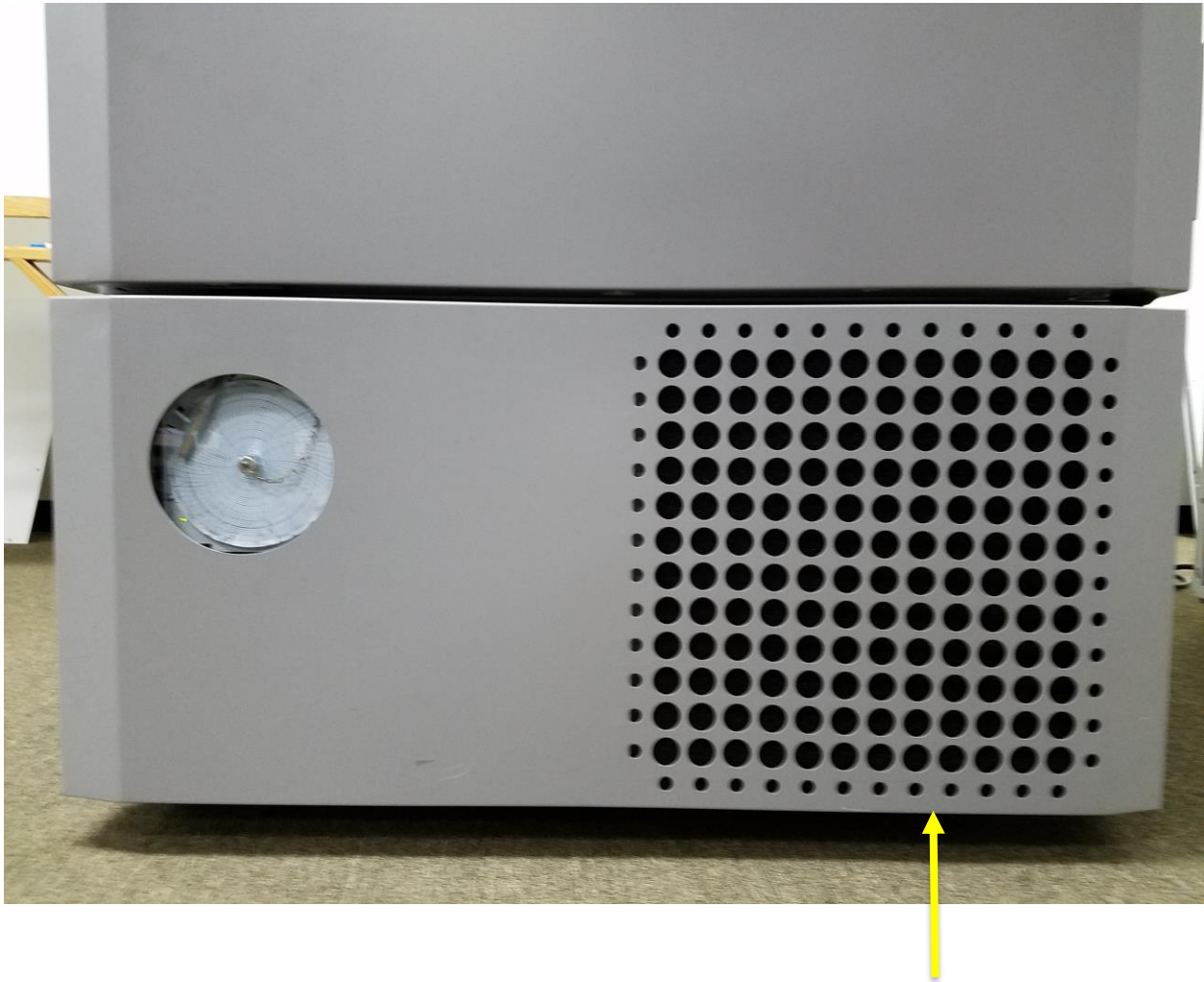

1. Open the door on the bottom of the unit.


## *How to Change battery on TSX Series ULT Freezers*

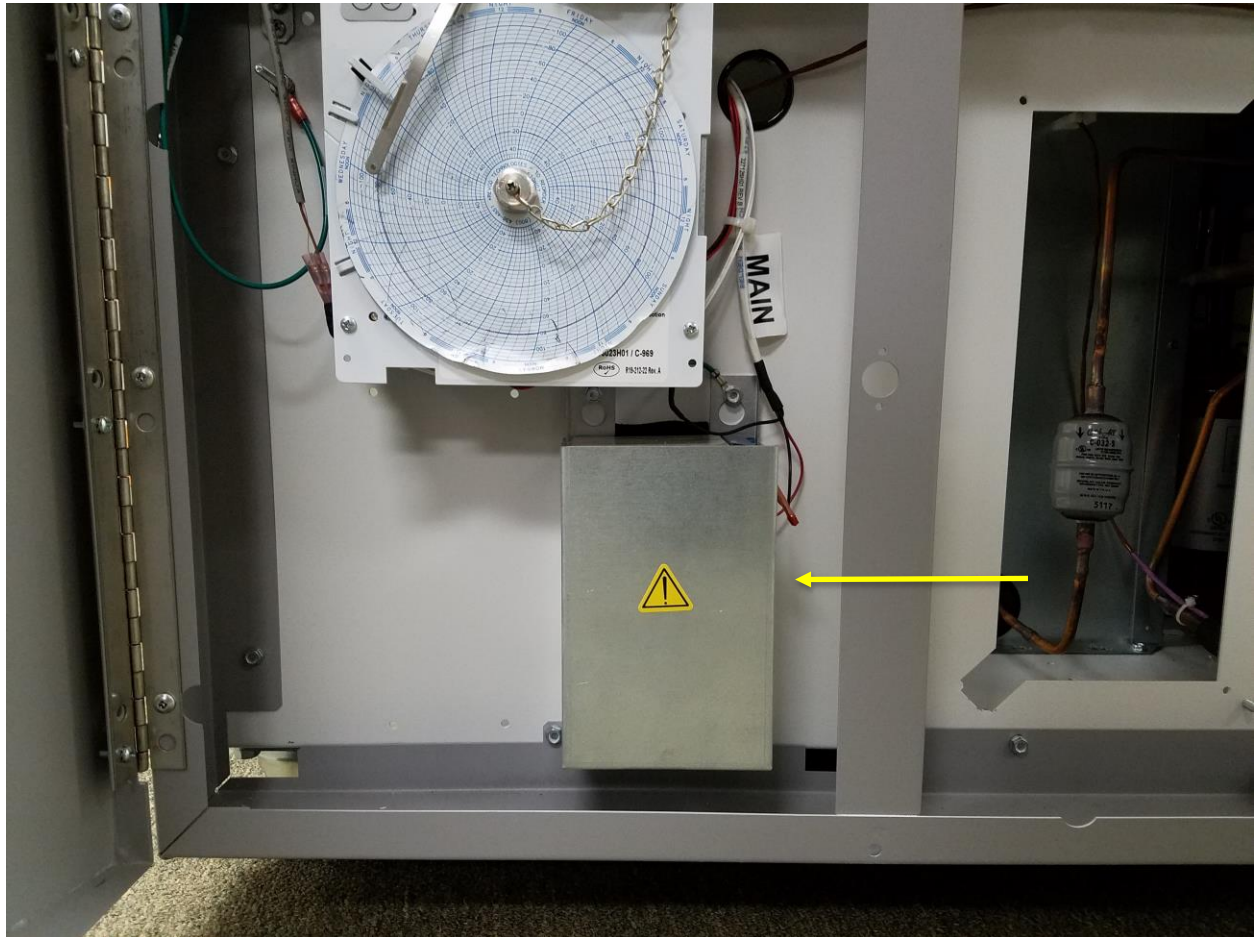

***The battery(s) are located at the front of the compressor deck, behind the bottom front panel door. The yellow arrow is pointing to bracket that holds battery in place.***

2. Label and disconnect the two red and black leads from the battery. Loosen the three bracket screws (do not remove screws) and remove the ground screw and lift bracket. Remove the expired battery and recycle or discard according to proper procedures.

*How to Change battery on TSX Series ULT Freezers*

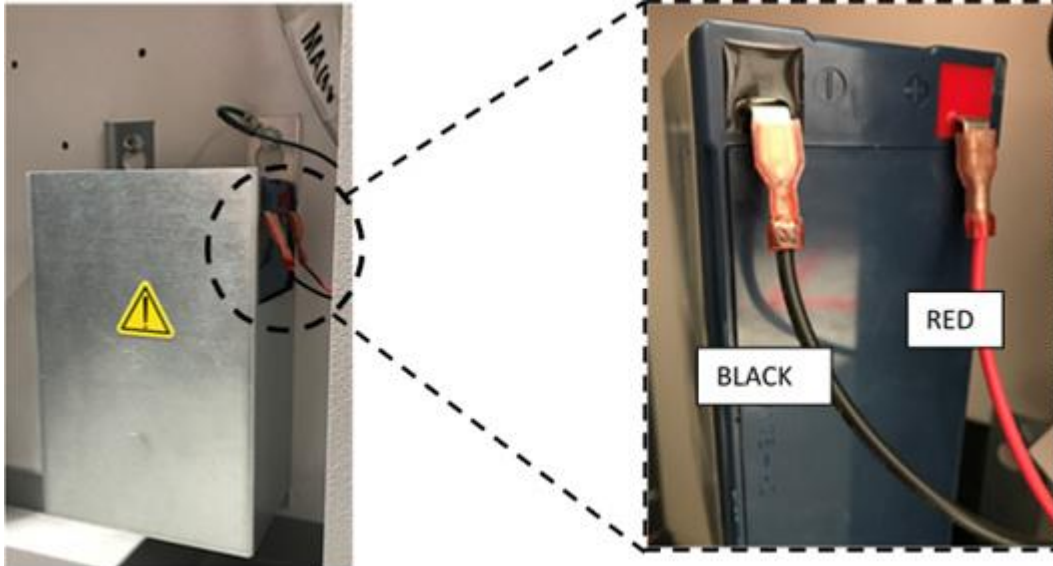


3. Install new battery (red is the positive lead and black is the negative) and tighten bracket screws.
4. Allow 4-6 days for the battery to fully charge. Monitor unit for proper operation.


*Reset Battery Clock*

Press buttons below in following order:

- Heart icon

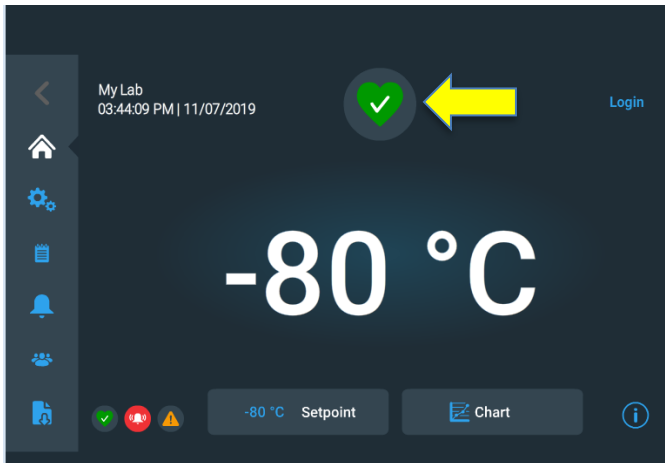

- Conditions Detail

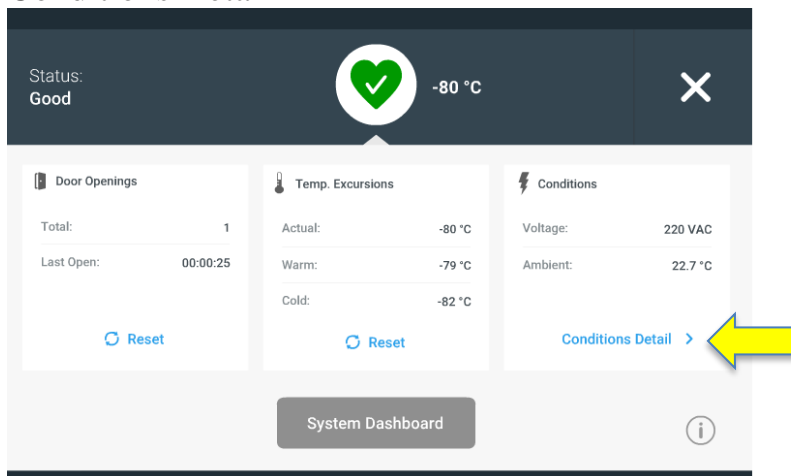

## How to Change battery on TSX Series ULT Freezers

- Reset (Once reset is pressed. The battery replace section will change to 12 months.)

