

KNOWLEDGE SHARING

Question/Issue/Symptom/Problem:


How to install sensors to unit thru 1-inch Access Port.

Resolution/Answer:

Follow Steps below:

There is a 1-inch port located on the rear lower right-hand side of High-Performance Refrigerator and Freezers +4 thru -30 units. Remove both exterior and interior plastic white caps. After this remove the batten insulation. Route independent sensors into unit. It is recommended to place by unit display sensor or as close as possible. It is important to seal up the access port after running sensors through. We suggest using a putty material - Permagem. This can be sourced locally or purchased through Thermo as part # 13015.

Note: It is recommended to thaw freezers before installation since access port insulation may be frozen.

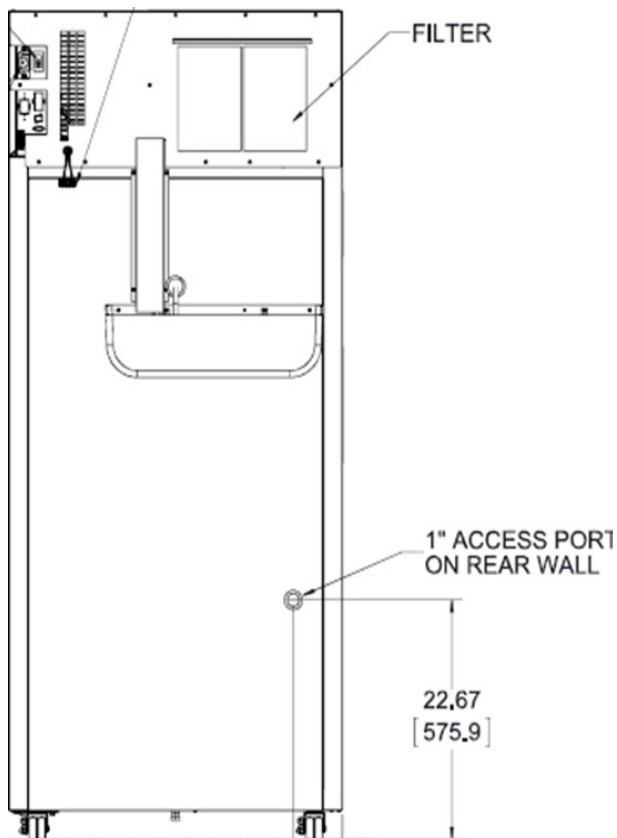

Interior plug example:



Exterior plug example

Title:

Instructions: TSX High Performance Refrigerators and Freezers - Installing Sensors thru Rear Access Port

Audience:

Public

Author:

Michael Warner
