


# KNOWLEDGE SHARING

## Question/Issue/Symptom/Problem:

How to Adjust Standard Performance Freezers Warm & Cold Alarm Setpoints.

## Resolution/Answer:

Follow Steps below:

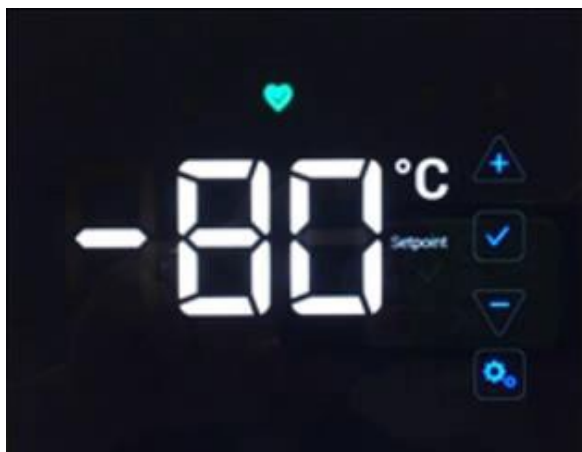

**Check Mark** – Saves a change to the selected value.



**Minus** – Decreases the value of the selected setting.



**Settings** – The settings icon represents the various settings including:

1. Press the settings button (gear icon on lower right corner).
2. Warm alarm will be displayed, use the + or – arrow buttons to adjust, setting will begin to flash.
3. Press the check mark button to save the change.
4. Press the setting button twice to view or adjust the Cold alarm setting. Use the + or – arrow to adjust, setting will begin to flash. Press the check mark button to save the change.

*Note: If set point security is active. Enter the 3 digit security code to access the settings screen. If you forget the setpoint security code, contact customer support.*

If wanting to adjust security code once in the setting screen, follow steps below.

- Press the settings button (Gear Icon) 4 times.
- The setpoint security code consists of 3 digits, each of which must be set in sequence from left to right.
- Use the plus or minus button to adjust each value, and the checkmark button to save each value of the 3- digit security code.

**Title:**

Instructions: FDE HDE RDE TDE IUE CAR VWR STP Series - How to Adjust Warm & Cold Alarm Setpoints

**Audience:**

Public

**Author:**

Michael Warner