

KNOWLEDGE SHARING


Question/Issue/Symptom/Problem:

How to Adjust Display Offset on a Carel Controller.

Resolution/Answer:

Follow Steps below:

Adjusting Display Offset (**Pharmacy Units Only**)

1. Press and hold the Program  Button for 3 seconds. You will see three dashes appear one at a time to represent the three seconds. Once they are all there you can release the program button.



2. Press the program button again to enter the parameter menu.



3. Press the up or down arrows to select "cr". Press the program button to select the parameter.



4. Press the up or down arrows to select the desired offset amount. Press the program button when done.



5. Press up or down arrows to select "ESC", and press program button. Repeat step until 3 dashes screen. At this point let the controller time out back to temperature display.



Title:

Instructions: TSG FBG CHG Series Pharmacy Refrigerators - How to Adjust Display Offset on Carel Controller

Audience:

Public

Author:

Ron Prickett

