

WX Plus User Management Reset

Direction:

1. Access to the Maintenance Screen for Service Engineer
TEMP→9→Hold→9→Hold→Enter.
2. Select “Initialization”
3. Select “User List / Schedule” to initialize the existing users list and schedules.
4. When there will be a pop-up screen saying, “Do you initialize it?”, please select “YES” to proceed.
5. After the initialization, you can see the below default user in the user list.
Username: Admin
Access level: Administrator
Password: 1111
6. Please follow the regular operation to register the new users.
Menu→Manager→User Management→Add User