

# KNOWLEDGE SHARING

## Question/Issue/Symptom/Problem:

How to Replace Battery on TSG205SA.

## Resolution/Answer:

Follow the Steps below:

1. Locate the battery compartment on the lower rear of the left side panel. See below.
2. Remove the two screws securing the pass-through port and access panel.
3. Remove the panel.
4. Remove and replace only with a CR123A or IEC16340 type battery.
5. Replace the panel and tighten the two screws.

*Note: We recommend that you keep spare batteries on hand. This battery will allow you to continue to monitor internal temperature in the event of a power outage.*



This battery will not power the refrigeration. You must ensure uninterrupted power to the refrigerator and monitoring system at all times.

**Title:**

Instructions – TSG205 Series - How to Replace Battery

**Audience:**

Public

**Author:**

Steven Thomas

---