

# KNOWLEDGE SHARING

## Question/Issue/Symptom/Problem:


Below are instructions on adjusting the temperature set point.

## Resolution/Answer:

Follow the Steps below:

1. Locate the control panel on the right front side of the unit as seen below.
2. Press the **Select/Control** button to display the **Main Menu**.
3. Use the **Scroll** buttons to move to the **Settings** option, press the **Select/Control** button to confirm/select **Settings**.
4. Scroll to the **Temp** option, press the **Select/Control** button to confirm/select **Temp**.
5. Scroll to select the **Set Point Temp** and press the **Select/Control** button to confirm.
6. **Temp Set** will display to confirm your selected **Set Point Temperature**.

## Control Panel

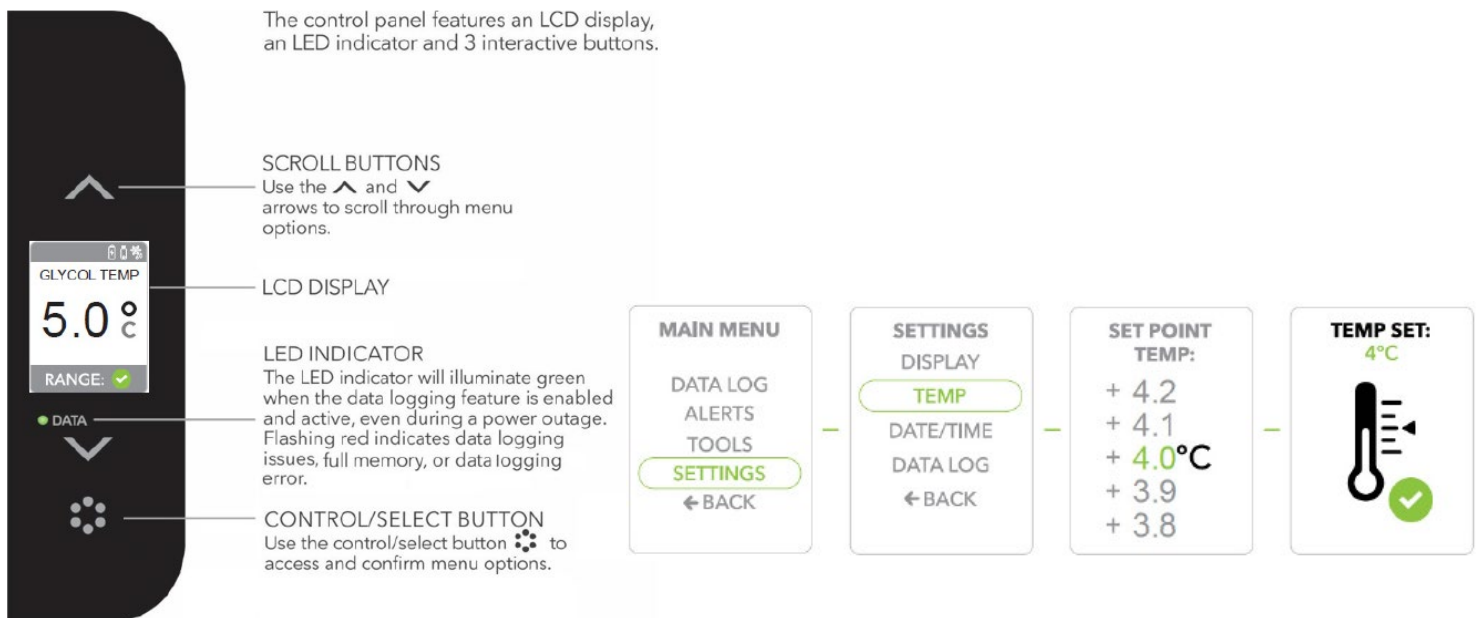

**Title:**

Instructions – TSG205 Series - How to Adjust Temperature Set Point

**Audience:**

Public

**Author:**

Steven Thomas

---