


KNOWLEDGE SHARING

Question/Issue/Symptom/Problem:

Adjusting setpoint security passcode on Standard Performance Freezers.

Resolution/Answer:

Follow the Steps below:

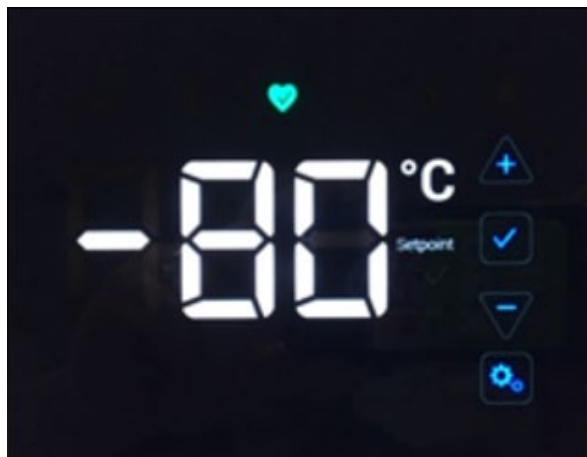

Check Mark – Saves a change to the selected value.



Minus – Decreases the value of the selected setting.



Settings – The settings icon represents the various settings including:

- If setpoint security is active. Enter 3-digit security code to access the freezer settings.
- To adjust setpoint security in the settings menu, press the settings Gear button 4 times.
- The Security code consists of 3 digits, each of which must be set in sequence.
- Use the plus and minus button to adjust each value, and the checkmark button to save each value of the 3-digit security code.



Note: If you forget the setpoint security code, contact customer support 1-866-984-3766.

Title:

Instructions – STP Freezer Series - How to Reset Setpoint Security Passcode

Audience:

Public

Author:

Steven Thomas