

## Release Notes

### Chromeleon 7 Chromatography Data System

Hotfix 111938 for SII For Empower 1.1 • Dec 2019

**Copyright**

Copyright © 2009-2019 Thermo Fisher Scientific Inc. All rights reserved.

**Trademarks**

Trademarks are property of Thermo Fisher Scientific Inc. and its subsidiaries.

Thermo Fisher Scientific Inc. provides this document to its customers with a product purchase to use in the product operation. The document is copyright protected; any reproduction of the whole or any part of this document is strictly prohibited, except with the written authorization of Thermo Fisher Scientific Inc.

This document is provided “as is”. The contents of this document are subject to being changed, without notice, in future revisions.

Thermo Fisher Scientific Inc. makes no representations that this document is complete, accurate, or error-free. Thermo Fisher Scientific Inc. assumes no responsibility and will not be liable for any errors, omissions, damage, or loss that might result from any use of this document, even if the information in the document is followed properly.

This document is not part of any sales contract between Thermo Fisher Scientific Inc. and a purchaser. This document shall in no way govern or modify any Terms and Conditions of Sale. The Terms and Conditions of Sale shall govern all conflicting information between the two documents.

**For Research Use Only. Not for use in diagnostic procedures.**

## Table of Contents

1	Introduction.....	4
2	Resolved Issues.....	5
3	Installation.....	6
3.1	Prerequisites .....	6
3.2	Installation Instructions .....	6
3.3	Compatibility.....	7
3.4	Verification.....	7

## 1 Introduction

This document describes the installation and compatibility of **Hotfix 111938 for SII For Empower**, which resolves an issue with SII For Empower 1.1.

## 2 Resolved Issues

ID	Description
<b>111938</b>	SII For Empower: Data points can be repeated if "Setup" is used prior to acquisition. In particular this can lead to problems when using Apex Tracking to evaluate data.

## 3 Installation

### 3.1 Prerequisites

Prior to installation, make sure you have the following:

- SII For Empower 1.1 installed
- Local administrative privileges
- Hotfix 111938 for SII For Empower.zip

### 3.2 Installation Instructions

1. Ensure that Chromeleon and Empower are both closed.
2. Create a folder on the Windows Desktop and unzip the contents of **Hotfix 111938 for SII For Empower.zip** into that folder.
3. Navigate to the SiiForEmpower folder found under the Empower installation e.g. **C:\Empower\Instruments\HTML\SiiForEmpower** and make a backup copy outside of the SiiForEmpower folder of the following file:
  - Thermo.SiiForEmpower.Driver.SiiInstrumentController.dll
4. Replace the Thermo.SiiForEmpower.Driver.SiiInstrumentController.dll file in the SiiForEmpower folder with the hotfix version from the unzipped folder.
5. The hotfix is now applied.
6. Reboot the PC.

### 3.3 Compatibility

Hotfix 111938 for SII For Empower is compatible with SII For Empower 1.1.

### 3.4 Verification

1. Performing a Station IQ will lead to the following error. This is expected and can be ignored by the user.



2. Navigate to the SiiForEmpower folder under the Empower installation folder:  
**C:\Empower\Instruments\HTML\SiiForEmpower**
3. Right mouse click on Thermo.SiiForEmpower.Driver.SiiInstrumentController.dll and select Properties
4. Select the Details tab of the File Properties dialog box.
5. Ensure that the File Version is shown as 1.1.0.28194:

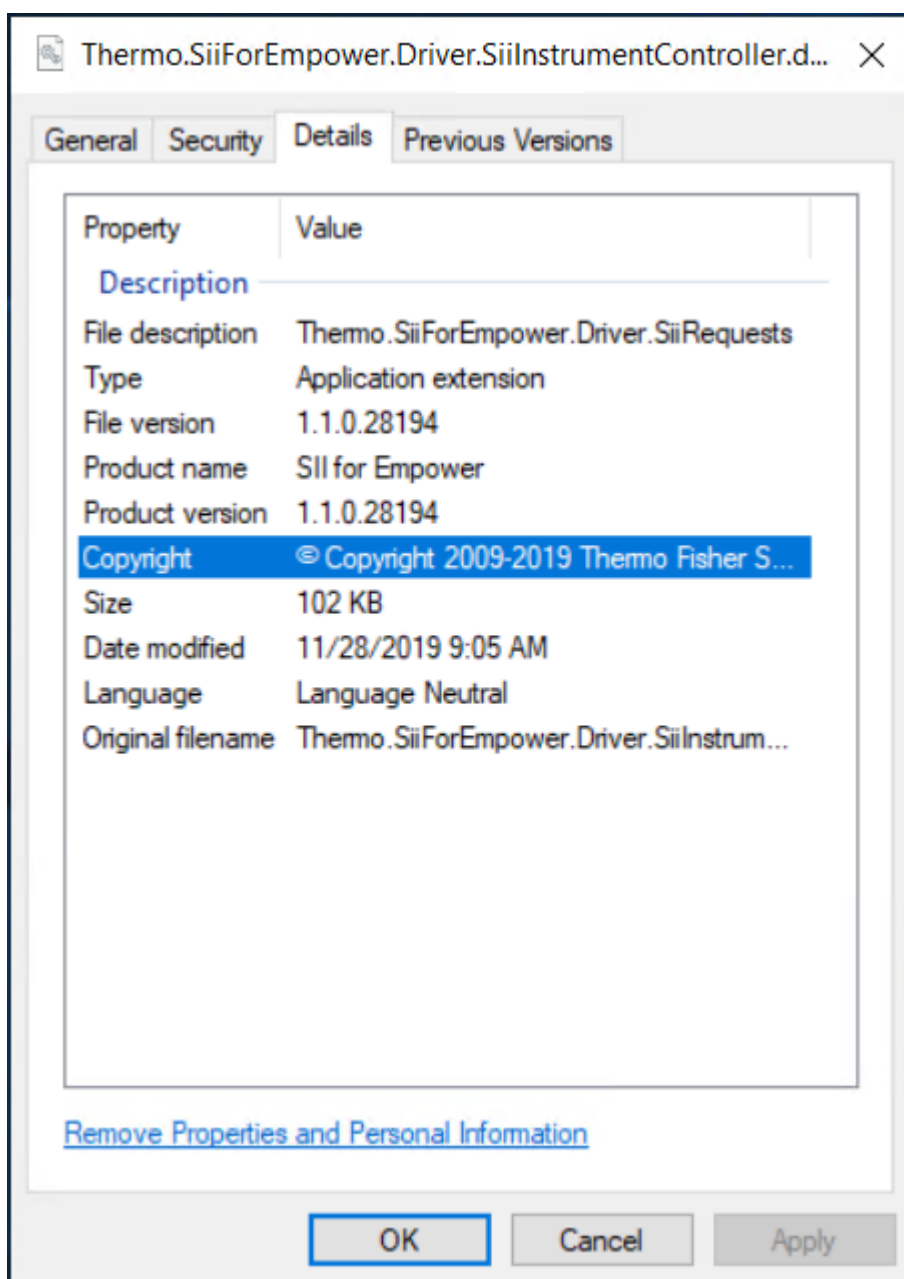


Figure 1: Thermo.SiiForEmpower.Driver.SiiInstrumentController.dll Properties dialog



[www.thermofisher.com](http://www.thermofisher.com)

---

© 2009-2019 Thermo Fisher Scientific Inc. All rights reserved.

**Thermo Fisher Scientific Inc.**  
168 Third Avenue  
Waltham, MA 02451  
USA

**ThermoFisher**  
S C I E N T I F I C